



MY11 Adapter for Pipettes calibration 11 ml



The drawings, photos and graphics used are for illustrative purposes only.

Overview

MY11 pipette calibration adapter enables calibration and/or verification of piston pipettes by means of a gravimetric method using the MYA microbalance.

Set content:

weighing pan weighing vessel evaporation ring glass chamber with an opening for dosing

Vessel capacity: 11 ml

Evaporation ring is an integral part of the adapter that increases relative humidity inside the weighing chamber. With this evaporation of the liquid during calibration/verification is stopped.

An integrated weighing pan enables centric loading of the vessel in the weighing chamber. The weighing pan and the vessel are housed in a glass chamber featuring an opening allowing the user to dispense liquid through it, and reducing air drafts.

Adapters for Pipettes Calibration

Pipette calibration adapters are intended for calibration and verification of piston pipettes using gravimetric

method. Use of the pipette calibration adapter reduces maximum capacity of the balance by the mass of the weighing vessel. The table below, which presents the minimum requirements in accordance with ISO 8655-6:2022 standard, enables to select suitable adapter for pipette calibration and respective balance version. The parameters of the balance have been selected so that the expanded uncertainty of measurement in use is less than ¼ of the maximum permissible error of the test object.

Nominal volume of apparatus under test
V

Resolution
(d) mg

Repeatability
(s)amg

Recommended balance Microbalances Analytical balances $0,5 \mu\text{l} \leq V < 20 \mu\text{l}$

0,001c 0,01d

0,006c, e 0,03d

MYA 21.5Y XA 6/21.5Y.M.A

XA 6/21.5Y.M XA 21.5Y.M.A

XA 21.5Y.M XA 21/52.5Y.M.A

XA 21/52.5Y.M XA 53.5Y.M.A

XA 53.5Y.M $20 \mu\text{l} \leq V < 200 \mu\text{l}$ 0,01 0,025 XA 52.5Y XA 110.5Y XA 82/220.5Y.A

XA 82/220.5Y XA 120/250.5Y.A

XA 120/250.5Y $200 \mu\text{l} \leq V \leq 10 \text{ ml}$ 0,1 0,2 Adapter for pipette calibration MY 11 XA 11 XA 11 XA 11 XA 11 XA

17 XA 17 XA 17 XA 100 XA 17 XA 100 Maximum vessel capacity 11 ml 11 ml 11 ml 11 ml 11 ml 17 ml 17 ml

17 ml 100 ml 17 ml 100 ml Weighing vessel mass ~ 10 g ~ 10 g ~ 10 g ~ 10 g ~ 10 g ~ 11 g ~ 11 g ~ 11 g ~

65 g ~ 11 g ~ 65 g

a The repeatability values, in this table, apply in the volumen determination of a single channel pipette. When a single-channel balance is used for volumen determination of single channel pipettes the repeatability values are double the values of this table.

c Single-channel balance.

d Multichannel balance, only valid for multi-channel pipettes. Multi-channel balances of 0,01 mg readability may be uses to test multi-channel pipettes with nominla volumes below $20 \mu\text{l}$ only if the expanded uncertainty in use is less than ¼ of the maximum permissible systematic error of the test object.

e For single-channel pipettes of nominal volumes of less than $2 \mu\text{l}$, a balance with repeatability better than the values in the table shall be used so that the expanded uncertainty in use is less than ¼ of the maximum permissible systematic error of the test object.


table {font-size: 12px} td {padding:10px; } tr {text-align:center;} @media (max-width: 800px) { table {font-size: 9px} td {padding: 1px 1px} }

Compatible with


MYA 5Y Microbalance

• MYA 21.5Y Microbalance [WL-109-0010]


Device dimensions

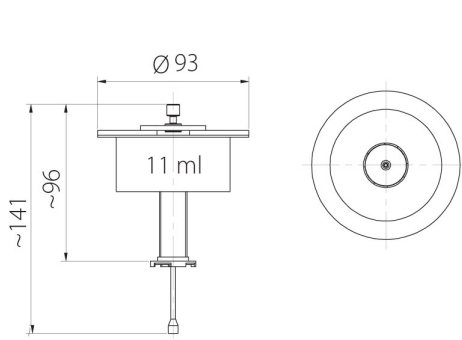


MY11

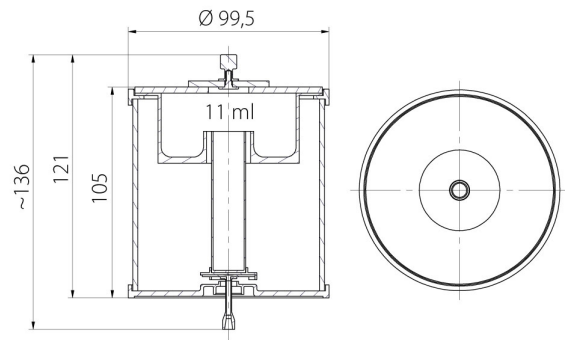

XA11

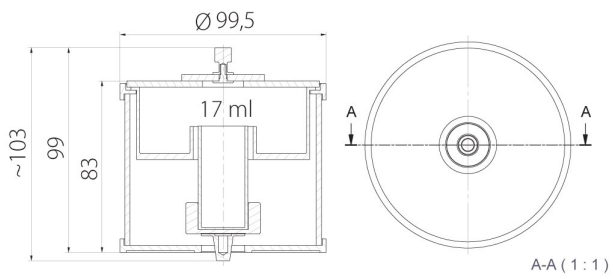

XA17

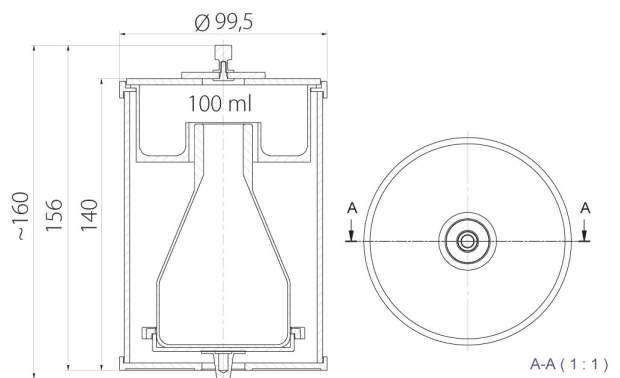

XA100