IMMB-12-02-05-22-EN-IT

Start up Guide

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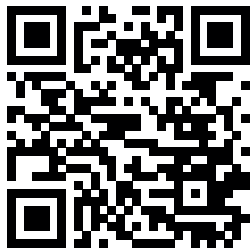
Moisture Analyzer

Termobilancia

- MA R

To view a complete User Manual go to the website or scan QR code:

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

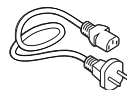
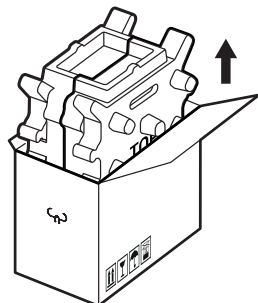
2

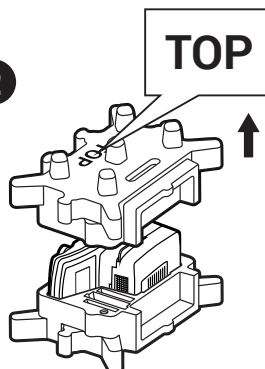3

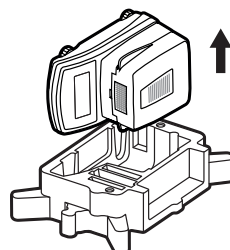
Keep the packaging should a warranty claim or service be required.

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3. RECOMMENDATIONS AND WARNINGS

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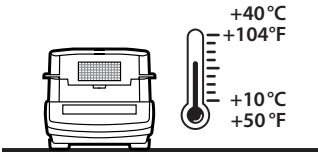
Operate the device in a room where the temperature ranges between 10–40 °C (50–104 °F) and where the relative humidity is below 80%.

Place the moisture analyzer on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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
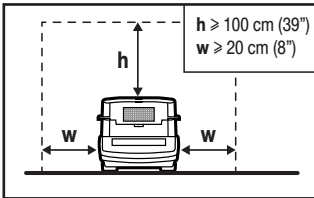
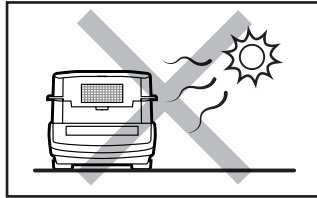
Place the device away from heat sources. Avoid exposing the device to the sunlight.

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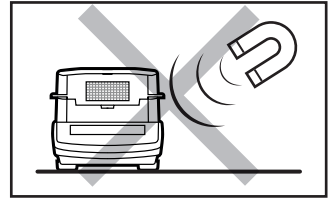
Avoid exposing the device to a magnetic field. Do not weigh magnetic substances.

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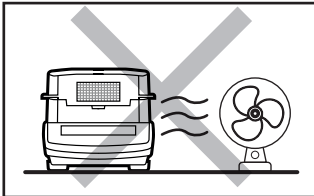
Avoid air drafts and air movements at the workstation.

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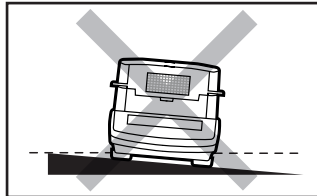
Make sure that the device is placed on an even surface.

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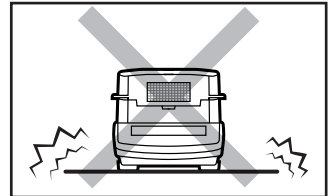
Do not place the device on an unstable surface exposed to shocks and vibrations.

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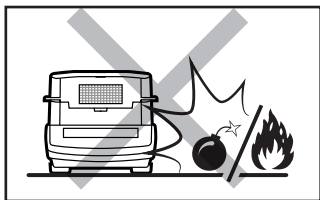
Do not place the device in a hazardous area. Do not weigh explosive or flammable materials.

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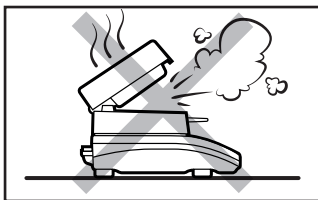Do not open in the course of operation since the heating components reach very high temperature.

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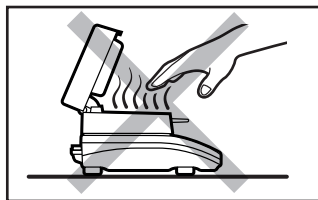Be extremely careful while taking the sample out. The sample and the moisture analyzer elements may be hot.

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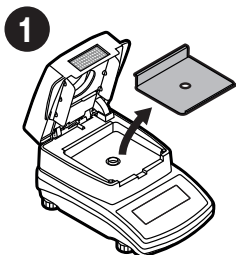
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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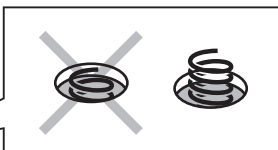
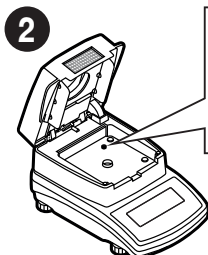
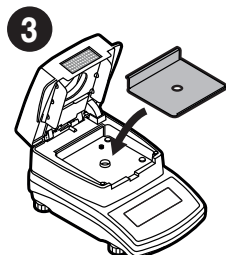
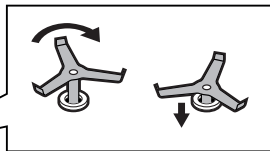
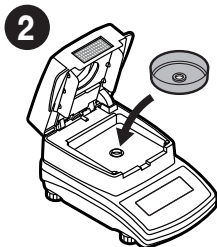
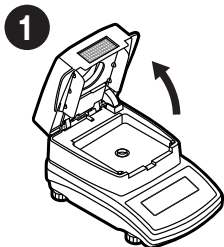
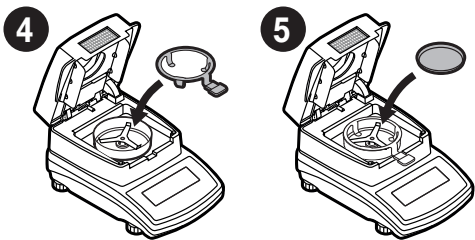
6. PREPARING FOR WORK

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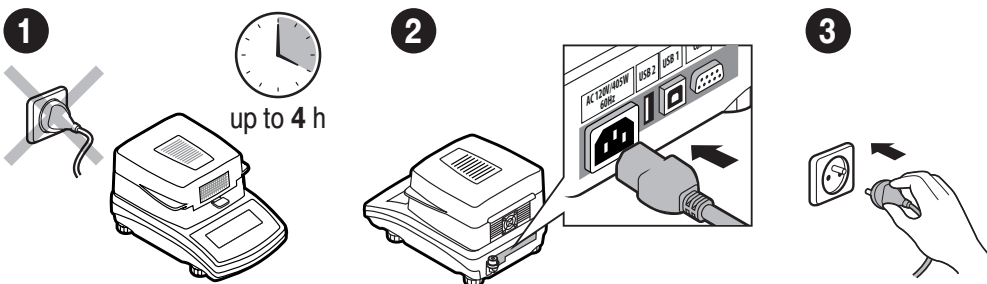
Place the moisture analyzer at the workstation, wait until it reaches room temperature.

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Prior to switching the device on let it reach room temperature (estimated stabilisation time: up to 4 hours).

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Connect the power cord to the AC socket located at the back of the device.

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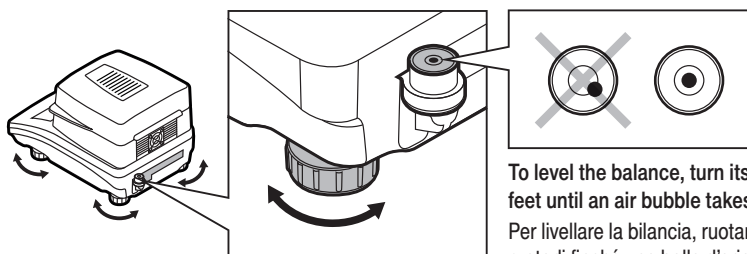
6.2. Levelling

Livellamento


To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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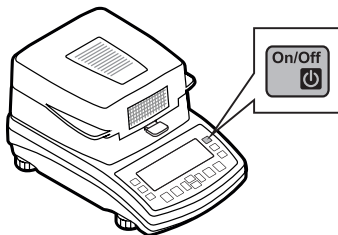Before using WEIGHING mode, carry out instrument adjustment: refer to section 8.1. (adjustment is not required for DRYING mode).

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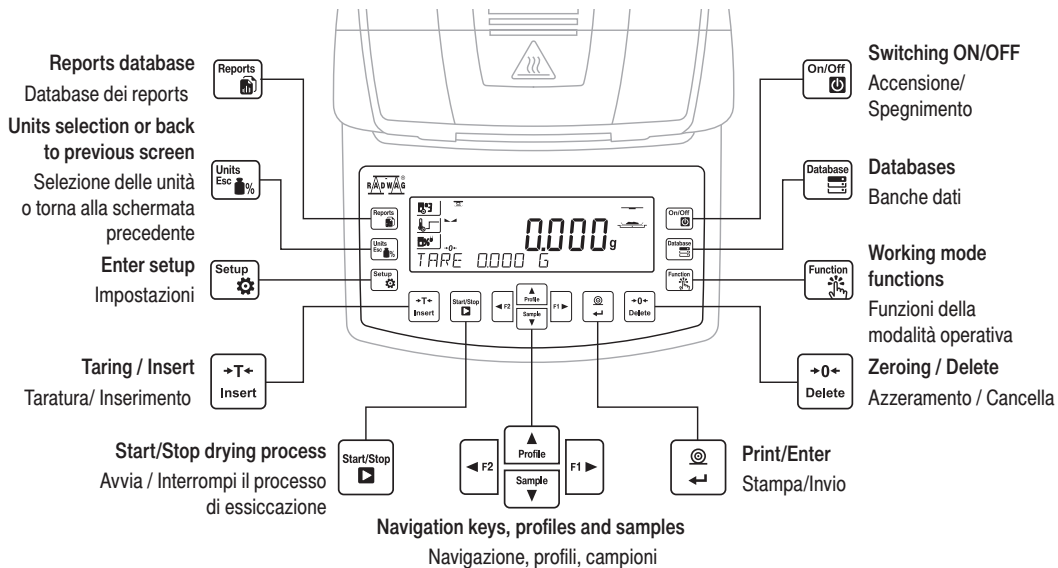
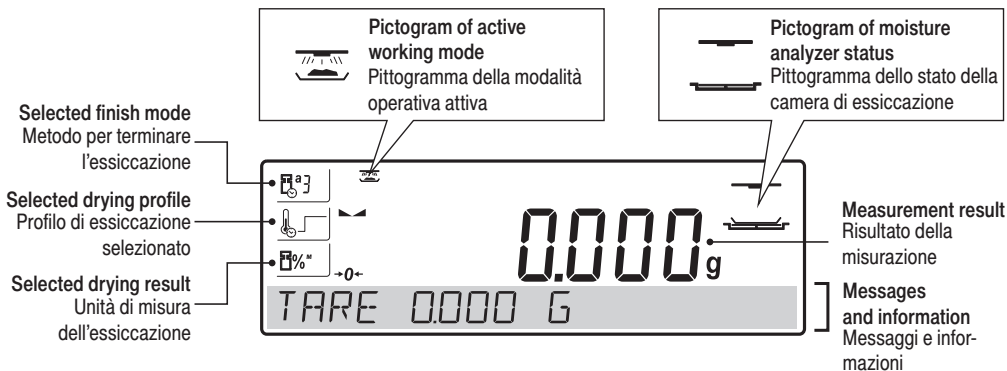
8.1. Adjustment

Regolazione

It is recommended to carry out adjustment operation prior to running the WEIGHING mode (adjustment is not required for DRYING). External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on device type and capacity.

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
The adjustment mass standard does not come with the moisture analyzer. Prior to adjustment prepare the right mass standard. You are recommended to use mass standard of class F1 or F2. View the table and select the mass standard you need. CAUTION! Make sure that the disposable weighing pan has been installed (read point 5.5.). Do not close the drying chamber in the course of adjustment.

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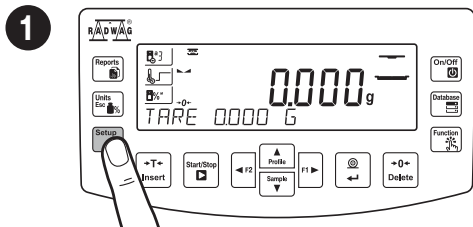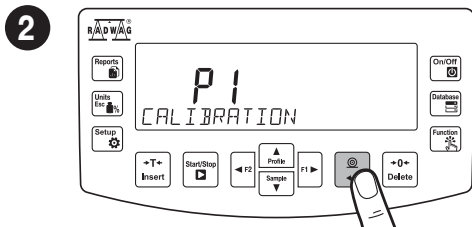
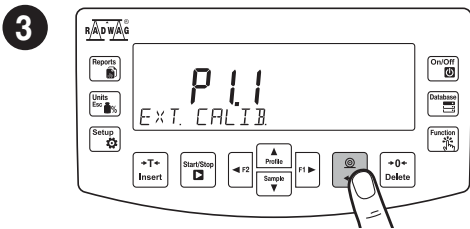
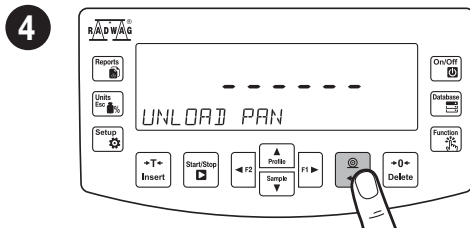
To proceed to calibration, press  key.

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
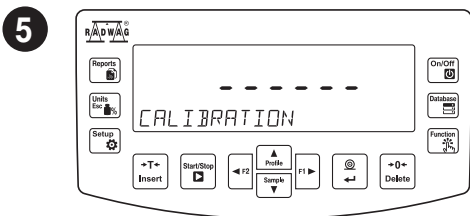
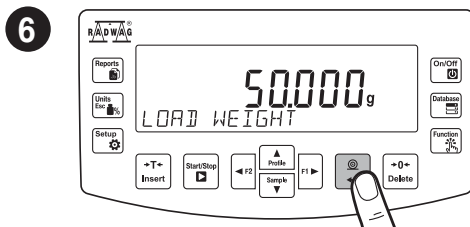
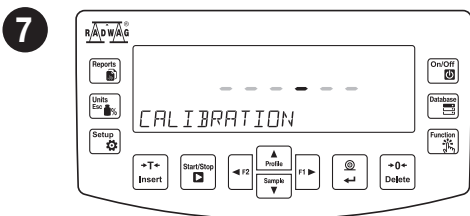
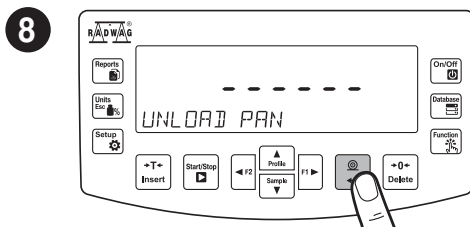
If "RANGE EXCEEDED" message (load too heavy) is displayed, remove the load and restart the adjustment process.

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
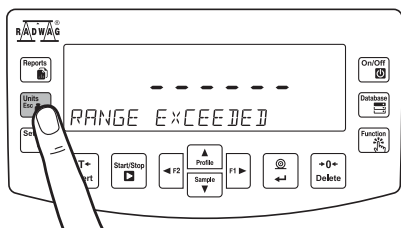
8.2. Zeroing

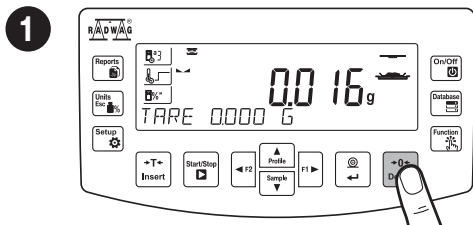Azzeramento

Make sure that the weighing pan is empty and press **+0-** key.

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
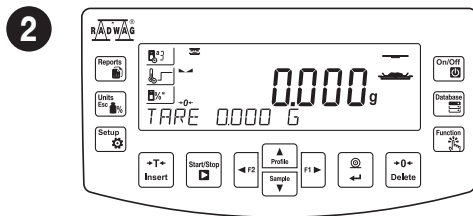
8.3. Taring

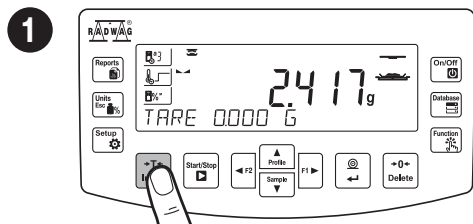Taratura

With loaded weighing pan: upon stabilisation of the weighing result, press **+T-** button.

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
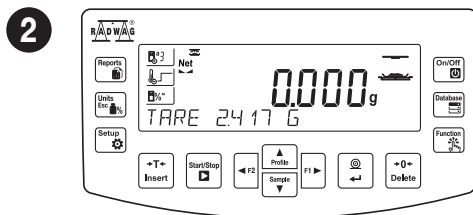
9.1. Settings for “Standard”, “Fast” and “Mild” profiles

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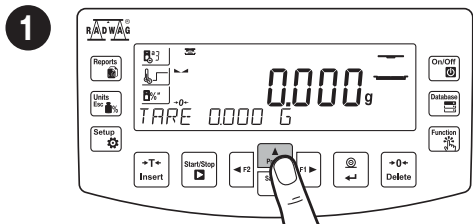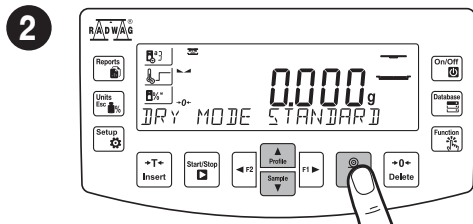
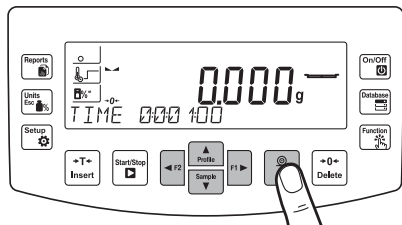
3

Using keys set drying time and press key to confirm.

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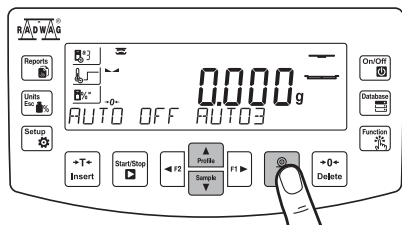
5

Using keys select drying switch-off option and press key to confirm.

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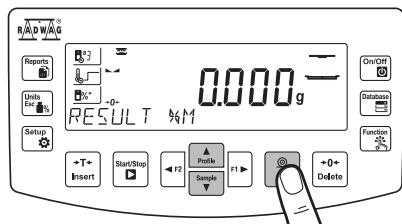
6

Using keys select measurement unit and press key to confirm.

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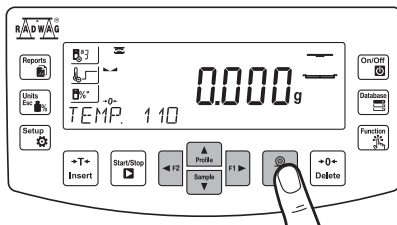4

Using keys select drying temperature and press key to confirm.

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
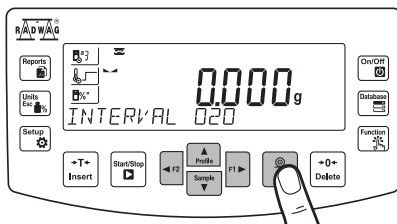
7

Using keys set interval for printouts of drying process results and press key to confirm.

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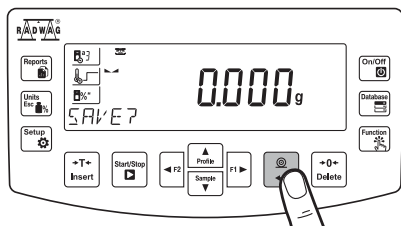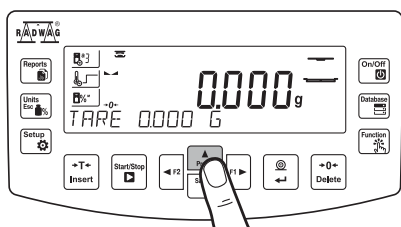
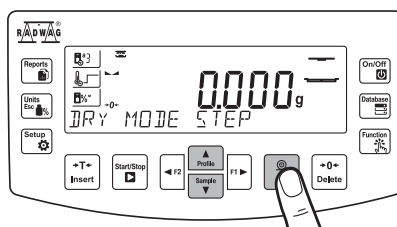
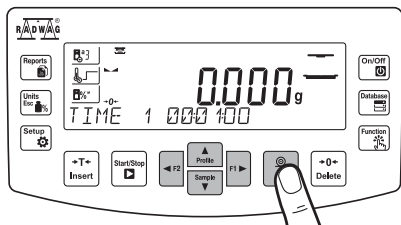
2

Using   keys select STEP drying profile and press  key to confirm.

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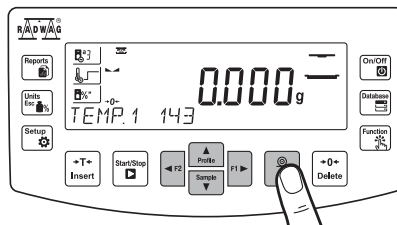
3

Using    keys set heating time for step 1 and press  key to confirm.

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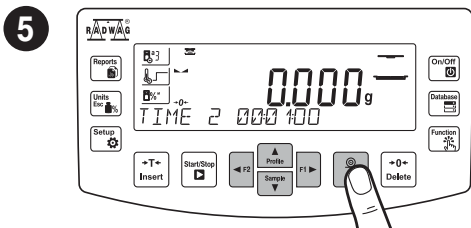
4


Using    keys set heating temperature for step 1 and press  key to confirm

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
Using keys set heating time for step 2 and press key to confirm.

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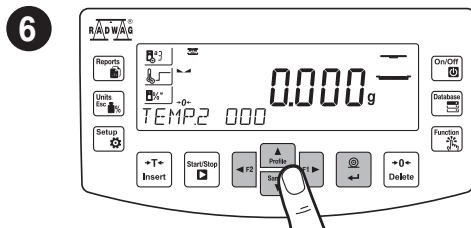
Using keys set heating temperature for step 2 and press key to confirm.

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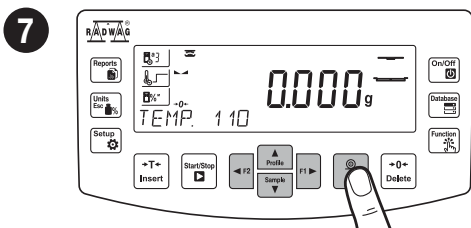
Using keys set drying temperature and press key to confirm.

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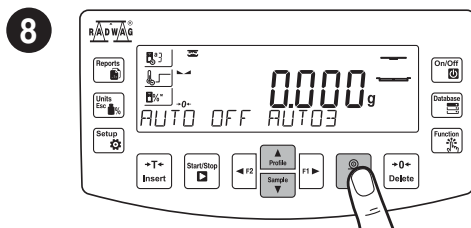
Using keys select finish mode option and press key to confirm.

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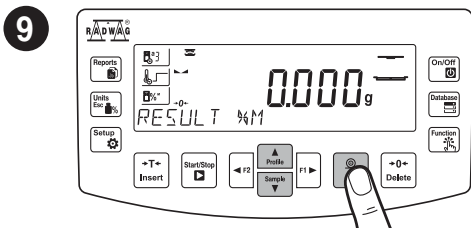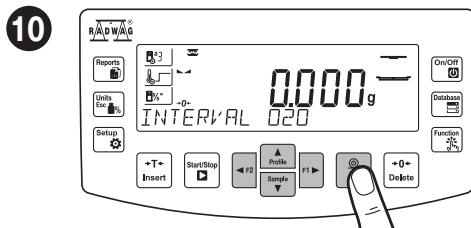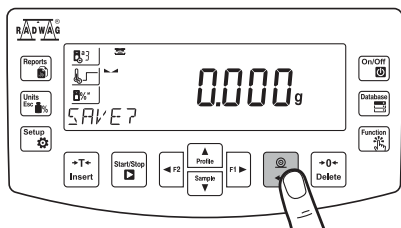
10. MOISTURE ANALYZER SETTINGS

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Some balance settings are accessible for Administrator exclusively.

Prior to balance parameters setup, log in as Administrator.

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
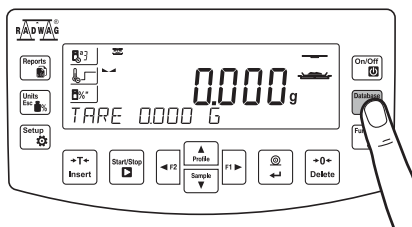
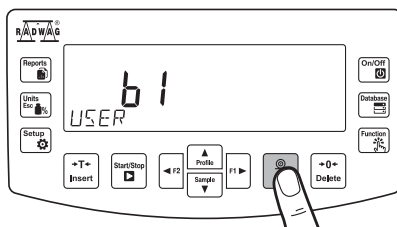
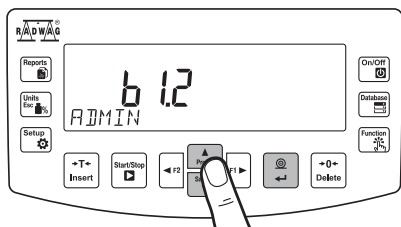
3

Using  keys select "Admin" and press  key to confirm.

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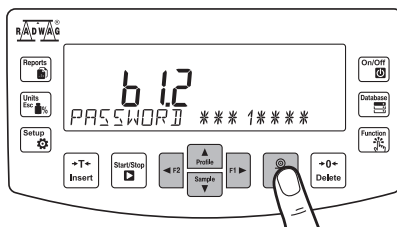
4

Using  keys select "1111" and press  key to confirm.

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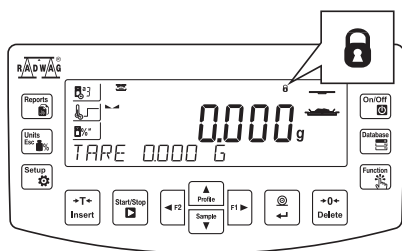10.2. User settings

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Prior to balance users setup, log in as Administrator (point 10.1).

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
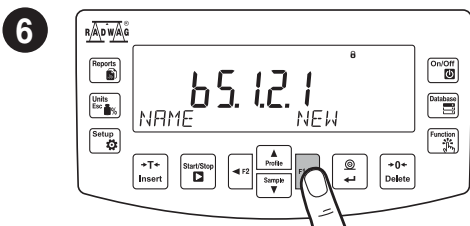
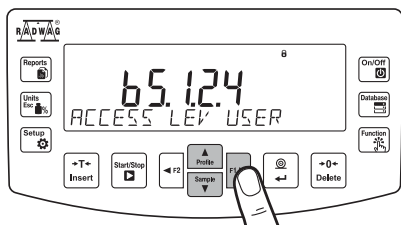
12

Using keys select “Access Level” and press key to confirm.

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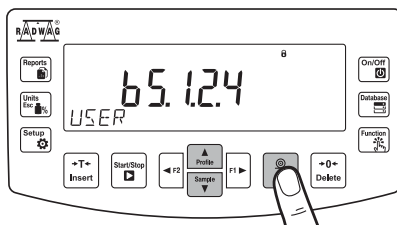
13

Using keys select permissions level and press key to confirm.

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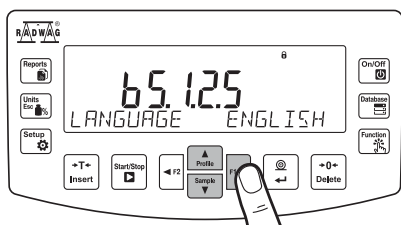
14

Using keys select “Language” and press key to confirm.

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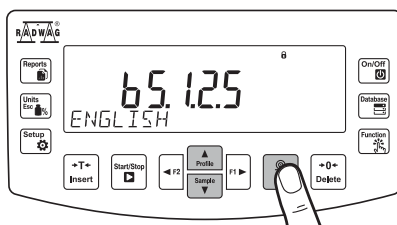15

Using keys select menu language and press key to confirm.

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11. SAMPLES AND DRYING METHODS

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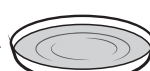
11.1. Preparing samples

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Optimal sample structure and size are crucial features influencing measurements speed, accuracy and repeatability.

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Solids:

- sample crushing necessary,
- sample of low weight required,
- large sample surface.

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- ampia superficie del campione.

11.2. Handling hazardous samples

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Some sample types require particular caution!

They may pose risk for people and objects. It is the moisture analyzer operator who is responsible for potential loss being a result of such samples use.

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Substances that in the course of heating may emit vapours causing corrosion (e.g. acids).

Apply small amount of samples, this helps to prevent risk of settling of vapour over cool housing components.

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Explosive and flammable materials and substances that in the course of heating emit flammable vapours.

Apply small amount of samples and lower drying temperature, this helps to prevent flames from rising and explosion. Use glasses and protective gloves.

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Toxic and corrosive substances and substances that in the course of heating emit toxic vapours.

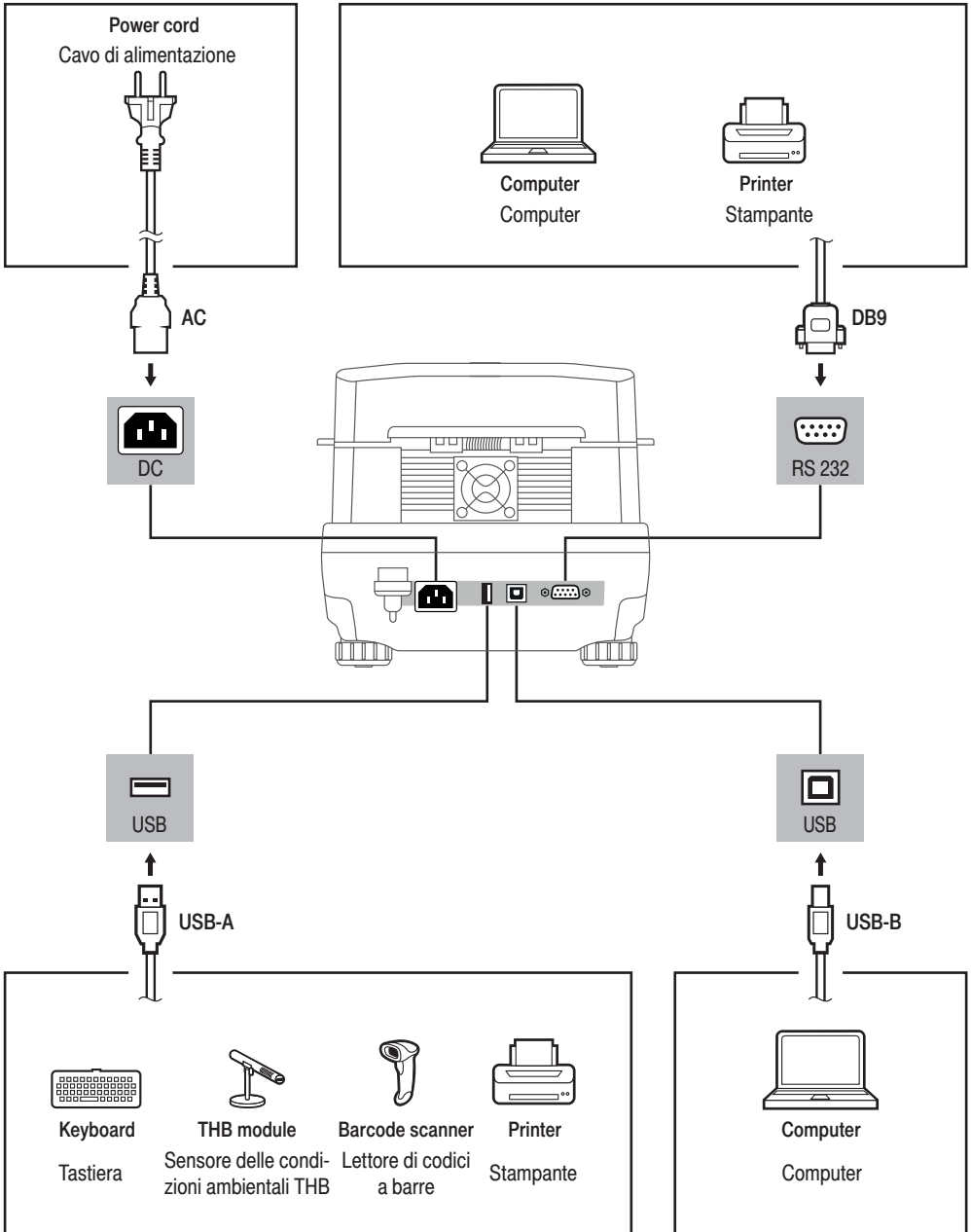
Due to risk of eyes, skin and lungs irritation, and due to risk of death, such samples must be dried in fume cupboard.

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
PULIZIA DEL DISPOSITIVO


13.1. Guidelines and precautions

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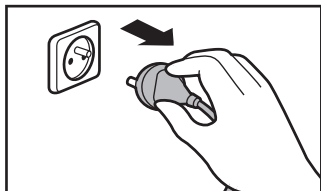
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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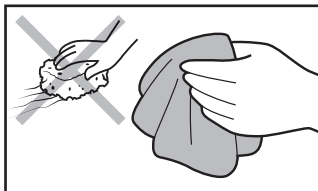Use soft cloth made of either natural or synthetic fiber exclusively. Avoid using cloth made of stiff/hard fabric, this could cause scratches on the surface!

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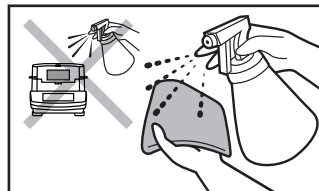Apply the cleanser onto the cloth. Mind not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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