IMMB-21-03-01-26-EN-PL

Start up Guide

Szybki start

Analytical and Precision Balances

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Analytical Balances:

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- AS X7

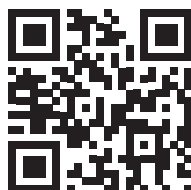
Precision Balances:

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- PS X7
- PS X7.M
- WLC X7

To view a complete User Manual go to the website or scan QR code:

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
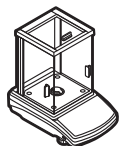
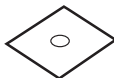
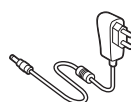
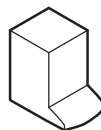
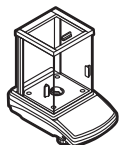
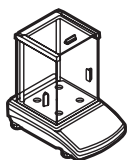
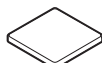
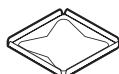
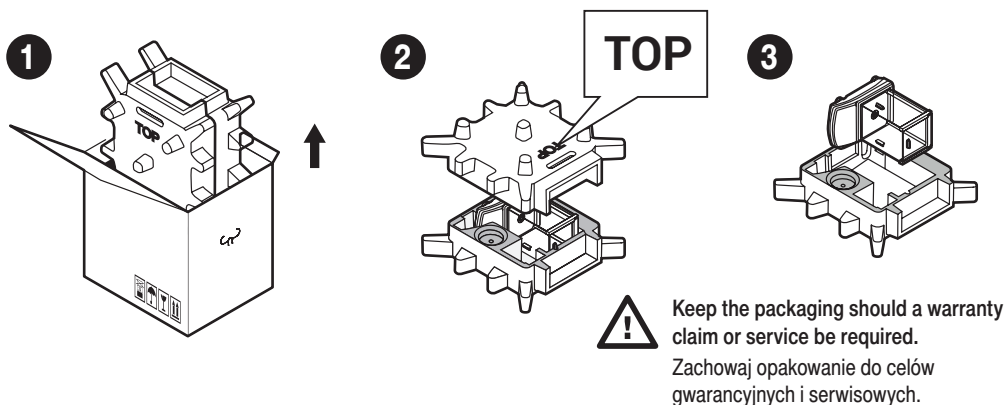
2. UNPACKING

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3. RECOMMENDATIONS AND WARNINGS

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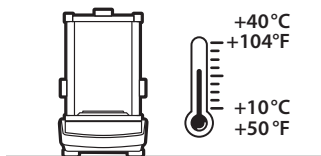
Operate the device in a room where the temperature ranges between 10–40 °C (50–104 °F) and where the relative humidity is below 80 %.

Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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
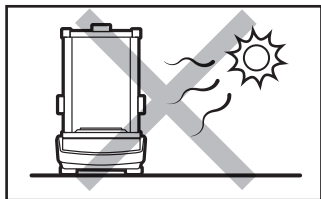
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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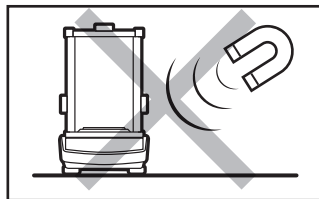
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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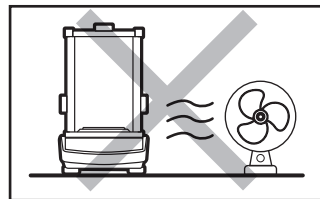
Avoid air drafts and air movements at the workstation.

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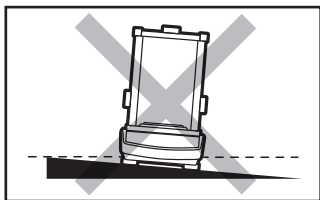
Make sure that the balance is placed on an even surface.

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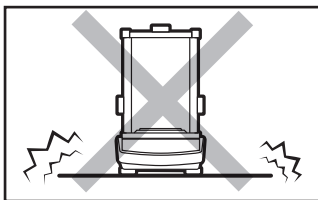Do not place the balance on an unstable ground exposed to shocks and vibrations.

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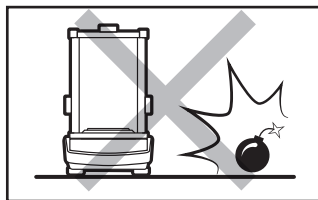Do not place the balance in a hazardous area. Do not weigh explosive materials.

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

4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Transport lock removal

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

AS X7 (d = 0,01 mg) | AS X7 (d = 0,1 mg)

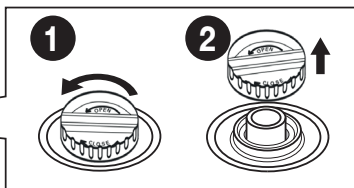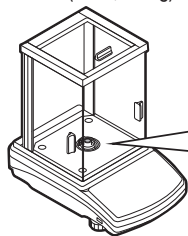Keep the transport lock should a warranty claim or service be required.

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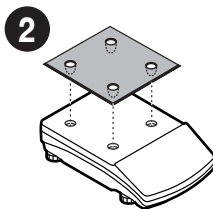
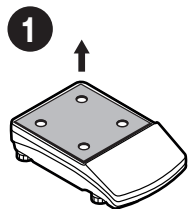
4.2. Grounding spring check

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AS X7 (d = 0,01 mg) | AS X7 (d = 0,1 mg) | WLC X7 (d = 1 mg)

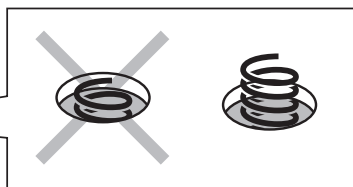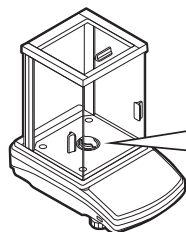Check whether the grounding spring takes its intended location.

Make sure that the grounding spring juts slightly out of the hole.

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
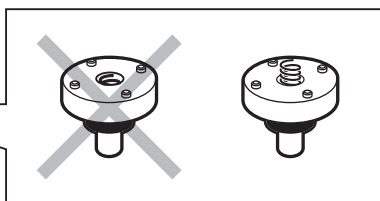
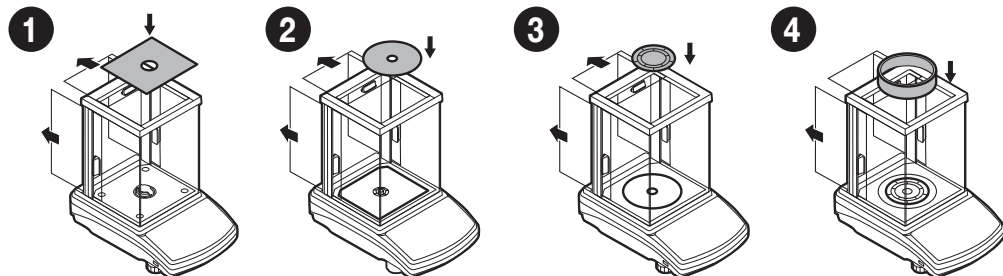
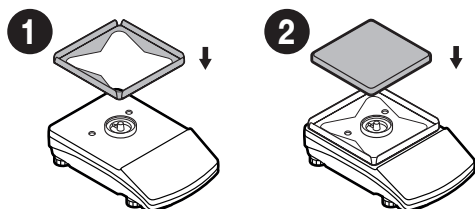
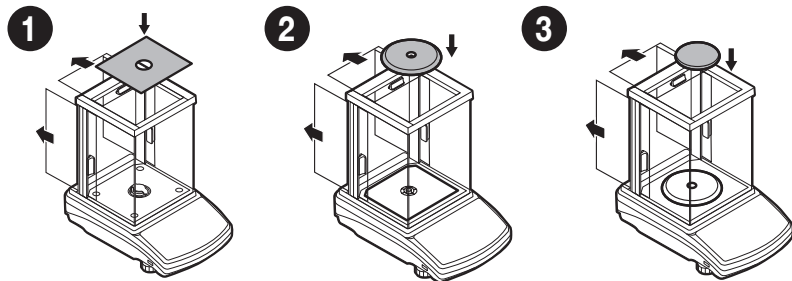
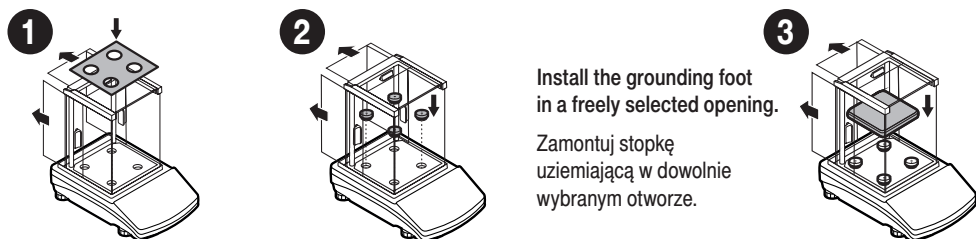
6. PREPARING FOR WORK

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6.1. Temperature stabilisation

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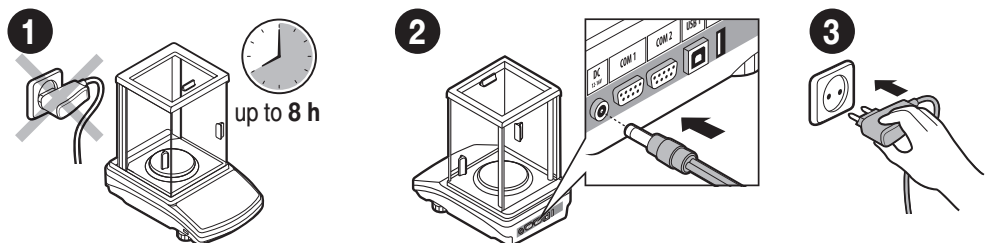Prior to switching the device on, let it reach room temperature (estimated stabilisation time: up to 8 hours).

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Connect the power adapter to DC connector.

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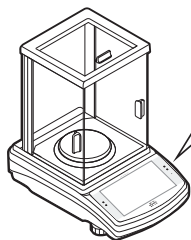
Connect the power adapter to the mains.

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6.2. Switching the balance on

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To turn on the device, double-tap the screen.

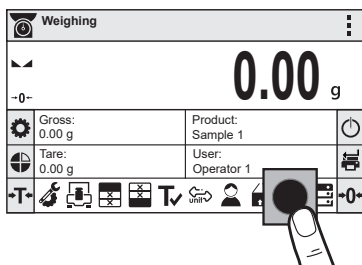
To switch the device off, press OFF key.

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
6.3. Levelling

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
Press leveling icon.

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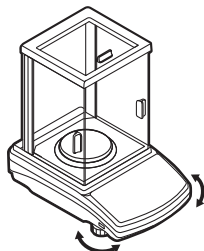
To level the balance turn its feet, keep turning the feet until the level indicator takes central position.

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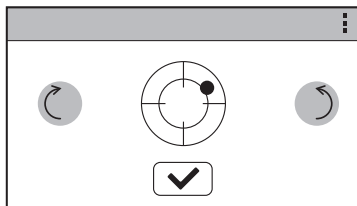Check the balance level state. The arrows shows how to move the balance feet to level the balance.

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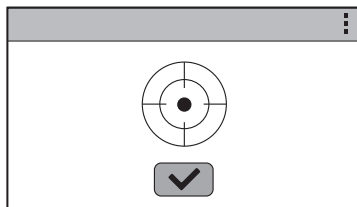Correct leveling screen. Press  button to end the operation.

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

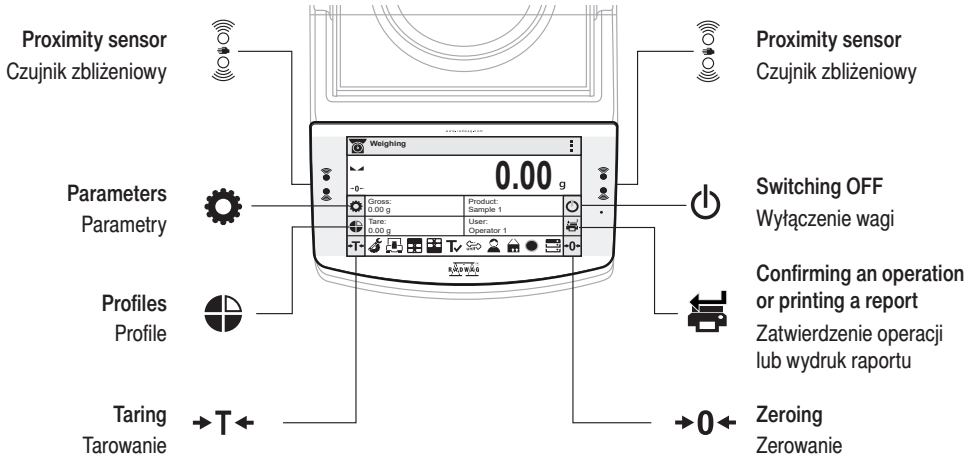
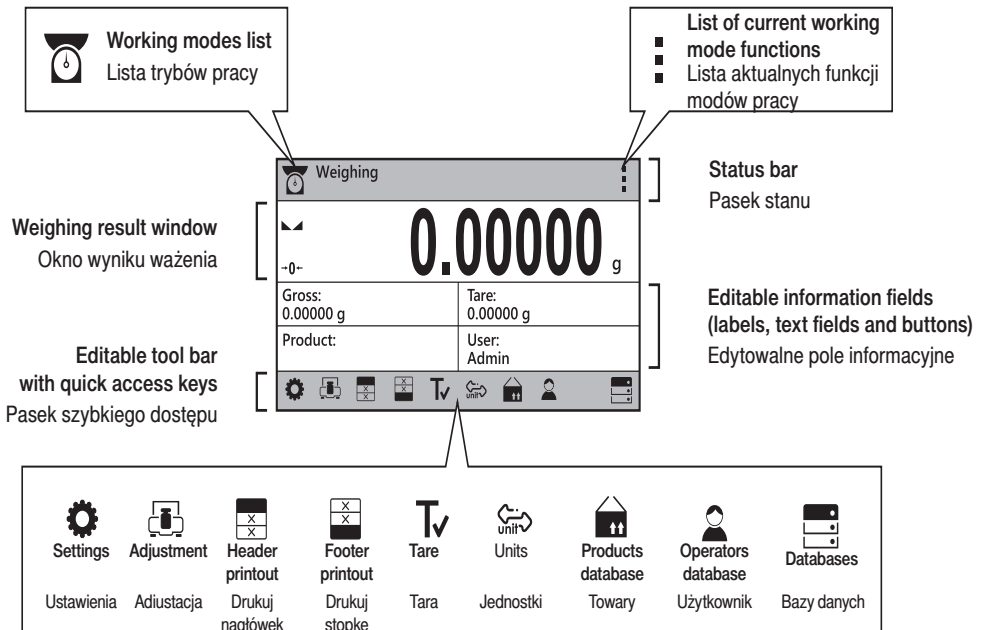
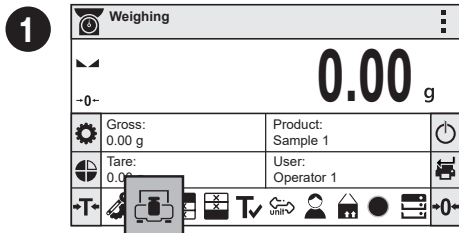
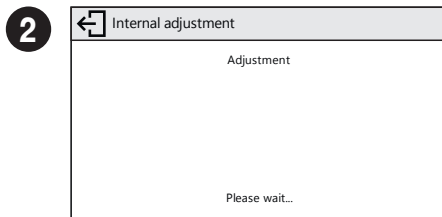
8.2. Zeroing

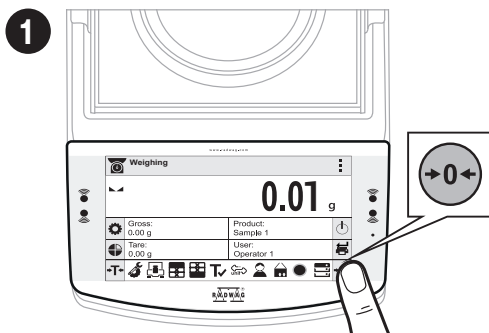Zerowanie

Make sure that the weighing pan is empty and press “Zeroing” button.

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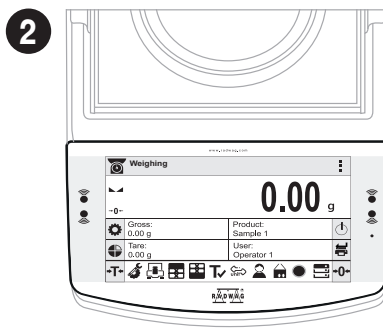
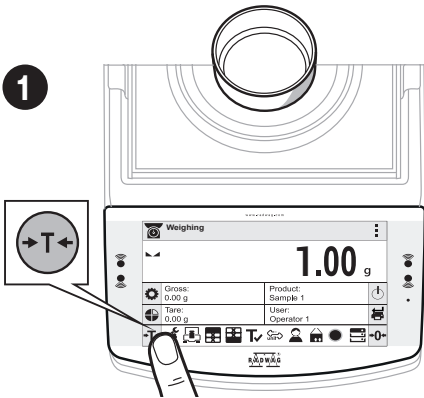
Tarowanie

2

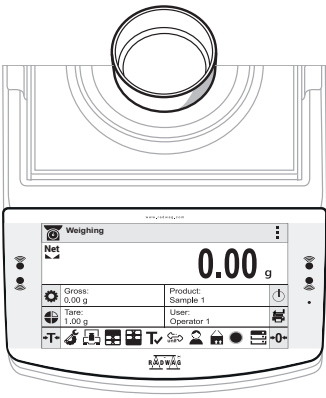With loaded weighing pan: upon stabilisation of the weighing result, press “Tare” button.

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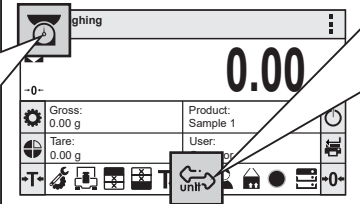
9. SETTINGS

USTAWIENIA

Some balance settings are accessible for Administrator exclusively.

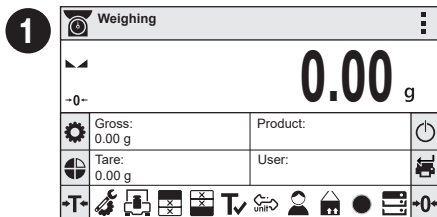
Prior balance parameters setup, log in as Administrator.

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
Logowanie administratora


Press "Operators" button.

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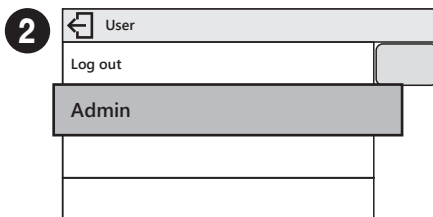Users list is displayed, select "Admin" option.

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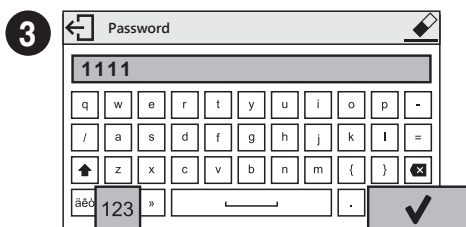
Press **123** button to activate numeric keyboard.

Enter „1111” password and press **✓** button to confirm.

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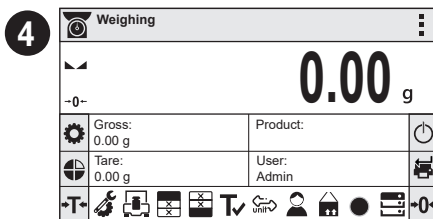You have logged as the Administrator.

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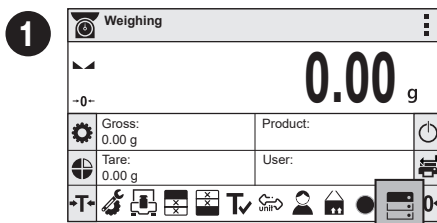
9.2. User settings

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
Prior to balance users setup, log in as Administrator (point 9.1.).

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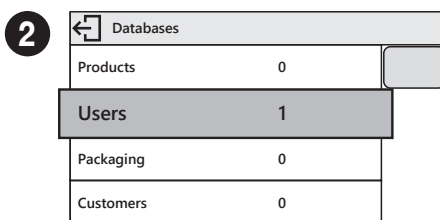4

New	
Name	New
Code	
Password	*****
Permissions	User

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6

New	
Name	John Smith
Code	
Password	*****
Permissions	User

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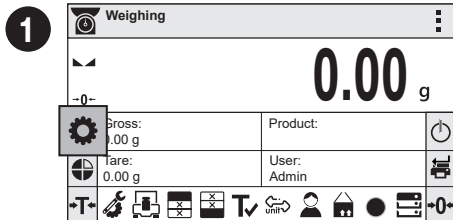
Prior balance proximity sensors setup, log in as Administrator (point 9.1.).

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
You can assign the proximity sensors with different functions for each working mode (the sensors will trigger different operation for each working mode). See the below procedure of sensors setup for the weighing mode.

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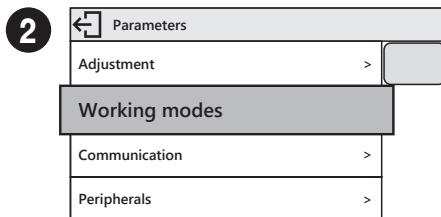
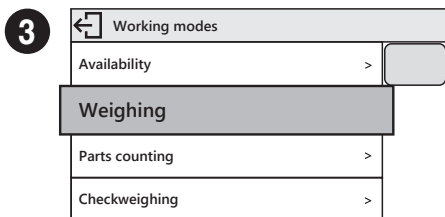
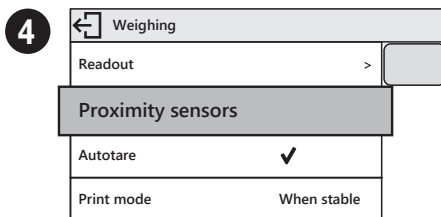
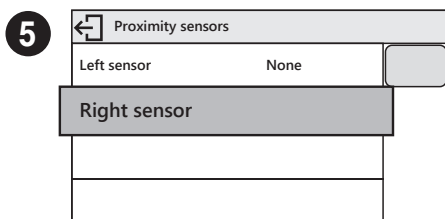
Select the appropriate sensor (right or left).

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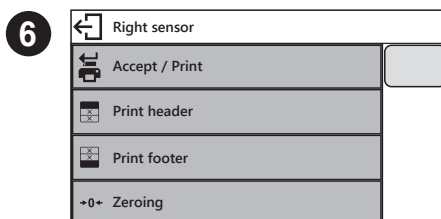Select function that is to be assigned to the chosen sensor.

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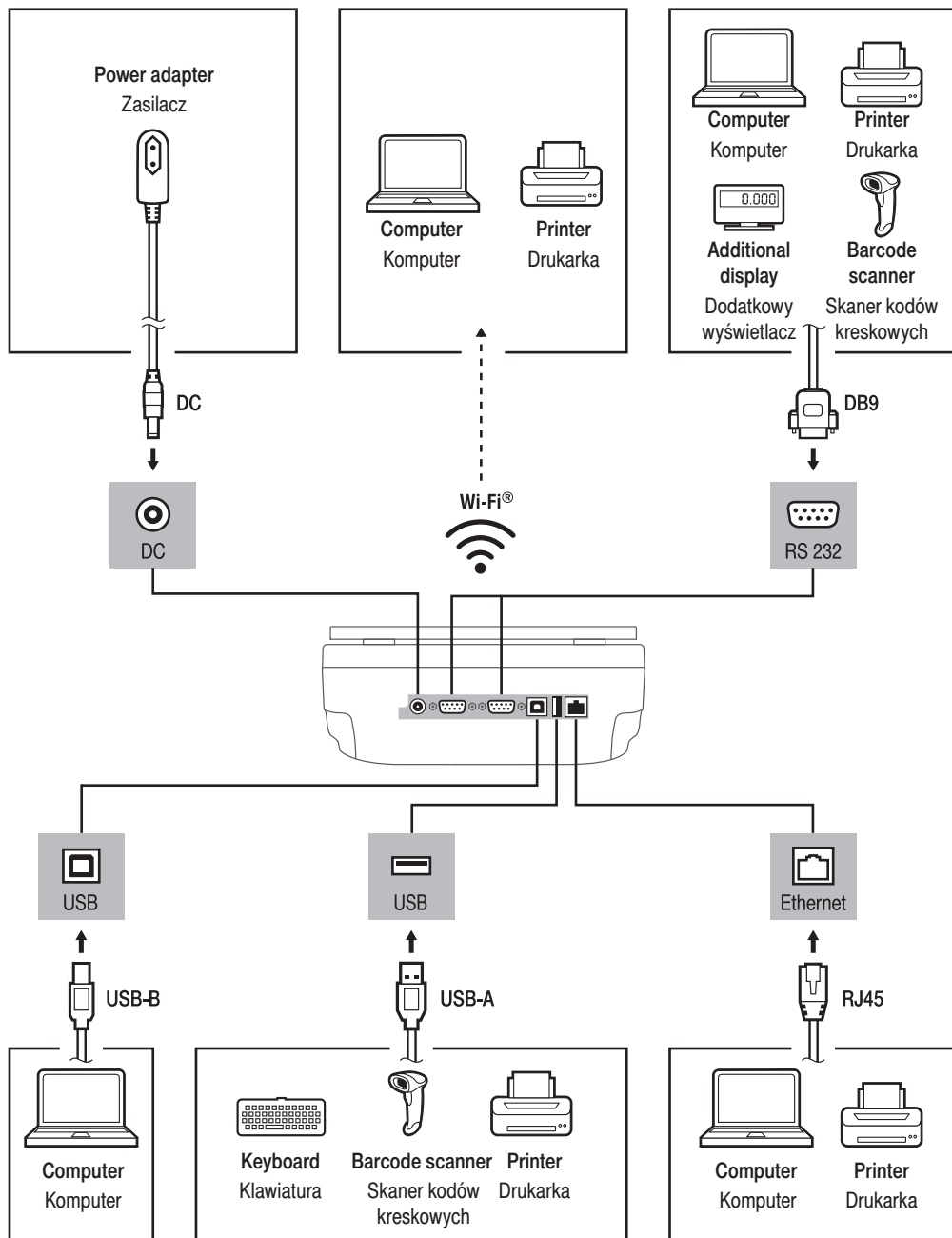
Following the above procedure you can configure sensors for other working modes.

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
11. DEVICE CLEANING

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
11.1. Guidelines and precautions

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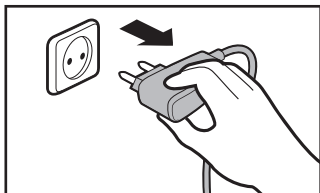
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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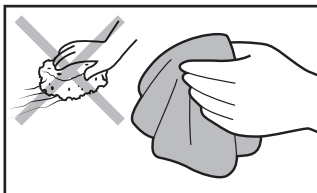Use soft cloth made of either natural or synthetic fiber exclusively. Avoid using cloth made of stiff/hard fabric, this could cause scratches on the surface!

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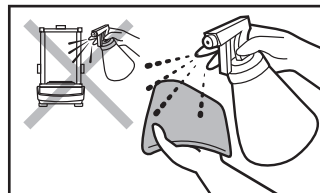Apply the cleanser onto the cloth. Mind not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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Shall it be necessary to disassemble the weighing chamber, follow the description provided in section 11.3.

- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either brush or handheld vacuum cleaner.
- Clean components can be dried using soft cloth or dust-free paper towel, with this the remaining moist will be absorbed.
- Install all the components providing they are completely dry.

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11.3. Weighing chamber elements disassembly

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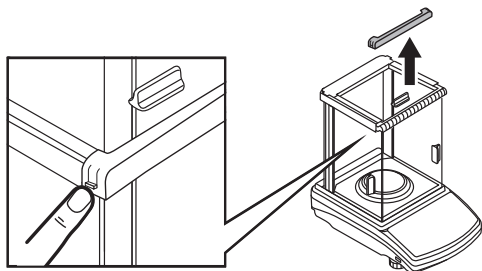
Prior disassembling weighing chamber elements make sure the balance is turned off and unplugged from the mains.

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
1

Press the moldings on the housing and lift the movable element up.

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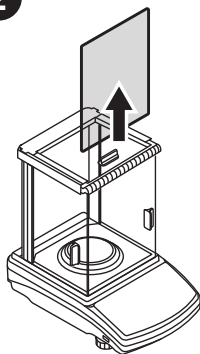
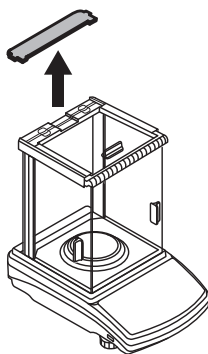
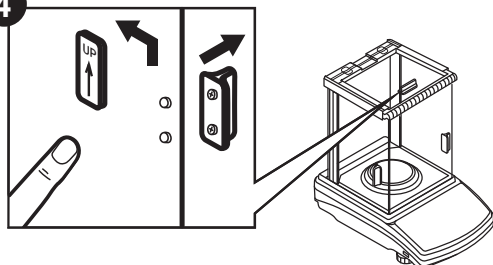
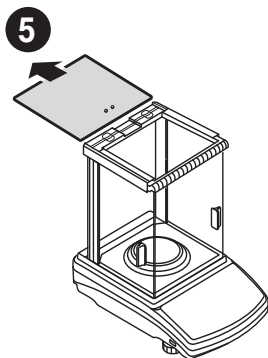
Now you can take out the upper glass.

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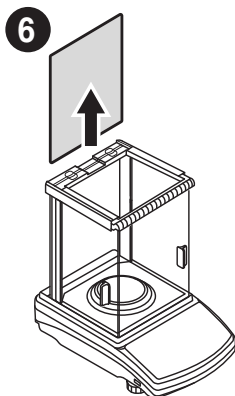
Take out the back glass.

The housing has recesses to facilitate removing the glass.

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
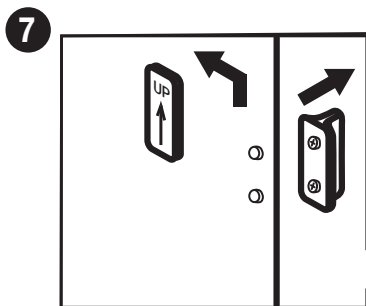
Remove the side door handles.

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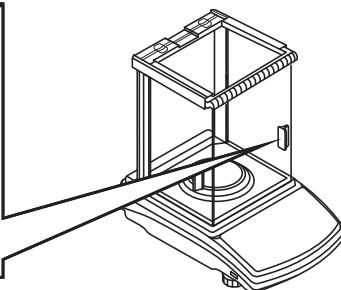Slide out the door.

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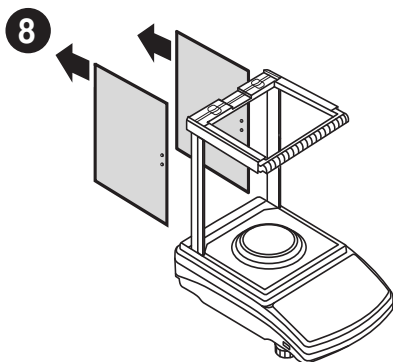All disassembled elements are prepared to be cleaned. Assembly all weighing chamber elements in an order reverse to the disassembly process.

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Wi-Fi® is a registered trademark of Wi-Fi Alliance.