



radwag.com

Scan the QR code to view additional scientific material that may be of interest to you.
There you will find more useful information in an accessible format!

Software manual

ITKP-09-03-12-25-EN

MODBUS

Communication protocol:

PUE 7.1 Weighing Terminal

PUE HY10 Weighing Terminal

WLY Precision Balances

WPY Multifunctional Scales

HY10 Multifunctional Scales

DECEMBER 2025

CONTENTS

1. GENERAL INFORMATION	4
2. IMPLEMENTED FUNCTIONS.....	4
3. MEMORY MAP	4
3.1. Input Address.....	4
3.2. Output Address.....	7

1. GENERAL INFORMATION

Modbus protocol implemented in the indicator can be applied when Ethernet are used (Modbus TCP).

2. IMPLEMENTED FUNCTIONS

Modbus communication is based on 3 functions:

- 03 (0x03) Read Holding Registers – output data readout.
- 04 (0x04) Read Input Registers – input data readout.
- 16 (0x10) Write Multiple Registers – output data record.

3. MEMORY MAP

3.1. Input Address

Input variables list:

Variable	Address	Length [WORD]	Data type
Platform 1 mass	0	2	float
Platform 1 tare	2	2	float
Platform 1 unit	4	1	word
Platform 1 status	5	1	word
Platform 1 LO threshold	6	2	float
Platform 2 mass	8	2	float
Platform 2 tare	10	2	float
Platform 2 unit	12	1	word
Platform 2 status	13	1	word
Platform 2 LO threshold	14	2	float
Process status (Stop, Start)	32	1	word
Inputs status	33	1	word
Min	34	2	float
Max	36	2	float
Lot number	42	2	dword
Operator	44	1	word
Product	45	1	word
Customer	46	1	word
Packaging	47	1	word
Source warehouse	48	1	word
Target warehouse	49	1	word
Formulation/Dosing	50	1	word

Platform mass – response: platform mass in current unit.

Platform tare – response: platform tare in an adjustment unit.

Platform unit – determines currently displayed mass unit of a platform.

Unit bits	
0	gram [g]
1	kilogram [kg]
2	carat [ct]
3	pound [lb]
4	ounce [oz]
5	Newton [N]

Example:

bit No.	B5	B4	B3	B2	B1	B0
value	0	0	0	0	1	0

The unit of the weighing instrument is kilogram [kg].

Platform status – determines status of a weighing platform.

Status bits	
0	Measurement correct (weighing instrument does not report an error).
1	Measurement stable.
2	Weighing instrument indicates zero.
3	Weighing instrument is tared.
4	Weighing instrument is in II weighing range.
5	Weighing instrument is in III weighing range.
6	Weighing instrument reports NULL error.
7	Weighing instrument reports LH error.
8	Weighing instrument reports FULL error.

Example:

bit No.	B8	B7	B6	B5	B4	B3	B2	B1	B0
value	0	0	0	0	1	0	0	1	1

The weighing instrument does not report error, stable measurement in II weighing range.

LO threshold – response: **LO** threshold value of a platform in an adjustment unit.

Process status – determines process status:

Decimal value	Process status	bit No.	
		B1	B0
0	process disabled	0	0
1	process start	0	1
2	process stop	1	0
3	process completed	1	1

Inputs status – response: status of set inputs:

Input No.	12	11	10	9	8	7	6	5	4	3	2	1
OFF	0	0	0	0	0	0	0	0	0	0	0	0
ON	1	1	1	1	1	1	1	1	1	1	1	1

Example:

Mask of set 2 and 4 inputs: 0000 0000 0000 1010

MIN – response: **MIN** threshold value (in the current unit selected for active working mode).

MAX – response: **MAX** threshold value (in the current unit selected for active working mode).

Lot number – response: lot number.

Operator – response: code of logged in operator.

Product – response: code of selected product.

Customer – response: code of selected customer.

Packaging – response: code of selected packaging.

Source warehouse – response: code of selected source warehouse.

Target warehouse – response: code of selected target warehouse.

Formulation – response: code of selected formulation.

3.2. Output Address

Output variables list:

Variable	Address	Length [word]	Data type
Command	0	1	word
Command with parameter	1	1	word
Platform	2	1	word
Tare	3	2	float
LO threshold	5	2	float
Outputs status	7	1	word
Min	8	2	float
Max	10	2	float
Lot number	16	2	dword
Operator	18	1	word
Product	19	1	word
Customer	20	1	word
Packaging	21	1	word
Source warehouse	22	1	word
Target warehouse	23	1	word
Formulation/Dosing	24	1	word

Basic command – setting respective value performs the task in accordance with the table:

Decimal value	Command
1	Zero the platform
2	Tare the platform
8	Delete statistics
16	Save/Print
32	Start
64	Stop (error)

Example:

0000 0000 0010 0000 – process start.

Complex command – setting respective value performs the task in accordance with the table:

Decimal value	Command
1	Setting tare value for a given platform
2	Setting LO threshold value for a given platform
3	Setting lot number
4	Setting outputs status
5	Operator selection
6	Product selection
7	Packaging selection
8	Setting MIN threshold value
9	Customer selection
10	Source warehouse selection
11	Target warehouse selection
12	Dosing selection
16	Setting MAX threshold value

	Complex command requires setting address of respective parameter (from 2 to 24 – refer to: "Complex command parameters").
---	--

Example:

0000 0000 0000 0010 – command sets LO threshold to the value set in LO parameter (address 5 – refer to: "Complex command parameters").

Platform – complex command parameter: weighing platform number.

Tare – complex command parameter: tare value (in an adjustment unit).

LO threshold – complex command parameter: LO threshold value (in an adjustment unit).

Outputs status – complex command parameter: determines status of weighing indicator outputs.

Output No.	12	11	10	9	8	7	6	5	4	3	2	1
OFF	0	0	0	0	0	0	0	0	0	0	0	0
ON	1	1	1	1	1	1	1	1	1	1	1	1

Example:

Mask of active 2 and 4 outputs: 0000 0000 0000 1010

MIN – complex command parameter: MIN threshold value (in the current unit selected for active working mode).

MAX – complex command parameter: MAX threshold value (in the current unit selected for active working mode).

Lot number – complex command parameter: lot number.

Operator – complex command parameter: code of logged in operator.

Product – complex command parameter: code of selected product.

Customer – complex command parameter: code of selected customer.

Packaging – complex command parameter: code of selected packaging.

Source warehouse – complex command parameter: code of selected source warehouse.

Target warehouse – complex command parameter: code of selected target warehouse.

Formulation – complex command parameter: code of selected formulation.



A command or a command with parameter is executed once when its bit setting is detected. If the command with the same bit is to be executed again, zero the bit.

Example:

Command	address 1	address 0
Tare	0000 0000	0000 0010
Zero command bits	0000 0000	0000 0000
Tare	0000 0000	0000 0010

