



radwag.com


Scan the QR code to view additional scientific material that may be of interest to you.
There you will find more useful information in an accessible format!

Software manual

ITKP-14-01-01-20-EN

PROFIBUS

Communication Protocol of PUE HY10 Terminal

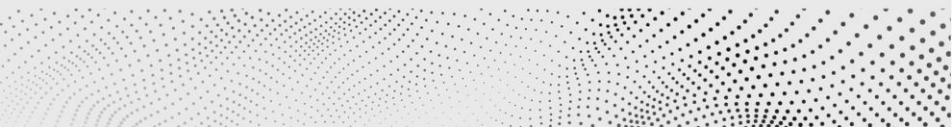

JANUARY 2020

CONTENTS

1. GENERAL INFORMATION	4
2. MEMORY MAP.....	5
2.1. Output address	5
2.2. Input address	6
3. DESCRIPTION OF VARIABLES.....	6
3.1. Output variables.....	6
3.2. Input variables.....	10

1. GENERAL INFORMATION

Profibus Communication Module ensures data exchange between a supervising controlling device (master) and a scale (Slave) in accordance with the Profibus DP protocol.

The supervising unit enables:

- Cyclic reading in input signals from an indicator PUE HY10 series,
- Cyclic saving outputs status to an indicator PUE HY10 series.

Profibus communication functionality with the indicator PUE HY10 enables:

- Operation of four weighing platforms,
- Tarring,
- Zeroing,
- Setting tare value,
- Setting the value of LO limit,
- Setting the value of Min threshold,
- Setting the value of Max threshold,
- Reading inputs status,
- Setting outputs,
- Selecting an operator,
- Selecting a product,
- Selecting a client,
- Selecting a packaging,
- Selecting a source warehouse,
- Selecting a destination warehouse,
- Selecting a formulation,
- Setting lot number,
- Process stopping,
- Process starting,
- Saving / Printing,
- Statistics zeroing.

2. MEMORY MAP

2.1. Output address

Address Offset	0	1	2	3	4	5	6	7	8	9
0	M 1	M 1	M 1	M 1	T 1	T 1	T 1	T 1	J 1	J 1
1	S 1	S 1	LO 1	LO 1	LO 1	LO 1	M 2	M 2	M 2	M 2
2	T 2	T 2	T 2	T 2	J 2	J 2	S 2	S 2	LO 2	LO 2
3	LO 2	LO 2	M 3	M 3	M 3	M 3	T 3	T 3	T 3	T 3
4	J 3	J 3	S 3	S 3	LO 3	LO 3	LO 3	LO 3	M 4	M 4
5	M 4	M 4	T 4	T 4	T 4	T 4	J 4	J 4	S 4	S 4
6	LO 4	LO 4	LO 4	LO 4	ST	ST	SW	SW	MIN	MIN
7	MIN	MIN	MAX	MAX	MAX	MAX	-	-	-	-
8	-	-	-	-	LOT	LOT	LOT	LOT	O	O
9	A	A	K	K	OK	OK	MZ	MZ	MD	MD
10	RC	RC	-	-	-	-	-	-	-	-

Where:

M	Mass of a weighing platform, 4 bytes, float
T	Tare of a weighing platform, 4 bytes, float
J	Measuring unit of a weighing platform, 2 bytes, word
S	Status of a weighing platform, 2 bytes, word
LO	Lo limit of a weighing platform, 4 bytes, float
MIN	MIN threshold, 4 bytes, float
MAX	MAX threshold, 4 bytes, float
LOT	Lot, 4 bytes, word
O	Operator, 2 bytes, word
A	Product, 2 bytes, word
K	Client, 2 bytes, word
OK	Packages, 2 bytes, word
MZ	Source warehouse, 2 bytes, word
MD	Destination warehouse, 2 bytes, word
RC	Formulation, 2 bytes, word

2.2. Input address

Address Offset	0	1	2	3	4	5	6	7	8	9
0	C	C	CP	CP	P	P	T	T	T	T
1	LO	LO	LO	LO	SW	SW	MIN	MIN	MIN	MIN
2	MAX	MAX	MAX	MAX	-	-	-	-	-	-
3	-	-	LOT	LOT	LOT	LOT	O	O	A	A
4	K	K	OK	OK	MZ	MZ	MD	MD	RC	RC

Where:

C	Command, 2 bytes, word
CP	Command with a parameter, 2 bytes, word
P	Active weighing platform, 2 bytes, word
T	Tare of a weighing platform, 4 bytes, float
LO	Lo limit of a weighing platform, 4 bytes, float
SW	Inputs/Outputs statuses, 2 bytes, word
MIN	MIN threshold, 4 bytes, float
MAX	MAX threshold, 4 bytes, float
LOT	Lot, 4 bytes, word
O	Operator, 2 bytes, word
A	Product, 2 bytes, word
K	Client, 2 bytes, word
OK	Packages, 2 bytes, word
MZ	Source warehouse, 2 bytes, word
MD	Destination warehouse, 2 bytes, word
RC	Formulation, 2 bytes, word

3. DESCRIPTION OF VARIABLES

3.1. Output variables

Reading the output variables enables obtaining data on device status.

List of output variables:

Variable	Address	Length [word]	Data type
Mass of platform 1	0	2	float
Tare of platform 1	4	2	float
Measuring unit of platform 1	8	1	word
Status of platform 1	10	1	word
Lo limit of platform 1	12	2	float
Mass of platform 2	16	2	float
Tare of platform 2	20	2	float
Measuring unit of platform 2	24	1	word
Status of platform 2	26	1	word
Lo limit of platform 2	28	2	float
Mass of platform 3	32	2	float
Tare of platform 3	36	2	float
Measuring unit of platform 3	40	1	word
Status of platform 3	42	1	word
Lo limit of platform 3	44	2	float
Mass of platform 4	48	2	float
Tare of platform 4	52	2	float
Measuring unit of platform 4	56	1	word
Status of platform 4	58	1	word
Lo limit of platform 4	60	2	float
Process status (Stop, Start)	64	1	word
Inputs status	66	1	word
Min	68	2	float
Max	72	2	float
Lot number	84	2	word
Operator	88	1	word
Product	90	1	word
Client	92	1	word
Packaging	94	1	word
Source warehouse	96	1	word
Destination warehouse	98	1	word
Formulation	100	1	word

Mass of platform – response is mass on a weighing platform in current measuring unit.

Tare of platform – response is the value of tare on a weighing platform in adjustment unit.

Measuring unit of a platform – determines current (displayed) measuring unit set for a weighing platform.

Measuring unit bits	
0	gram [g]
1	kilogram [kg]
2	carat [ct]
3	pound [lb]
4	ounce [oz]
5	Newton [N]

Example:

Bit no.	B5	B4	B3	B2	B1	B0
Value	0	0	0	0	1	0

The scale measures with a unit: kilogram [kg].

Status of a platform – determines status of a weighing platform.

Bity statusu	
0	correct measurement (the scale does not report an error)
1	stable measurement
2	scale in precise zero
3	scale tarred
4	scale in 2nd measuring range
5	scale in 3rd measuring range
6	scale reports NULL error
7	scale reports LH error
8	scale reports FULL error

Example:

Bit no.	B8	B7	B6	B5	B4	B3	B2	B1	B0
Value	0	0	0	0	1	0	0	1	1

The scale does not report an error, the measurement is stabilized in the 2nd measuring range.

LO – response is the value of **LO** limit in an adjustment unit of a given weighing platform.

Process status – determines status of a process:

Decimal value of a variable	Process status	Bit no.	
		B1	B0
0	Process inactive	0	0
1	Process start	0	1
2	Process stop	1	0
3	Process end	1	1

Inputs status – response is the status of set inputs:

Input no.	12	11	10	9	8	7	6	5	4	3	2	1
OFF	0	0	0	0	0	0	0	0	0	0	0	0
ON	1	1	1	1	1	1	1	1	1	1	1	1

Example:

Mask of set inputs 2 and 4: 0000 0000 0000 1010

MIN – response is the value of set **MIN** threshold (in a measuring unit of an enabled working mode).

MAX – response is the value of set **MAX** threshold (in a measuring unit of an enabled working mode).

Lot number – response is the value of lot number.

Operator – response is the value of a logged operator.

Product – response is the value of a selected product.

Client – response is the value of a code of a selected client.

Packaging – response is the value of a code of a selected packaging.

Source warehouse – response is the value of a code of a source warehouse.

Destination warehouse – response is the value of a code of a destination warehouse.

Formulation – response is the value of a code of a selected formulation.

3.2. Input variables

Saving input variables in an indicator PUE HY10 series enables influencing its operation.

List of input variables:

Variable	Address	Length [word]	Data type
Command	0	1	word
Command with a parameter	2	1	word

List of parameters of a complex command:

Parameter	Address	Length [word]	Data type
Platform	4	1	word
Tare	6	2	float
LO limit	10	2	float
Output status	14	1	word
Min	16	2	float
Max	20	2	float
Lot number	32	2	word
Operator	36	1	word
Product	38	1	word
Client	40	1	word
Packaging	42	1	word
Source warehouse	44	1	word
Destination warehouse	46	1	word
Formulation	48	1	word

Basic command – Setting a bit causes carrying out a task as specified in a below table:

Command bit	Command
0	Zero platform
1	Tare platform
3	Clear statistics
4	Save / Print
5	Start
6	Stop

Example:

0000 0000 0010 0000 – the command carries out process start.

Complex command – Setting a bit causes carrying out a task as specified in a below table:

Command bit	Command
0	Setting tare value of a weighing platform
1	Setting the value of LO limit of a weighing platform
2	Setting outputs status
3	Setting the value of MIN threshold
4	Setting the value of MAX threshold

	<i>A complex command requires setting an appropriate parameter (addresses from 4 to 48. – see table “List of parameters in a complex command”).</i>
-----------------------------------------------------------------------------------	------------------------------------------------------------------------------------------------------------------------------------------------------------

Example:

0000 0000 0000 0010 – a command carries out setting of the LO limit for a value given in parameter LO (address 10 – see table “List of parameters in a complex command”).

Platform – complex command parameter: number of a weighing platform.

Tare – complex command parameter: tare value (in an adjustment unit).

LO – complex command parameter: the value of LO limit (in an adjustment unit).

Outputs status – complex command parameter: determines outputs status of an indicator.

Output no.	12	11	10	9	8	7	6	5	4	3	2	1
OFF	0	0	0	0	0	0	0	0	0	0	0	0
ON	1	1	1	1	1	1	1	1	1	1	1	1

Example:

Mask of active outputs 2 and 4: 0000 0000 0000 1010

MIN – complex command parameter: the value of MIN threshold (in a measuring unit of an enabled working mode).

MAX – complex command parameter: the value of MAX threshold (in a measuring unit of an enabled working mode).

Lot number – complex command parameter: the value of lot number.

Operator – complex command parameter: the value of a code of a logged operator.

Product – complex command parameter: the value of a code of a selected product.

Client – complex command parameter: the value of a code of a selected client.

Packaging – complex command parameter: the value of a code of a selected packaging.

Source warehouse – complex command parameter: the value of a code of a selected source warehouse.

Destination warehouse – complex command parameter: the value of a code of a selected destination warehouse.

Formulation – complex command parameter: the value of a code of a selected formulation.

	<p><i>A command or a command with a parameter is carried out once on detecting the setting of a corresponding bit. If it is necessary to repeat a command with the same bit, then first it has to be zeroed.</i></p>
------------------------------------------------------------------------------------	-----------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------

Example:

Command	address 1	address 0
Tarring	0000 0000	0000 0010
Zeroing command bits	0000 0000	0000 0000
Tarring	0000 0000	0000 0010

