IMMB-03-05-01-26-EN-PL

Start up Guide

Szybki start

Analytical and Precision Balances

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Analytical balances:

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- AS R1 PLUS
- AS R2 PLUS

Precision balances:

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- PS R1
- PS R1.M
- PS R2
- PS R2.M
- PS R2.H
- PS R2.M.H

To view a complete User Manual go to the website or scan QR code:

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
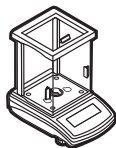
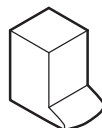
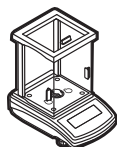
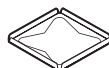
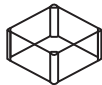
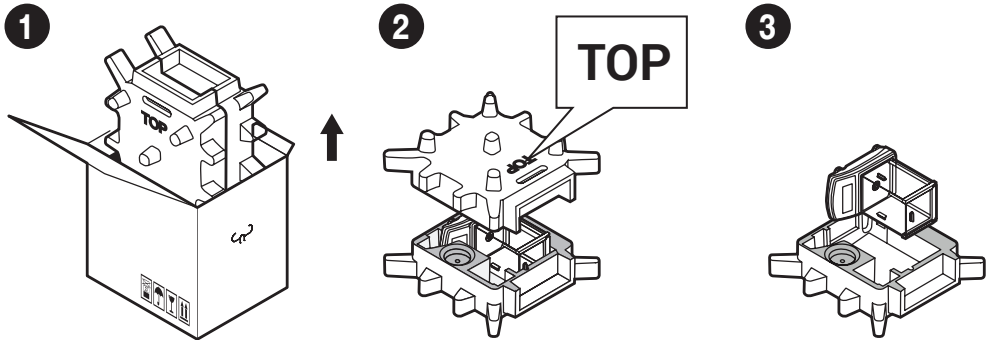
3. RECOMMENDATIONS AND WARNINGS

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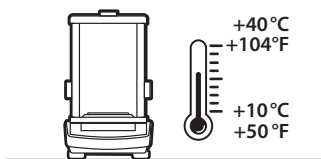
Operate the device in a room where the temperature ranges between 10–40 °C (50–104 °F) and where the relative humidity is below 80%.


Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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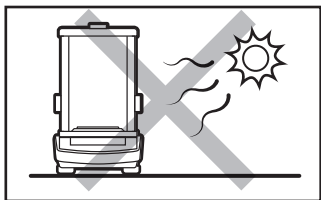
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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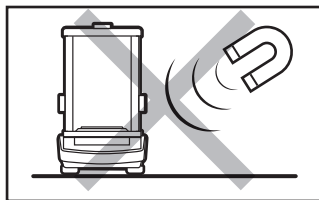
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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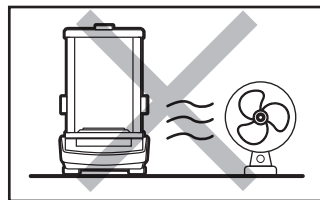
Avoid air drafts and air movements at the workstation.

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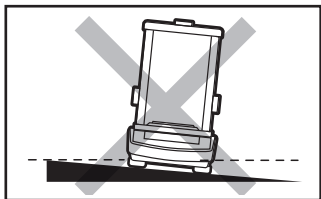
Make sure that the balance is placed on an even surface.

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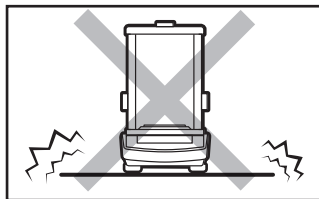Do not place the balance on an unstable ground exposed to shocks and vibrations.

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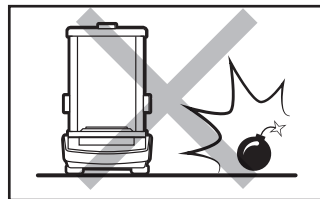Do not place the balance in a hazardous area. Do not weigh explosive materials.

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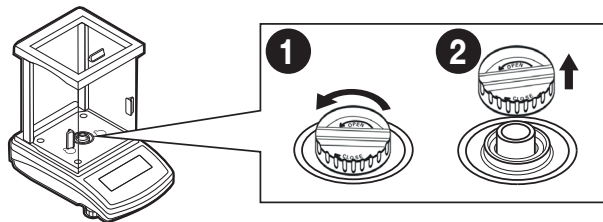
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Transport lock removal

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
AS R Series

Keep the transport lock should a warranty claim or service be required.

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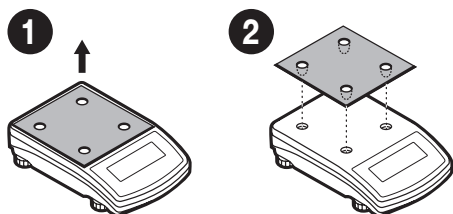
4.2. Grounding spring check

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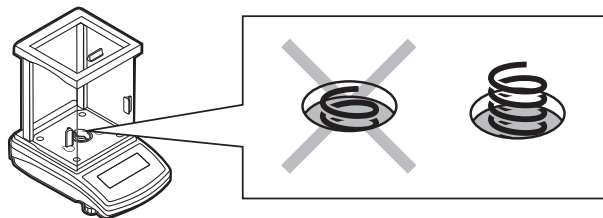AS R Series


Check whether the grounding spring takes its intended location. Make sure that the grounding spring juts slightly out of the hole.

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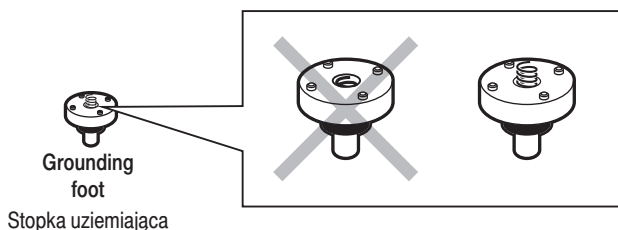
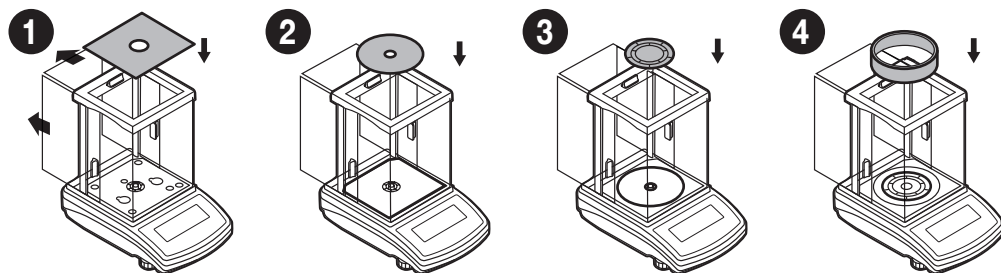
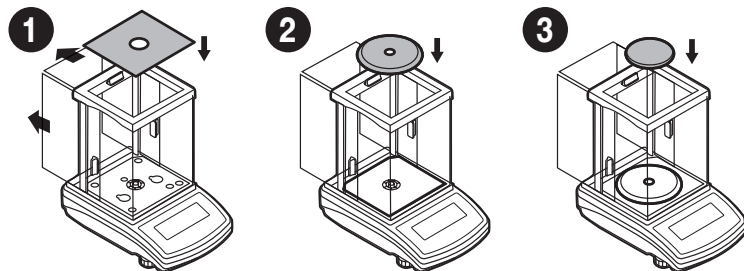
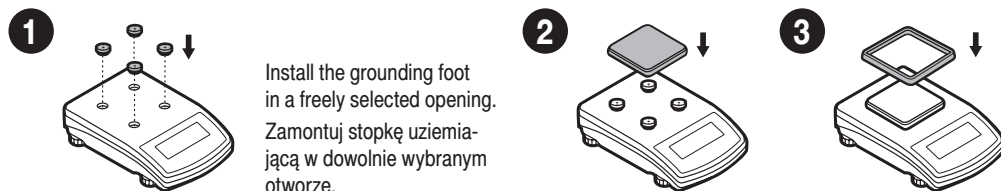
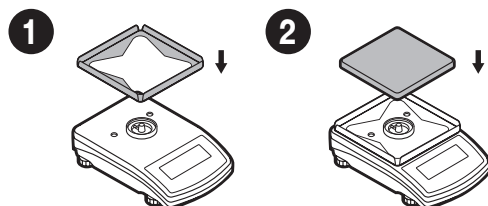
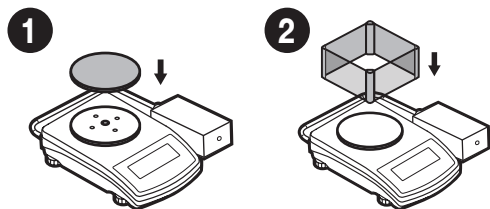
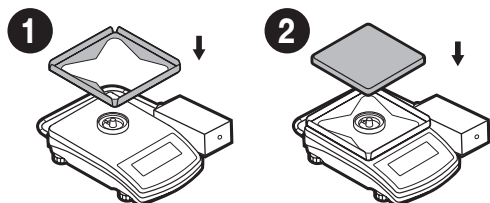
6. PREPARING FOR WORK

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6.1. Temperature stabilization

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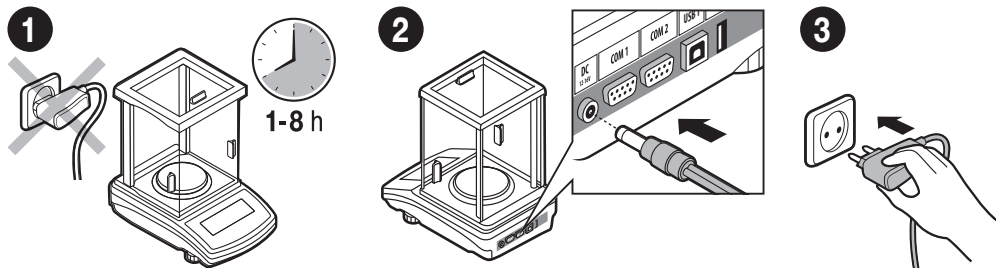Prior to switching the device on, let it reach room temperature (estimated stabilisation time: 1-8 hours).

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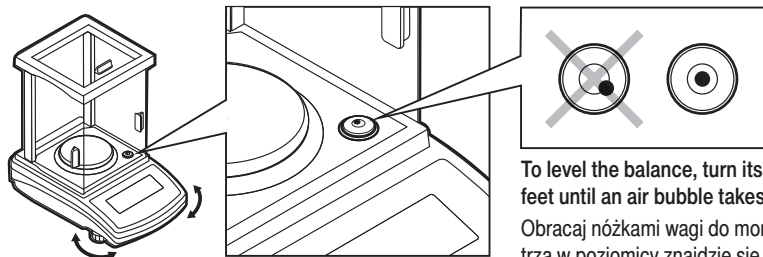
6.2. Leveling

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
To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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6.3. Switch the balance on

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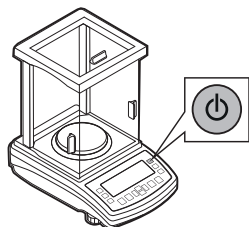Your balance is ready to use!

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To switch the device on, press ON / OFF key.

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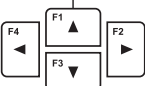
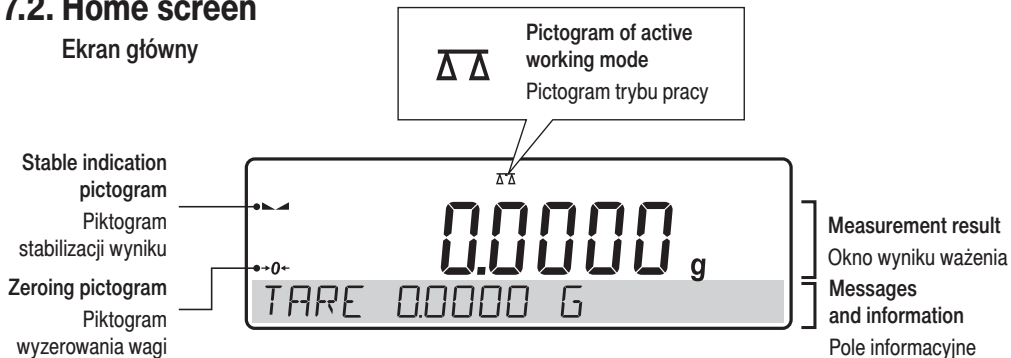
Ekran główny


8. BASIC OPERATIONS

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8.1.1. Adjustment (for models with an internal adjustment only)

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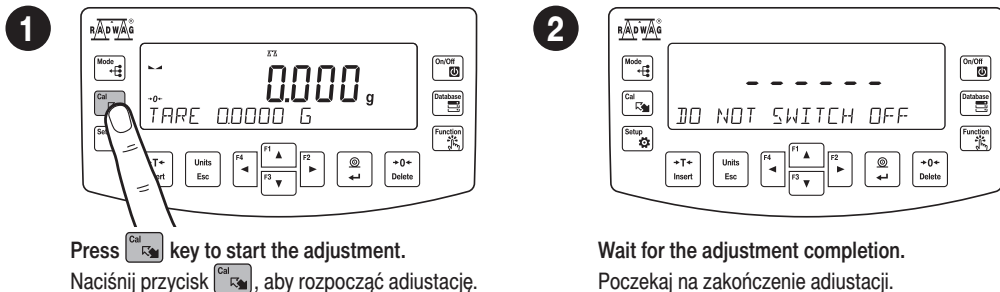8.1.2. Adjustment (for models with an external adjustment only)

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It is recommended to carry out adjustment operation prior to running the **WEIGHING** mode. External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on device type and capacity.

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The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use mass standard of class F1 or F2. View the table and select the mass standard you need.

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

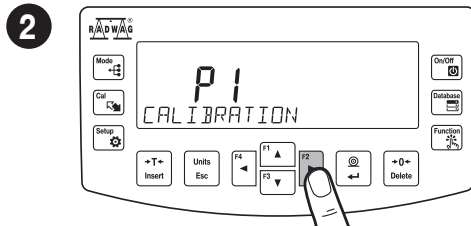
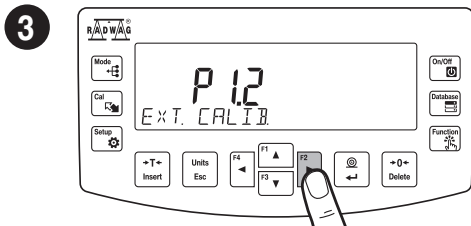
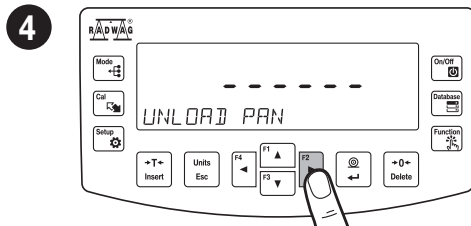
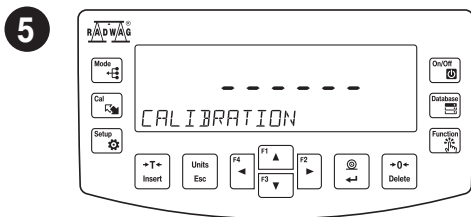
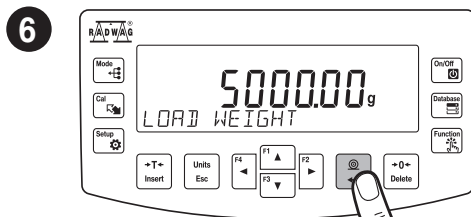
8.2. Zeroing

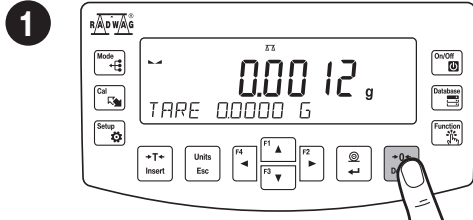Zerowanie

Make sure that the weighing pan is empty and press **0** key.

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
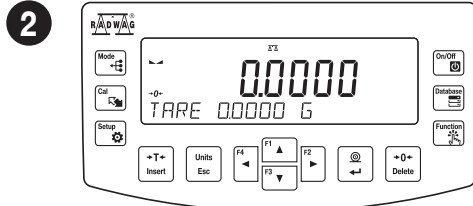
8.3. Taring

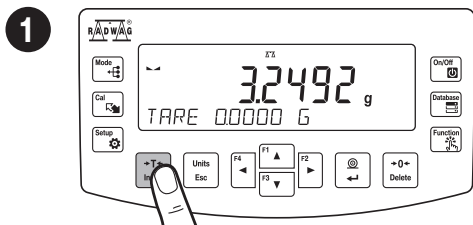Tarowanie

With loaded weighing pan: upon stabilisation of the weighing result, press **T** key.

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
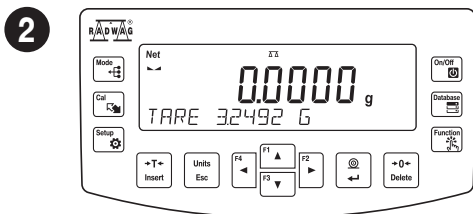
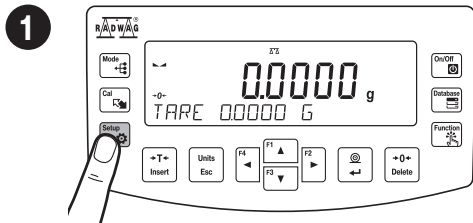
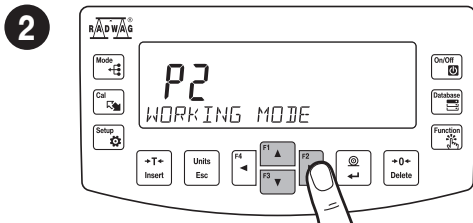
Using **F2** keys select "Working mode" and press **F2** key.

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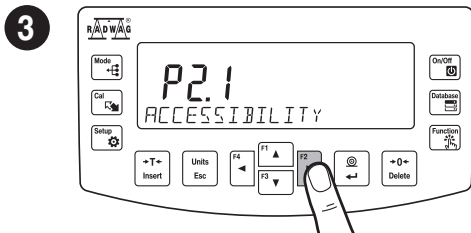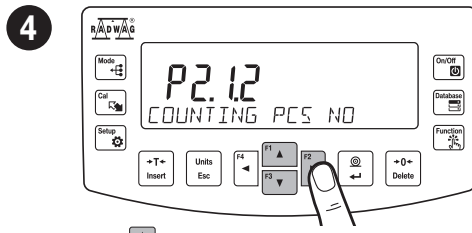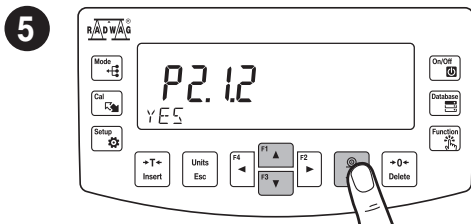
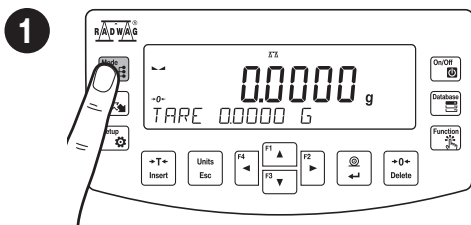
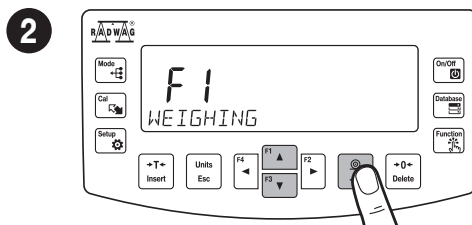
The menu features the following working modes:

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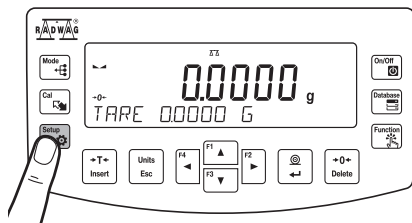
Detailed overview of the settings of the working modes is to be found in balance's main user manual.

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
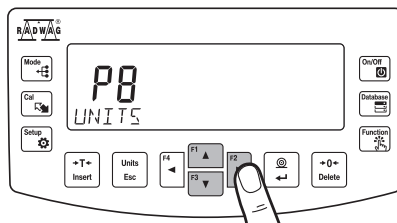
2

Using keys select "Units" and press key.

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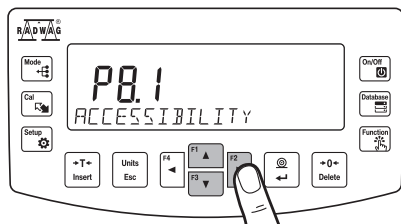
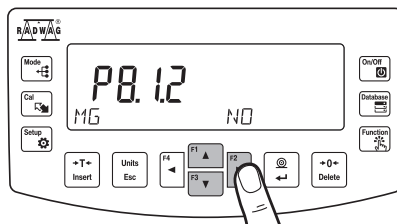
4

Using keys select unit for which accessibility is to be set and press key.

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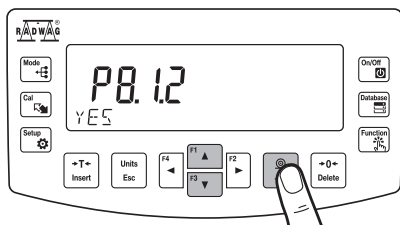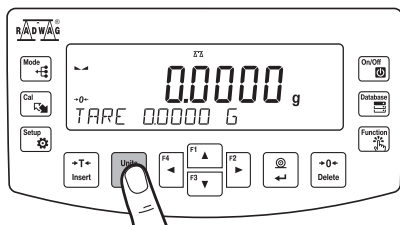
The balance enables selecting the following weighing units:

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* Units available for verified balances exclusively

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9. SETTINGS

USTAWIENIA

Some balance settings are accessible for Administrator exclusively.

Prior balance parameters setup, log in as Administrator.

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
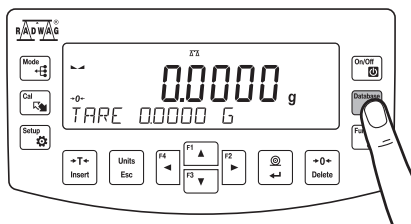
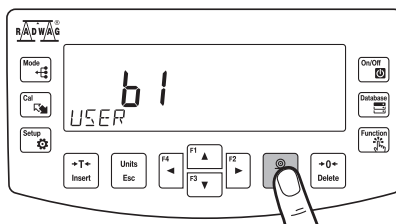
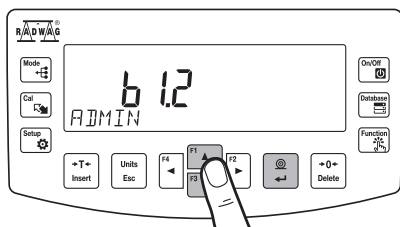
3

Using keys select "Admin" and press key to confirm.

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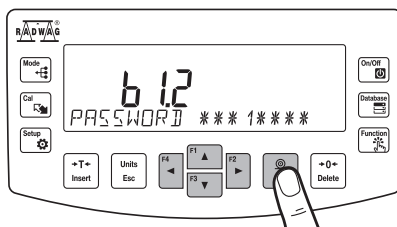
4


Using keys select "1111" and press key to confirm.

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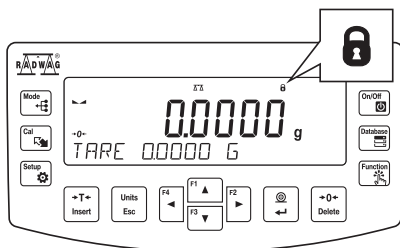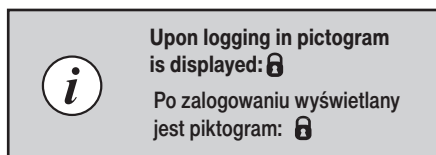9.2. User settings

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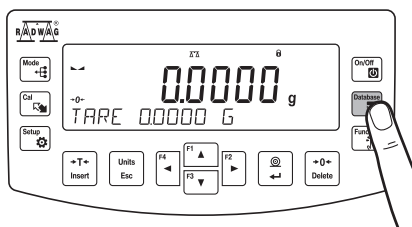
Prior to balance users setup, log in as Administrator (point 9.1).

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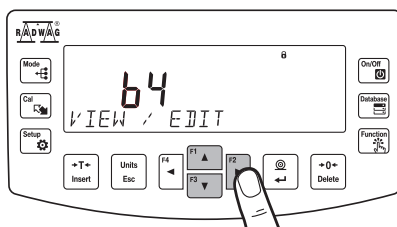
2

Using keys select "View/Edit" option and press key to confirm.

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
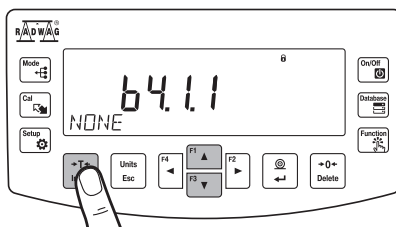
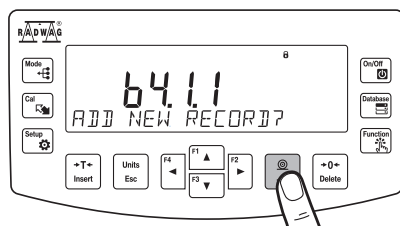
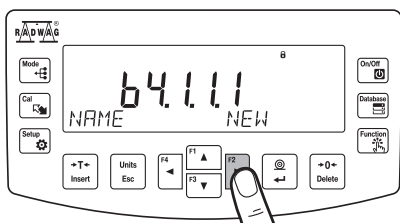
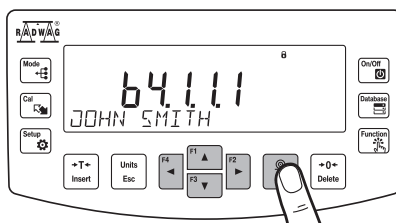
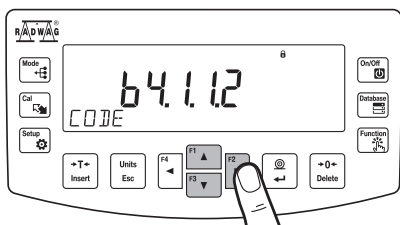
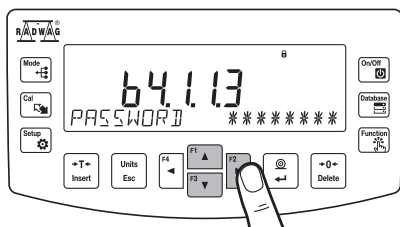
10

Using keys select "Password" and press key to confirm.

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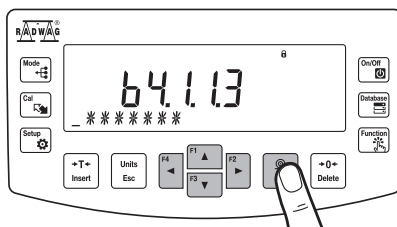
11

Using keys enter user password and press key to confirm.

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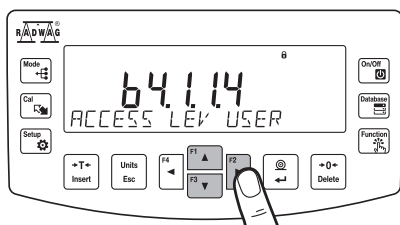
12

Using keys select "Access Level" and press key to confirm.

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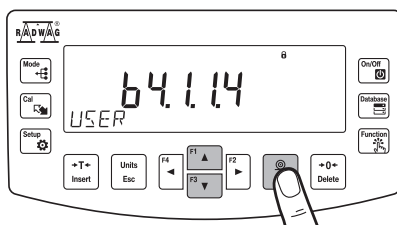
13

Using keys select permissions level and press key to confirm.

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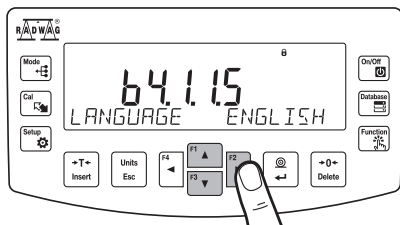
14

Using keys select "Language" and press key to confirm.

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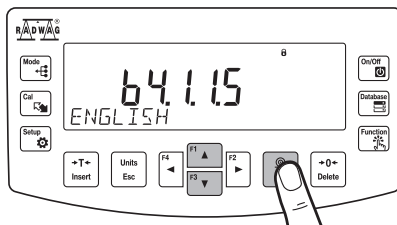15


Using keys select menu language and press key to confirm.

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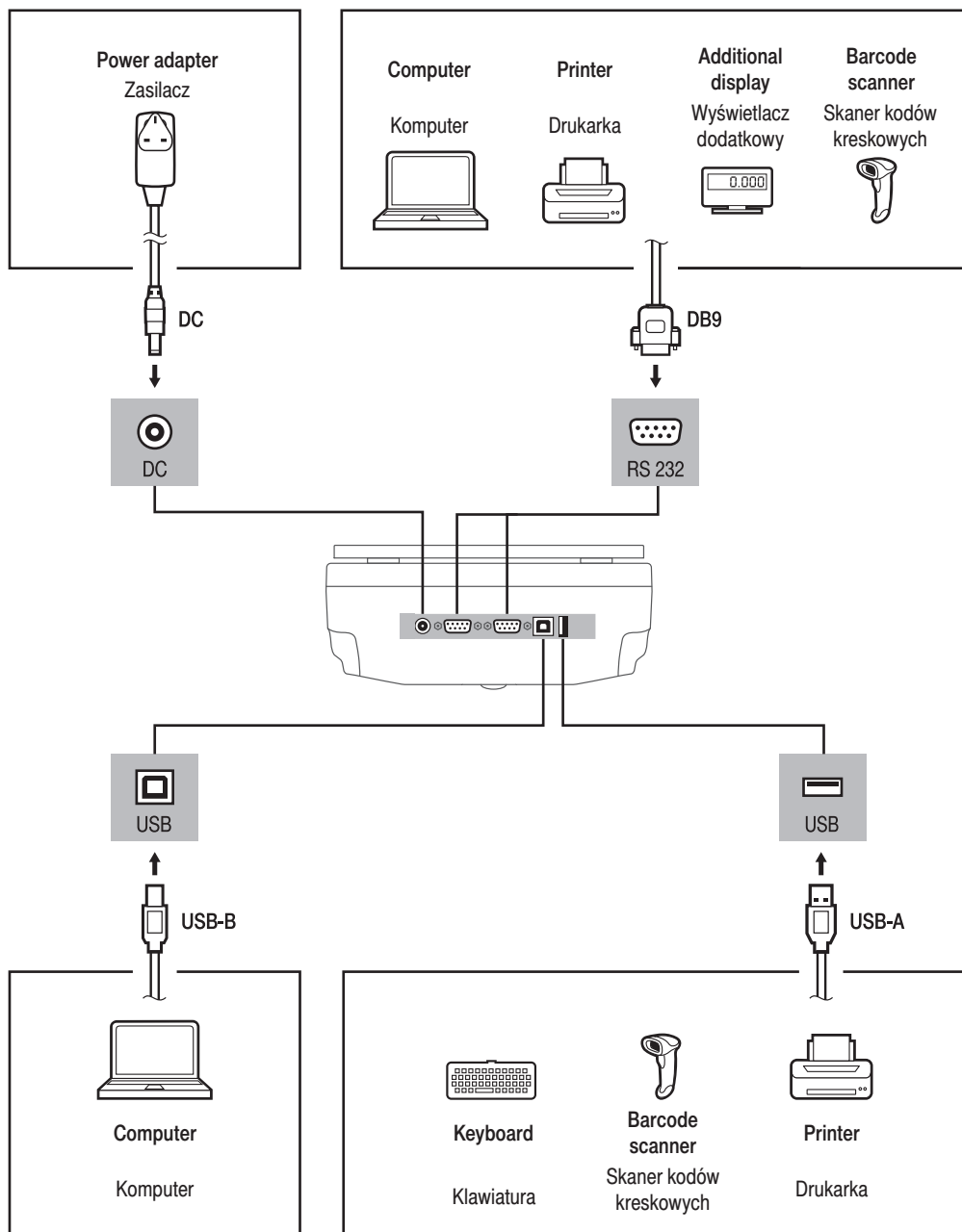
11. DEVICE CLEANING

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
11.1. Guidelines and precautions

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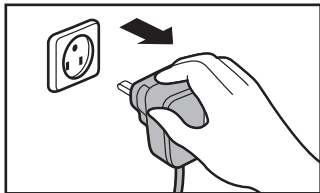
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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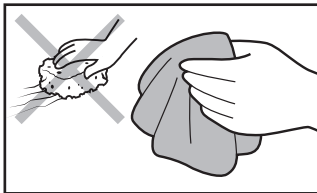Use soft cloth made of either natural or synthetic fiber exclusively. Avoid using cloth made of stiff/hard fabric, this could cause scratches on the surface!

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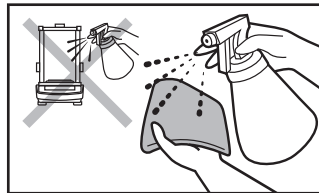Apply the cleanser onto the cloth. Mind not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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Shall it be necessary to disassemble the weighing chamber, follow the description provided in section 11.3.

- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either brush or handheld vacuum cleaner.
- Clean components can be dried using soft cloth or dust-free paper towel, with this the remaining moist will be absorbed.
- Install all the components providing they are completely dry.

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11.3. Weighing chamber elements disassembly

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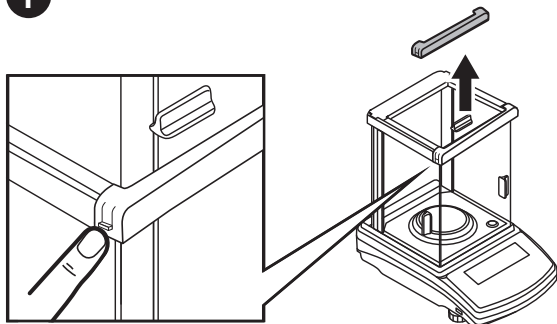
Prior disassembling weighing chamber elements make sure the balance is turned off and unplugged from the mains.

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
1

Press the moldings on the housing and lift the movable element up.

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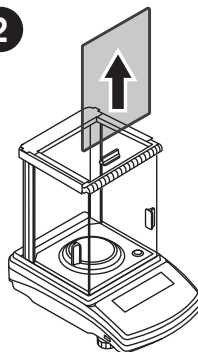
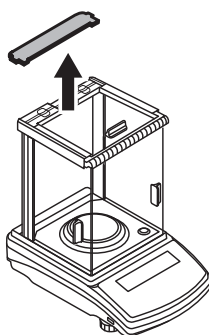
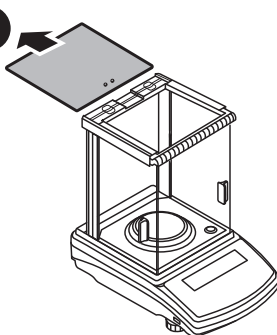
Now you can take out the upper glass.

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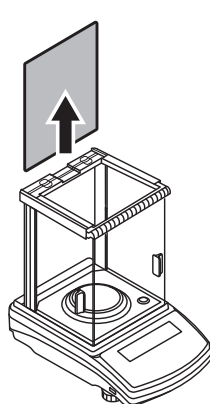
5

Take out the back glass. The housing has recesses to facilitate removing the glass.

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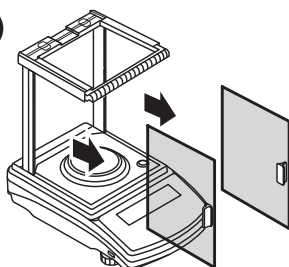6

Slide out the door.

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After cleaning the weighing chamber elements, they must be assembled in the reverse order.

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