IMMB-08-03-06-24-EN-PL

**Start up Guide**

**Szybki start**

# **Precision balances**

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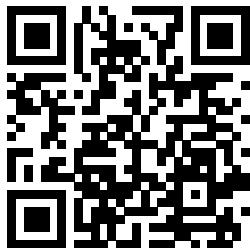
---

- WTC

To view a complete User Manual go to the website or scan QR code:

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
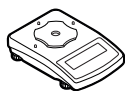
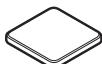
## 2. UNPACKING

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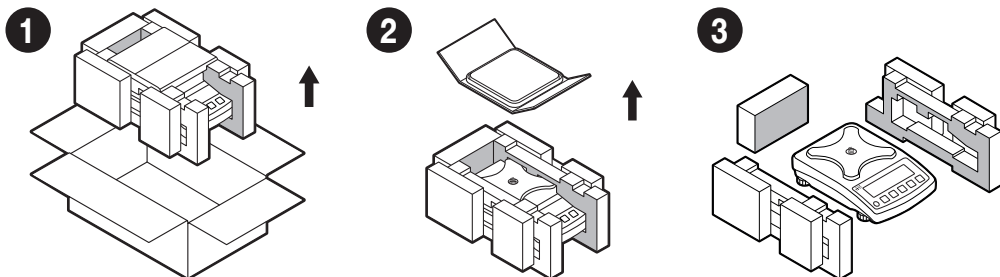
Keep the packaging should a warranty claim or service be required.

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
## 3. RECOMMENDATIONS AND WARNINGS

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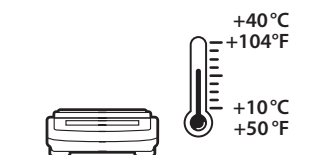
Operate the device in a room where the temperature ranges between 10 – 40 °C (50–104 °F) and where the relative humidity is below 80%. Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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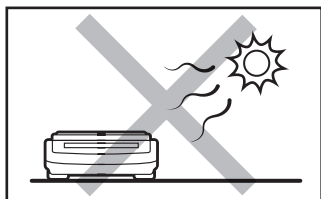
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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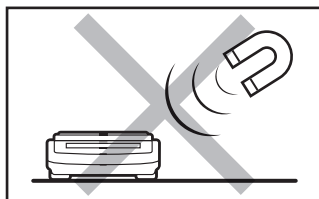
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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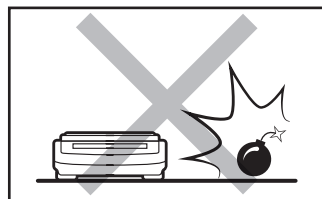
Do not place the balance in a hazardous area. Do not weigh explosive materials.

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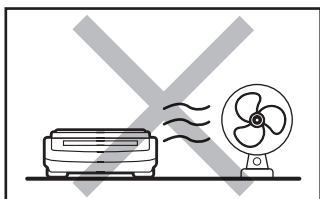
Avoid air drafts and air movements at the workstation.

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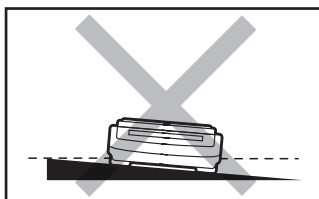Make sure that the balance is placed on an even surface.

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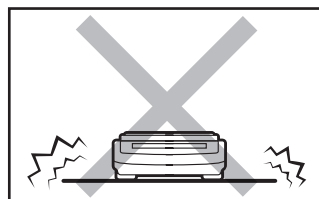Do not place the balance on an unstable ground exposed to shocks and vibrations.

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
## 4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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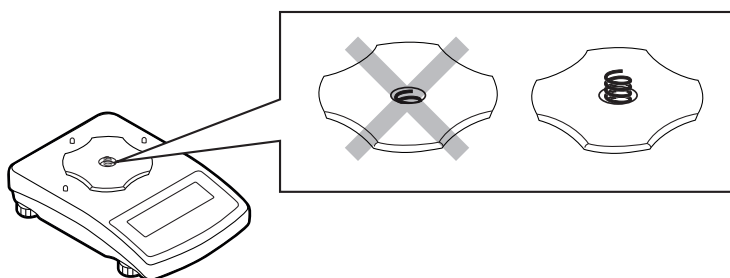### 4.1. Grounding spring check

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

Check whether the grounding spring takes its intended location. Make sure that the grounding spring juts slightly out of the hole.

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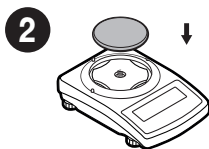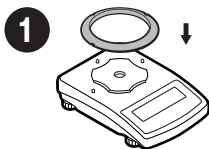
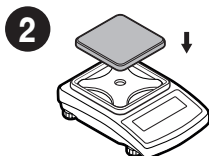
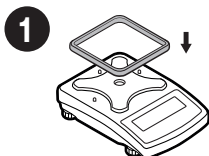
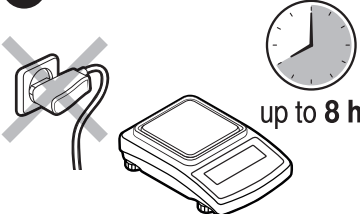
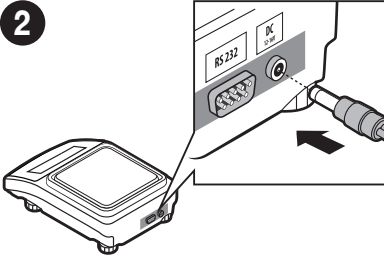
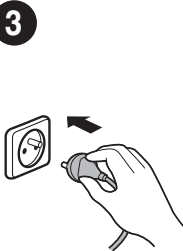
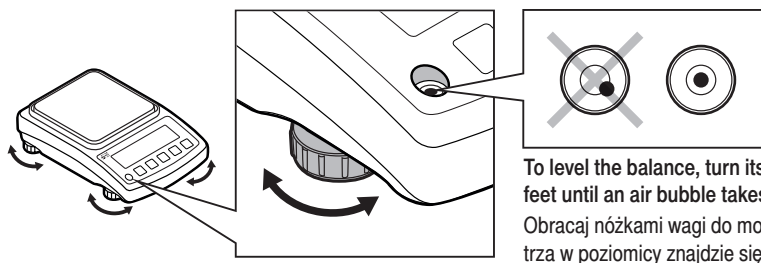
#### 6.3. Switching the balance on

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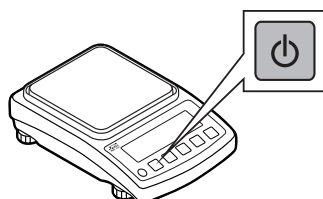Your balance is ready to use!

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To switch the device on, press ON / OFF key.

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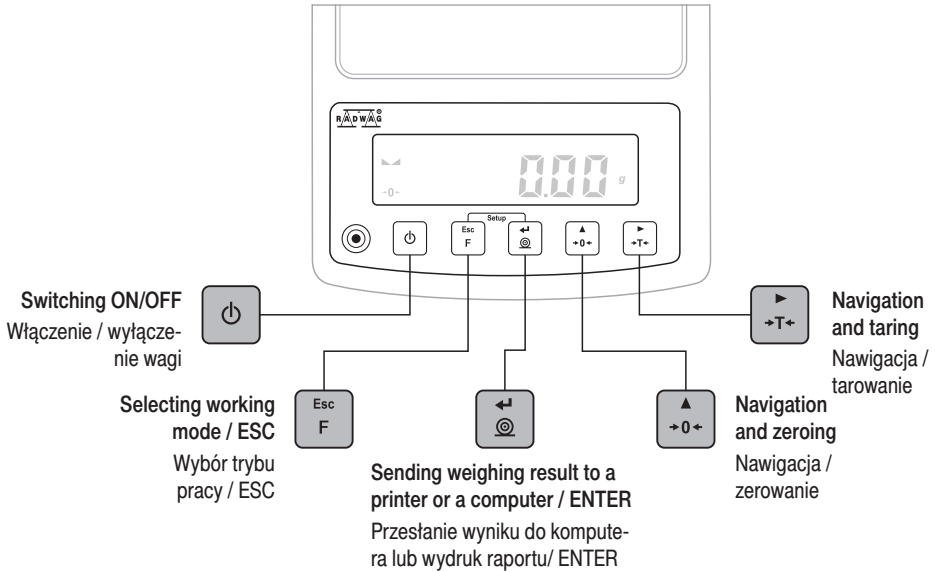
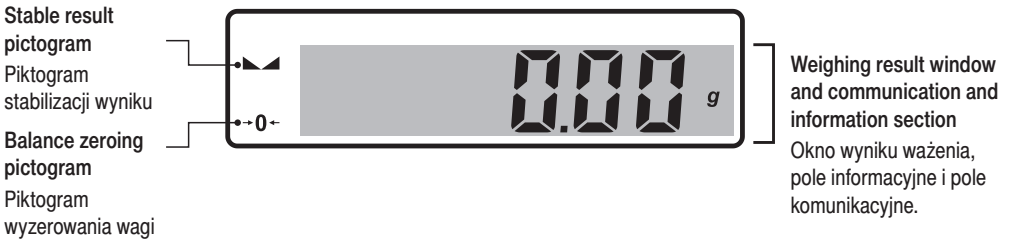
## 8.1. Adjustment

### Adiustacja

External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on balance type and capacity.

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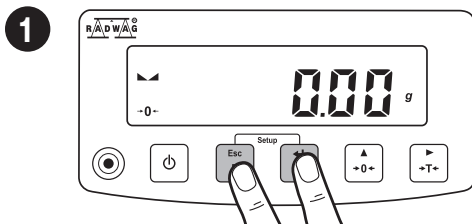
**The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use a mass standard of class F1 or F2. View the table and select the mass standard you need.**

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

Simultaneously press + keys to enter home screen.

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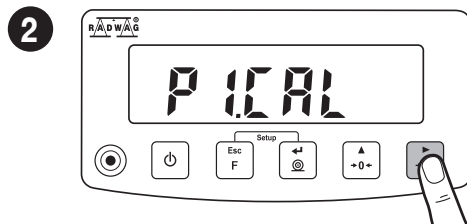
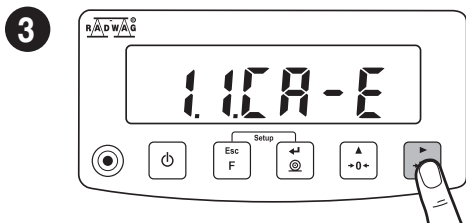
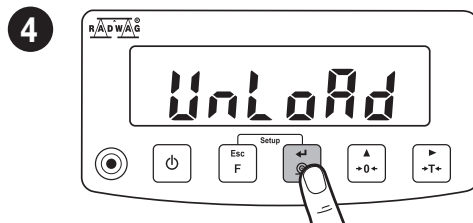
Unload the weighing pan and press key.

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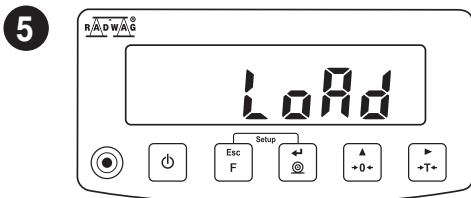
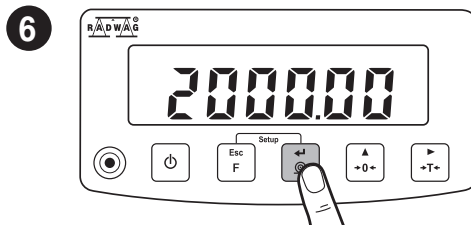
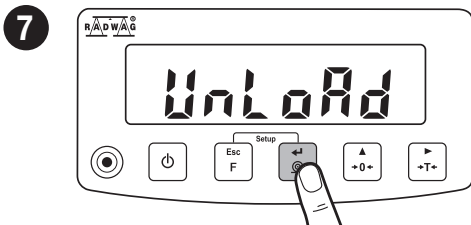
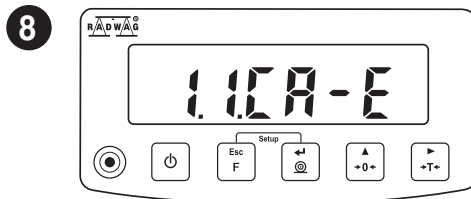
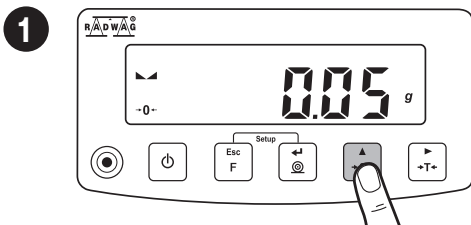
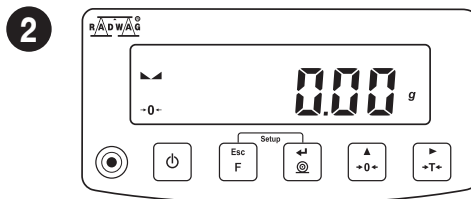
## 8.3. Taring
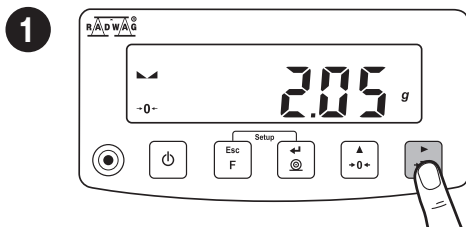
### Tarowanie

With loaded weighing pan, upon result stabilization, press key.

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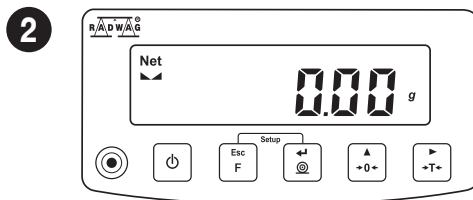
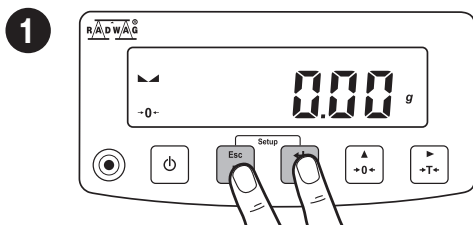
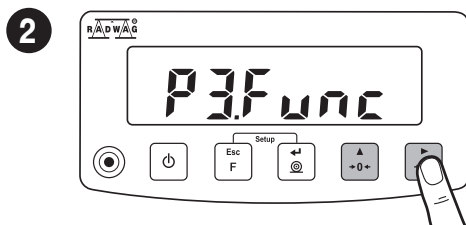
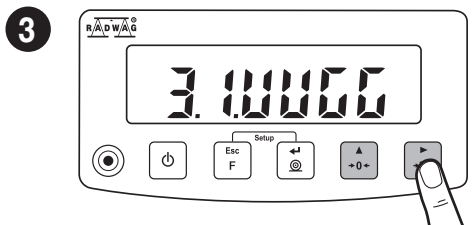
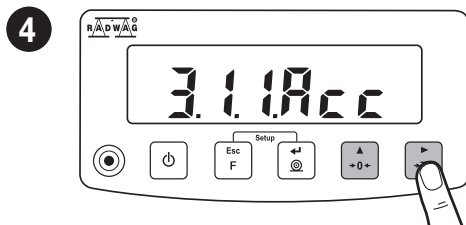
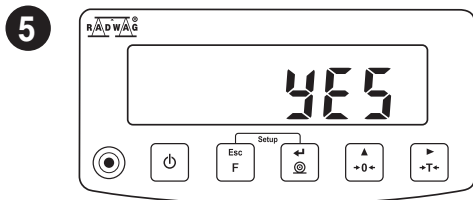
Using key select "YES" value and press key to confirm.

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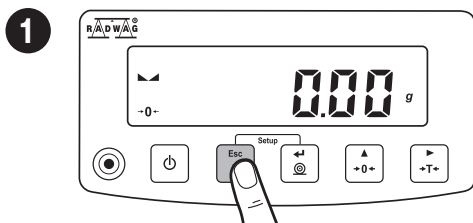
### Tryby pracy

Using key enter working modes menu.

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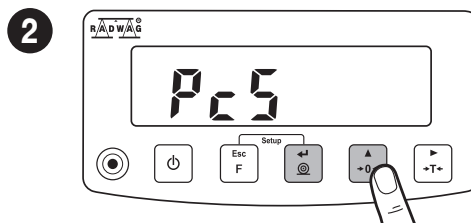Using key select working mode and press .

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### The menu features the following working modes:

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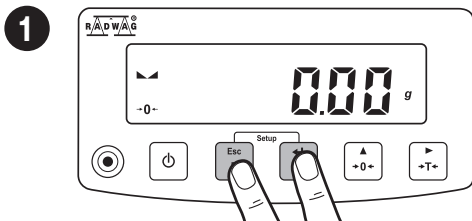
## 8.6. Weighing unit selection

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
When the main unit is g (gram), the user has the following units: g (gram), ct (karat), lb (pound).\*

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
1 Simultaneously press keys to enter home screen.

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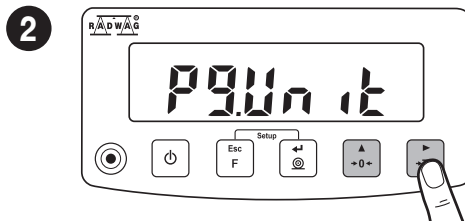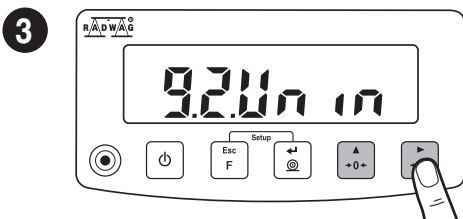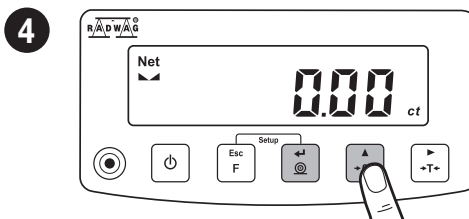
4 Using key select weighing unit and press key to confirm.

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
\*(lb) pound unit unavailable for verified balances.

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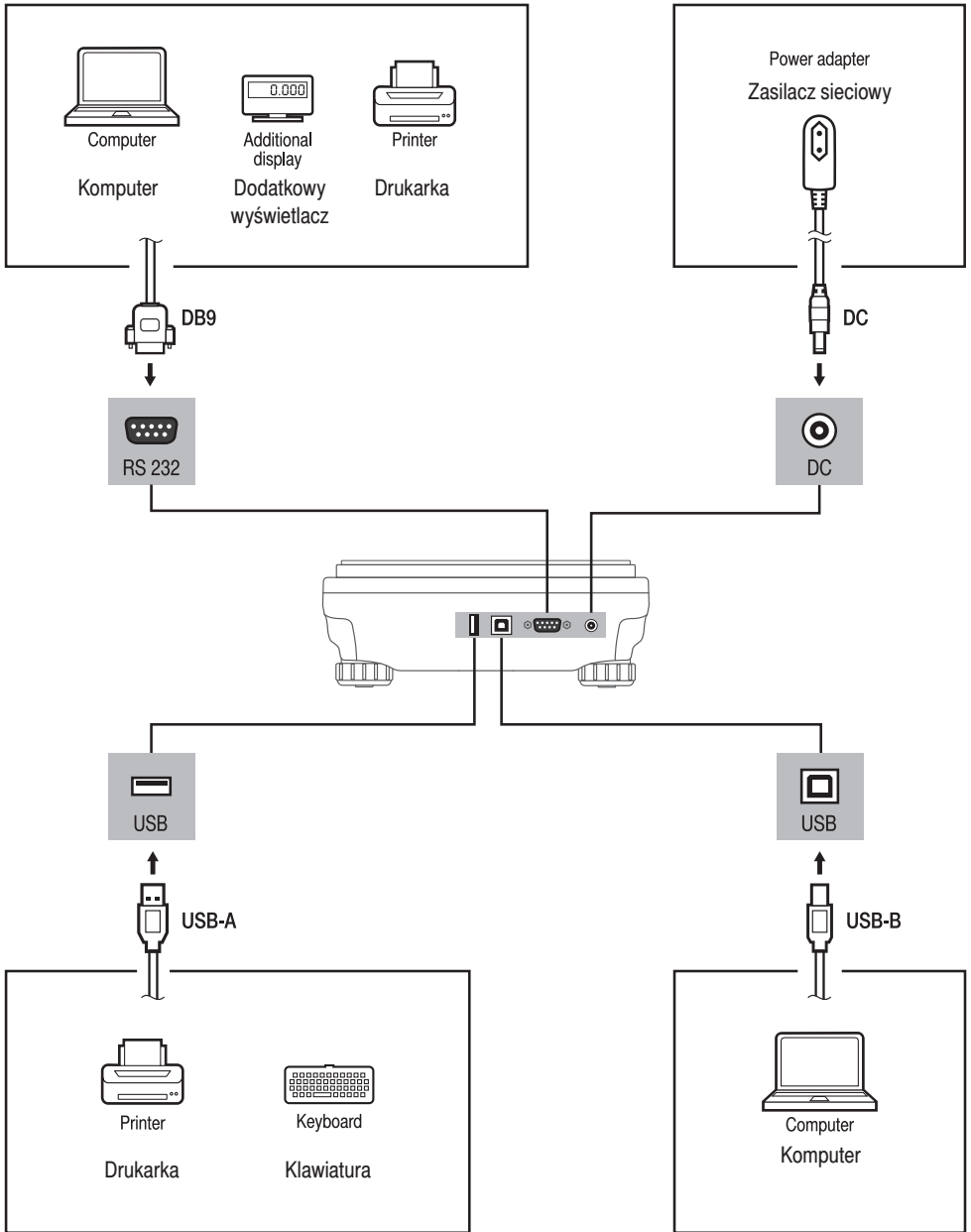
## 10. DEVICE CLEANING

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
#### 10.1. General recommendations

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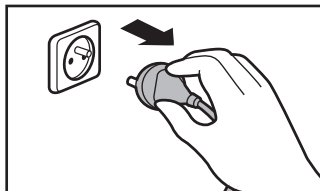
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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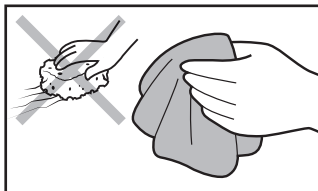Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface.

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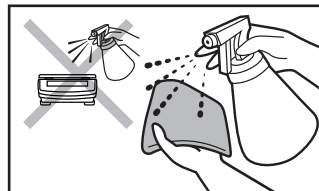Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device.

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**Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).**

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**Neither use substances of sharp structure nor intended for scrubbing.**

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**Make sure that neither dust nor liquid gets inside the weighing system (device inside).**

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