IMMB-17-01-05-23-EN-PL

Start Up Guide

Szybki start

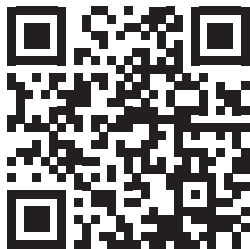
Automatic Device for Multi-channel Pipette Calibration

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- AP-12.5Y
- AP-12.1.5Y

To view the complete User Manual, go to the website or scan the QR code:

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
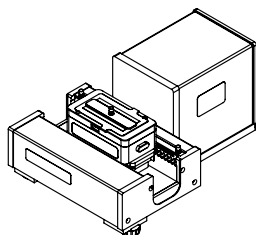
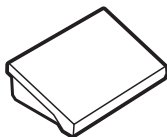
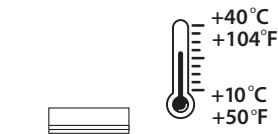
2. RECOMMENDATIONS AND WARNINGS

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Operate the device in a room where the temperature ranges between 10 – 40 °C (50 – 104 °F) and where the relative humidity is below 80%. The temperature change must occur gradually and slowly. The use of an anti-vibration table is recommended for obtaining stable and repeatable weighing results.

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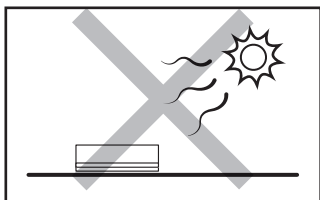
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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
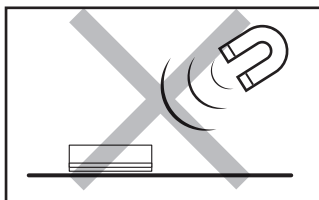
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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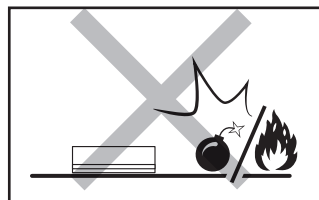
Do not place the balance in a hazardous area. Do not weigh explosive or flammable materials.

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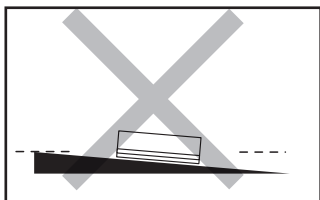
Make sure that the balance is placed on an even surface.

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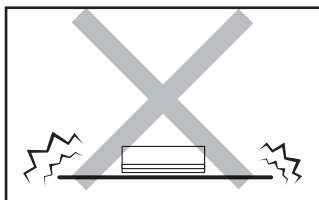Do not place the balance on an unstable surface exposed to shocks and vibrations.

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
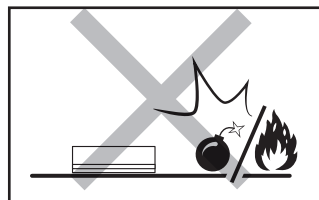
AP-12.5Y | AP-12.15Y

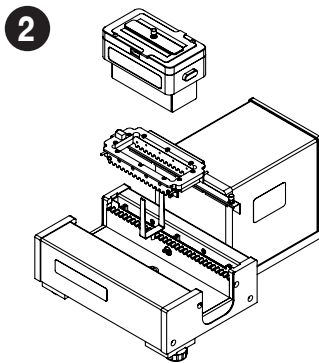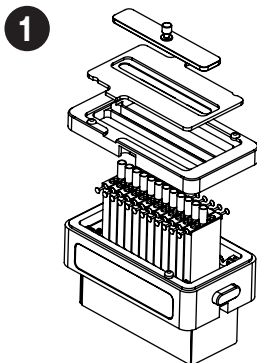4. PREPARING FOR WORK

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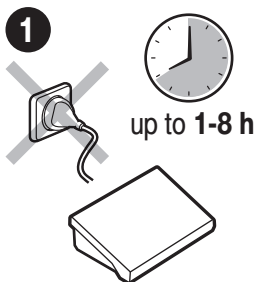
Place the device at the workstation, wait until it reaches room temperature.

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
4.1. Temperature Stabilization

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Prior to switching the device on, let it reach room temperature (estimated stabilization time: up to 8 hours).

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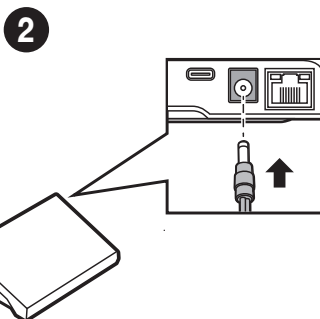
Connect the power supply cable to the DC socket on the back of the balance or on the back of the terminal.

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Do not connect the balance and the terminal to the power supply at the same time.

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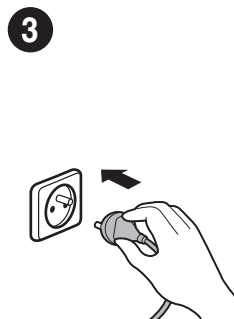
Connect the power supply to the mains.

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
4.2. Switching the Balance On

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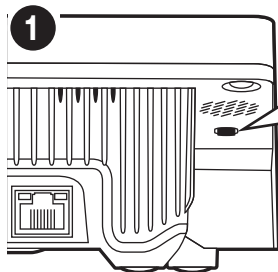
The balance switches on automatically when connected to the mains. If the balance is manually switched off using the top menu, it must be switched on using the button underneath the display housing on the left.

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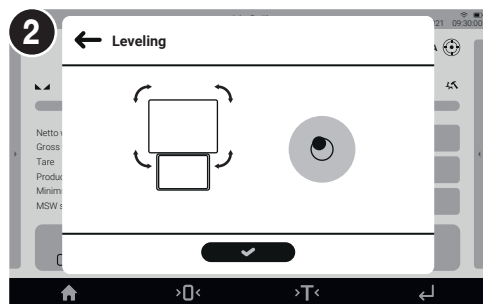The balance with automatic levelling starts the program and the levelling process runs automatically.

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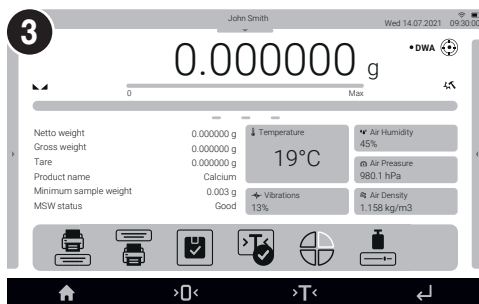
When autolevelling is complete, the home screen is displayed.

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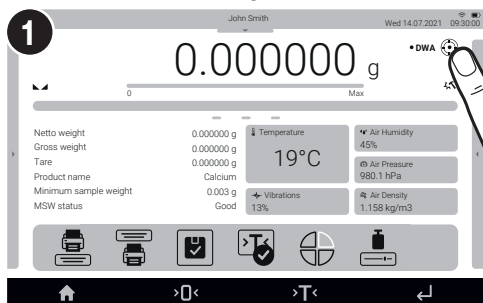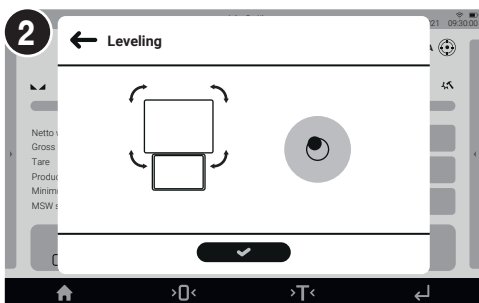
The control panel of the levelling function is displayed. To level the balance, turn its feet; keep turning until the pictogram on the screen is in the central position. Press  to return to the home screen.

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5. HOME SCREEN

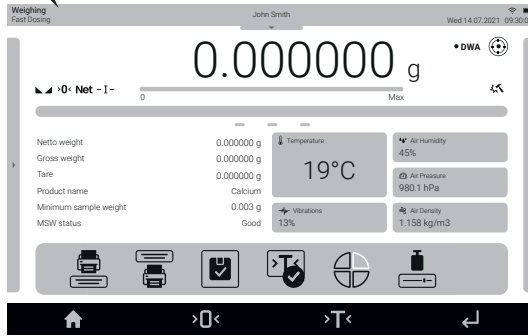
EKRAN GŁÓWNY

information about the currently used working mode, logged-in operator, date, time, active connection with a computer.

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
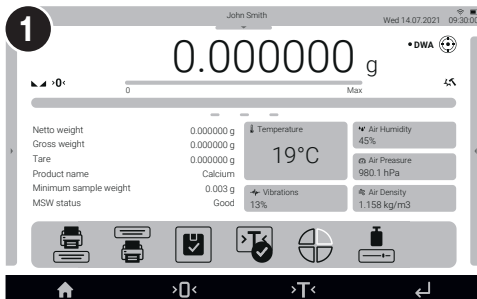
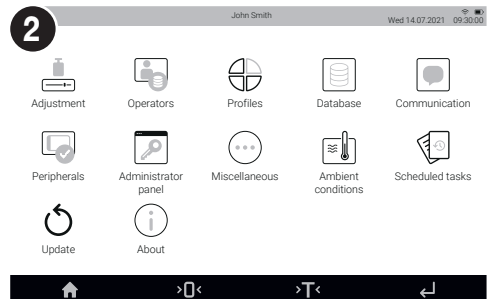
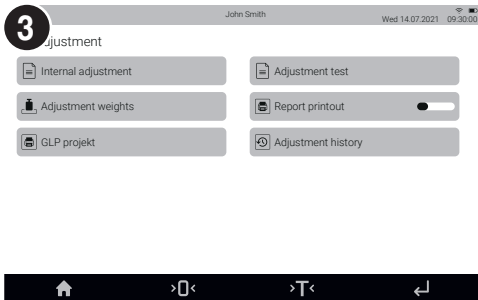
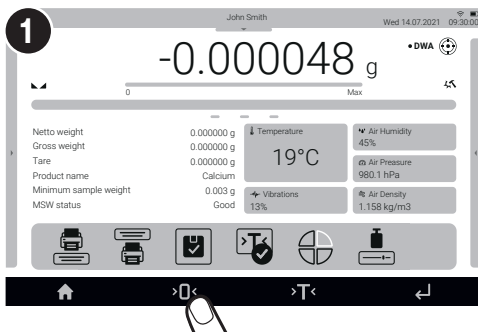
6.2. Zeroing Zerowanie


Make sure that the weighing pan is empty and press “Zeroing” button.

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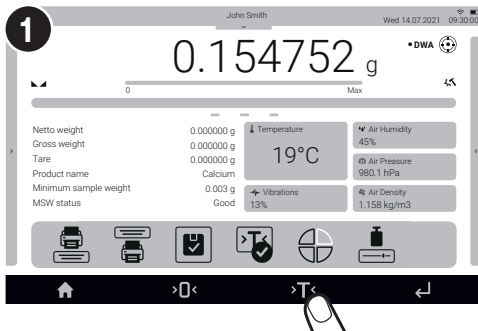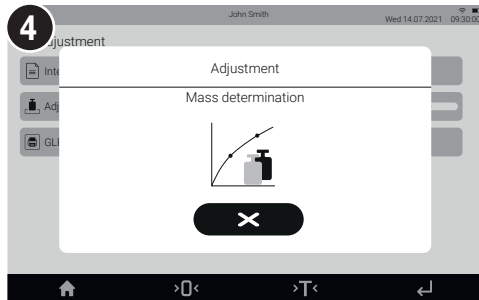
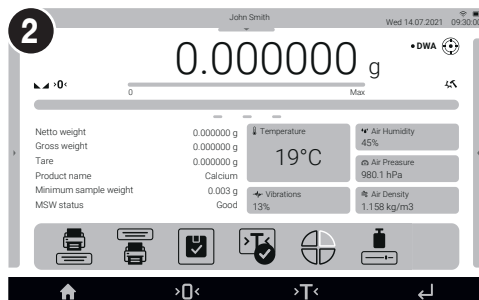
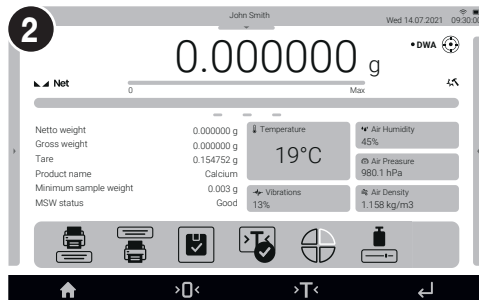
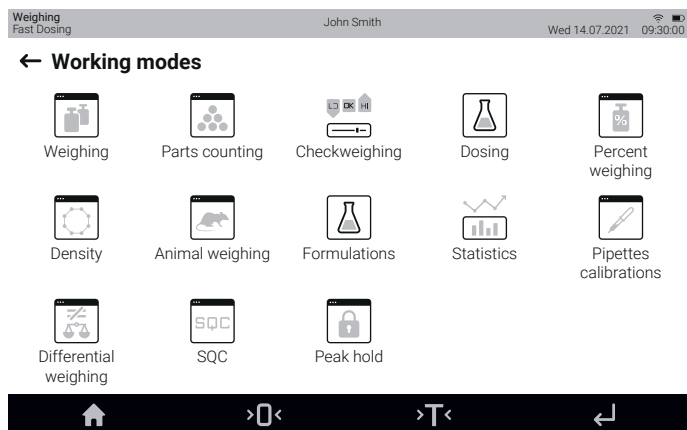
6.5. Unit Selection

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Press the current mass unit pictogram to display the list of available units.

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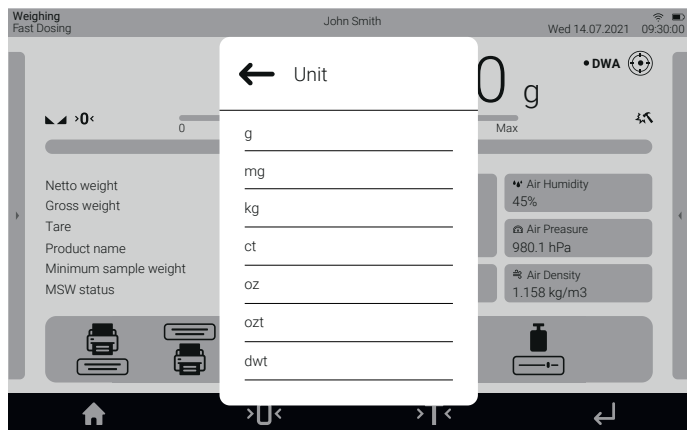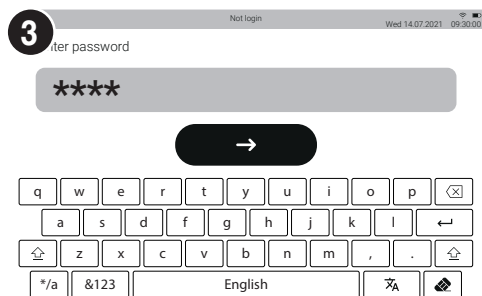
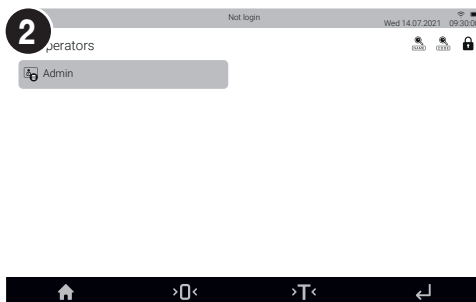
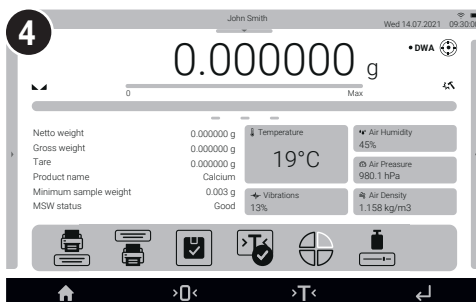
7.2. Operator Settings

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Prior to operators configuration, log in as the Administrator (see section 9.1).

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
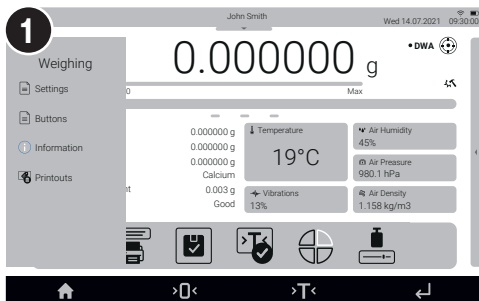
You can assign a function of your choice to each proximity sensor. In addition, you can assign various functions to the same sensor in different working modes. See the exemplary procedure for sensors configuration for the „Weighing” mode.

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Slide out the menu.

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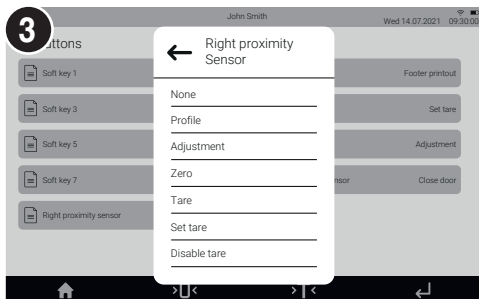
Select a respective sensor (right or left) from the list.

Select a function to be assigned to the sensor.

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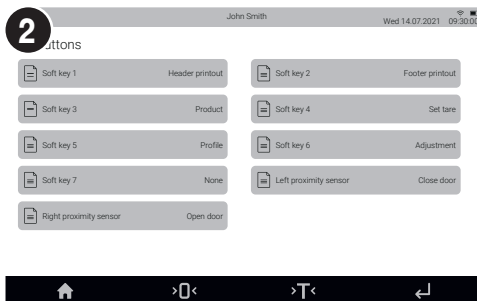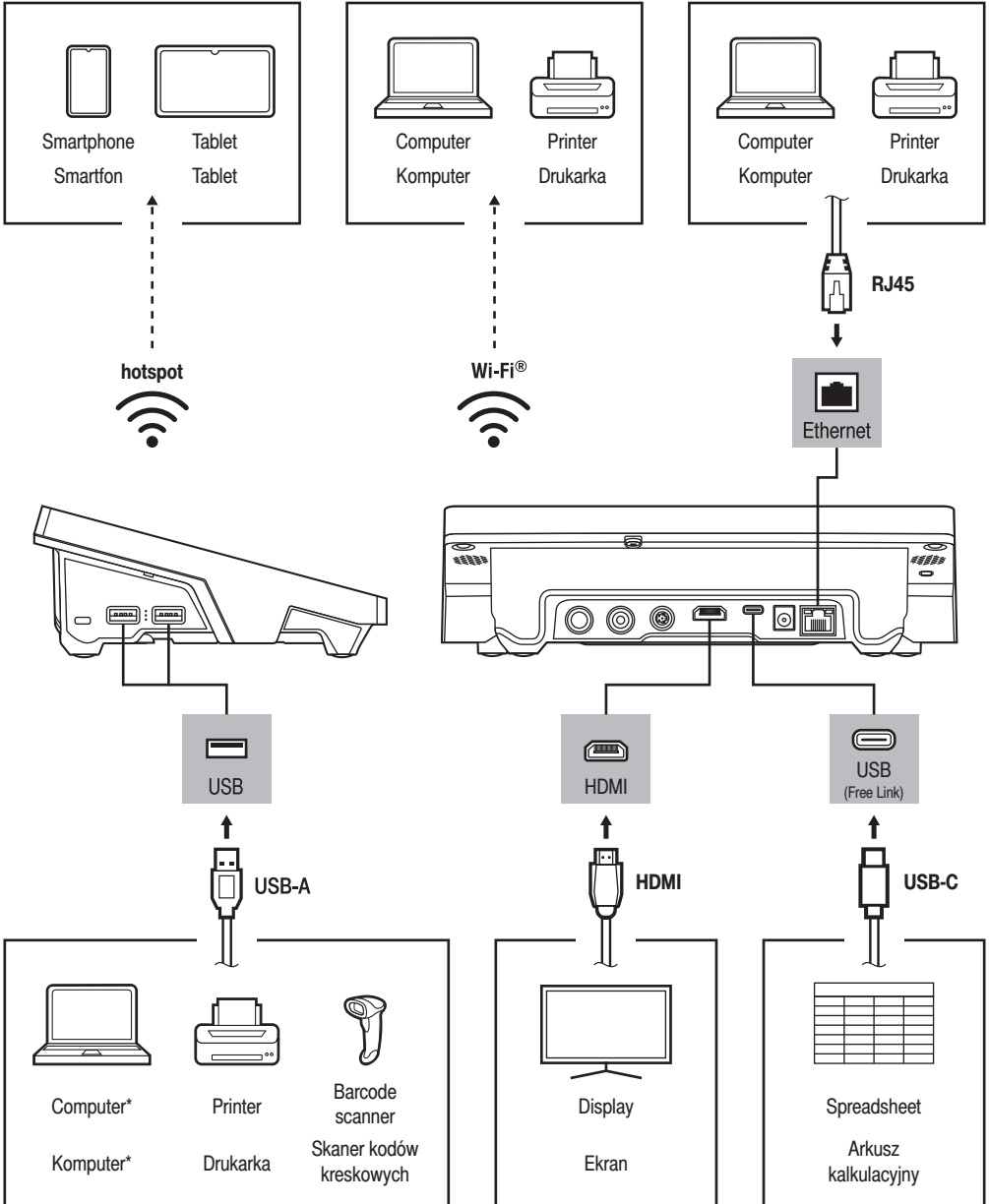
Wi-Fi® is a registered trademark of Wi-Fi Alliance.

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*A USB to RS232 converter must be used for proper connection.

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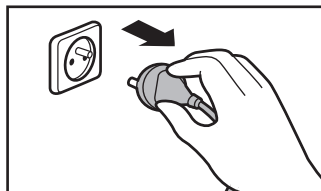
9. CLEANING AND MAINTENANCE

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
9.1. Guidelines and Precautions

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
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet and disconnect all cables.

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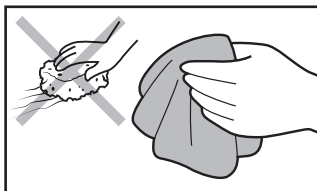Use soft cloth made of either natural or synthetic fibers exclusively. Avoid using stiff or hard fabric, as this may cause scratches on the surface!

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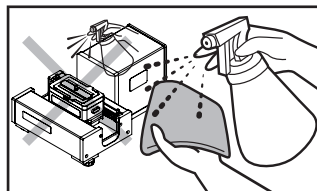Apply the cleanser to the cloth. Remember not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Use neither sharp-structured nor scrubbing-purposed substances.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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If it is necessary to disassemble the weighing chamber, follow the instructions in section 11.3.

- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either a brush or a handheld vacuum cleaner.
- Clean components can be dried using a soft cloth or a dust-free paper towel. With this, the remaining moisture will be absorbed.
- **Install all the components, providing they are completely dry.**

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