IMMB-05-02-05-22-EN-PL

Start up Guide

Szybki start

Precision Balances

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- PM 4Y
- WLY

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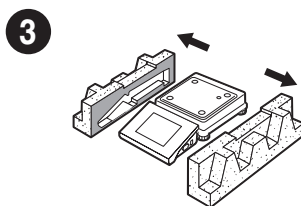
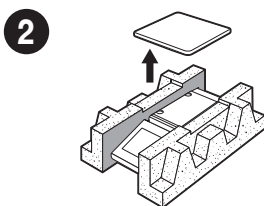
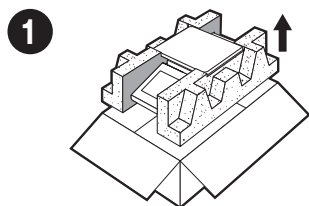
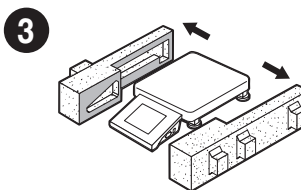
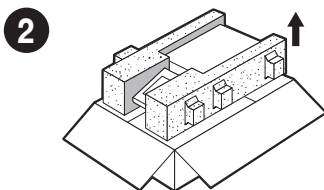
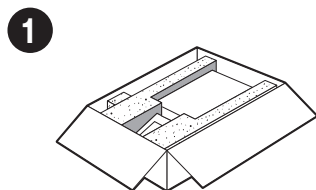
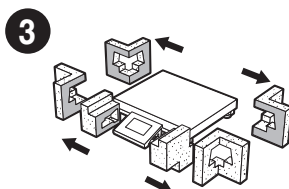
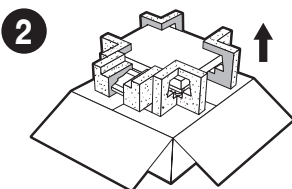
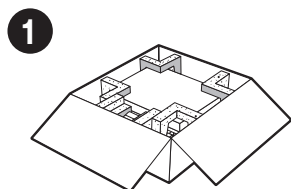
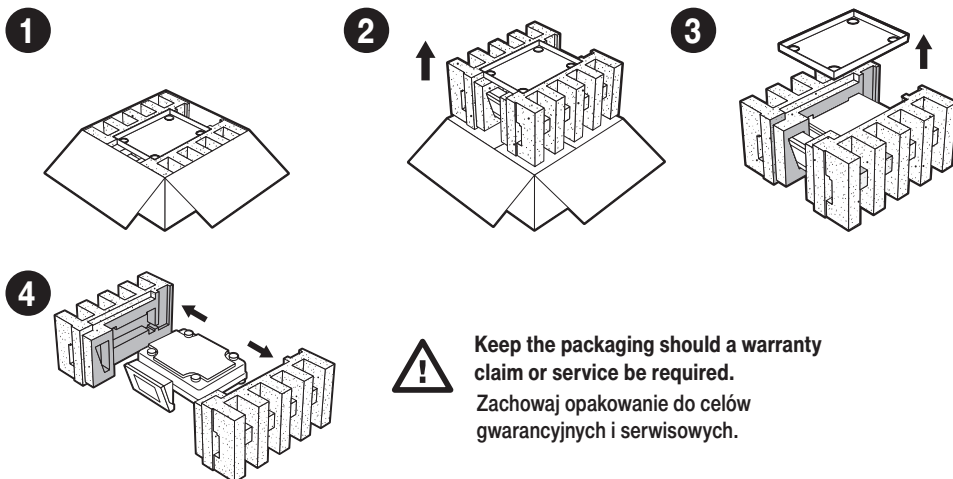
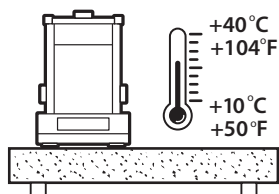
PM 4Y ($d \leq 0,1 \text{ g}$)

3. RECOMMENDATIONS AND WARNINGS

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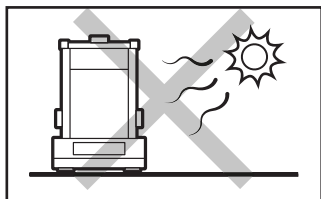
Operate the device in a room where the temperature ranges between 10–40 °C (50 – 104 °F) and where the relative humidity is below 80 %. Temperature must change slowly and gradually.

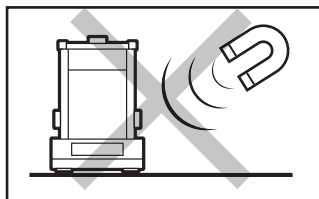
Using the anti-vibration table is recommended in order to enable the most stable and repeatable measurement.

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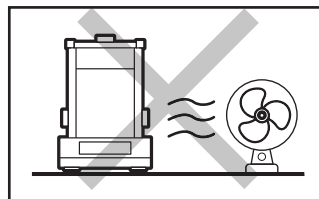
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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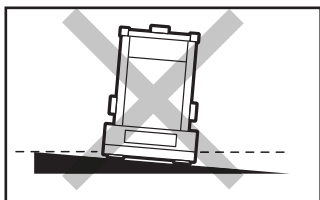
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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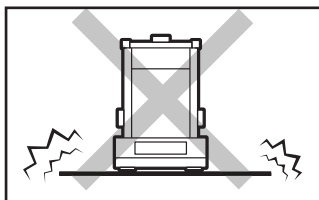
Avoid air drafts and air movements at the workstation.

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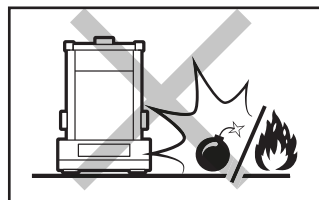
Make sure that the balance is placed on an even surface.

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Do not place the balance on an unstable surface exposed to shocks and vibrations.

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Not intended for usage within EX zones. Not intended for weighing explosive or flammable substances.

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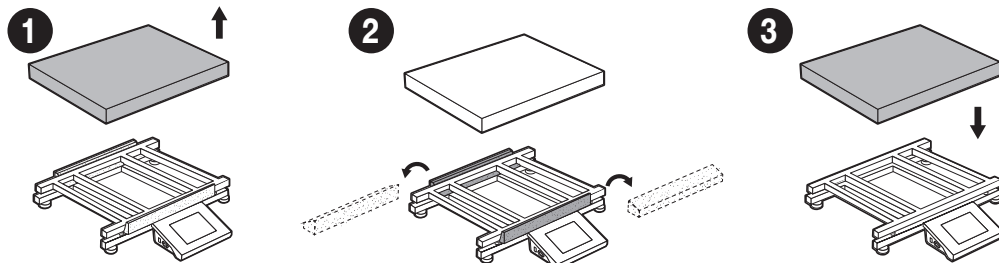
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Transport lock removal

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WLY F1 | WLY C2 | PM 4Y ($d \geq 0,5$ g)

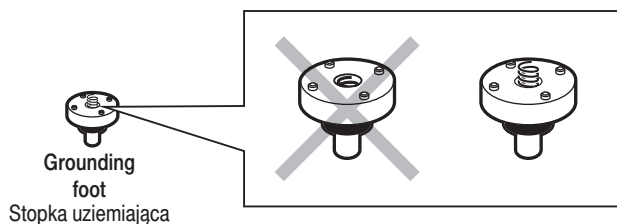
Keep the transport lock should a warranty claim or service be required.

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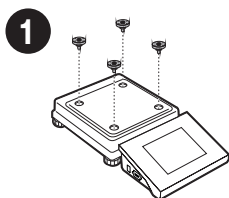
5. COMPONENTS ASSEMBLY

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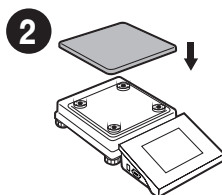
WLY D2

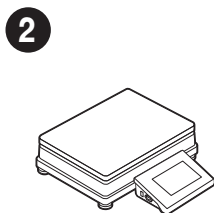
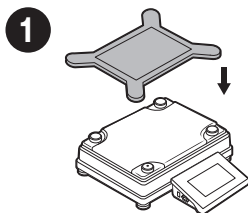
Install the grounding foot in a freely selected opening.

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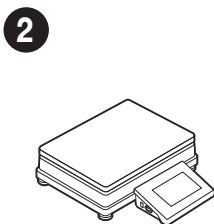
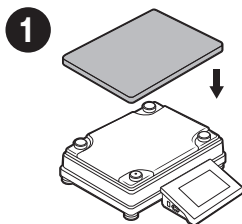
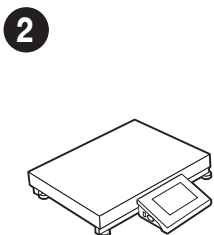
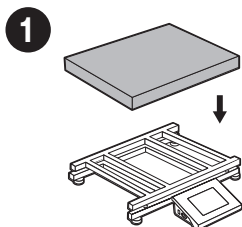
WLY F1 | WLY C2 | PM 4Y (d ≥ 0,5 g)

6. PREPARING FOR WORK

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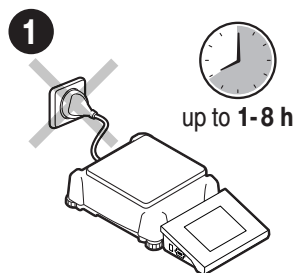
Place the device at the workstation, wait until it reaches room temperature.

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6.1. Temperature stabilization

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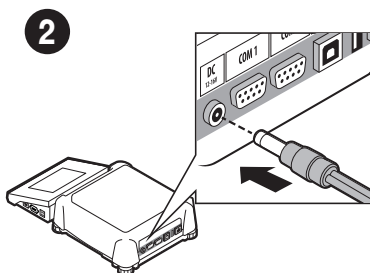
WLY D2

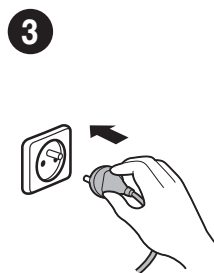
Prior switching the device on let it reach room temperature (estimated stabilization time: up to 8 hours).

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Connect the power adapter to DC connector.

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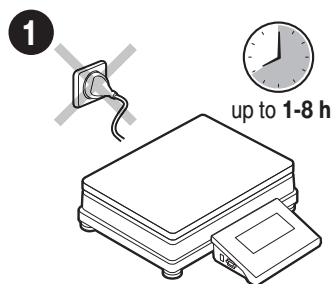
Connect the power adapter to the mains.

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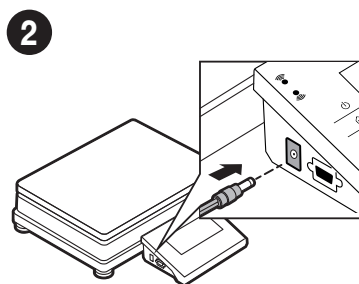
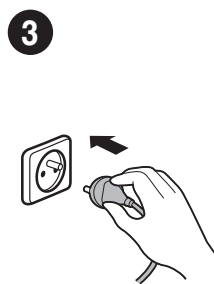
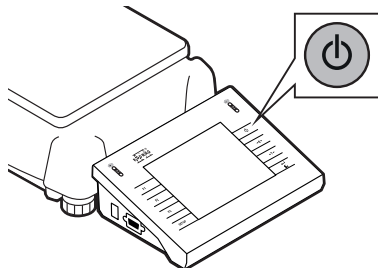
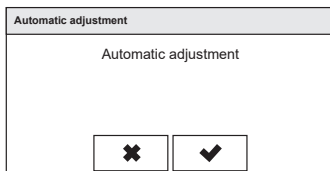
Confirm or cancel the start of the automatic adjustment process.

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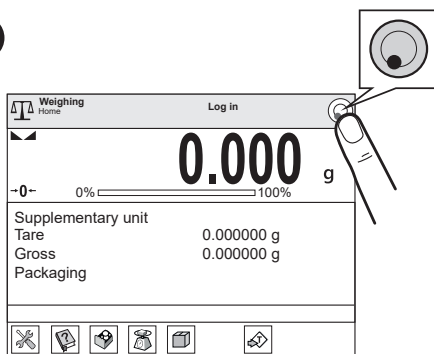
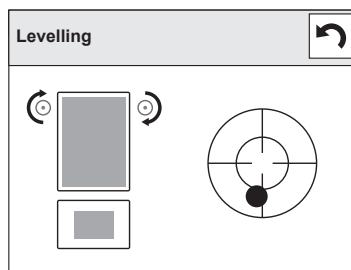
Press "Levelling" icon.

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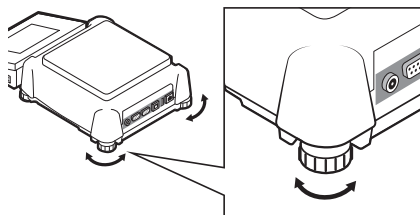
2

Check the balance level state. The arrows shows how to move the balance feet to level the balance.

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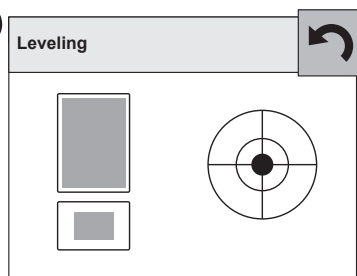
3

To level the balance turn its feet, keep turning the feet until the level indicator takes central position.

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4

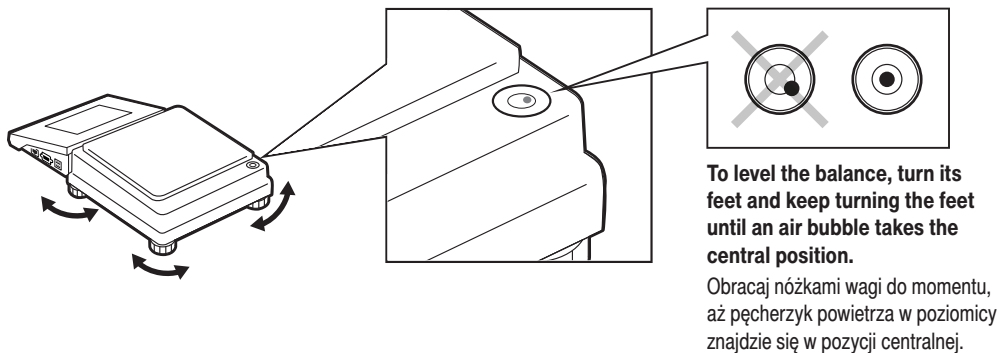
Balance has been levelled. To close the levelling window press  button.

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WLY D2

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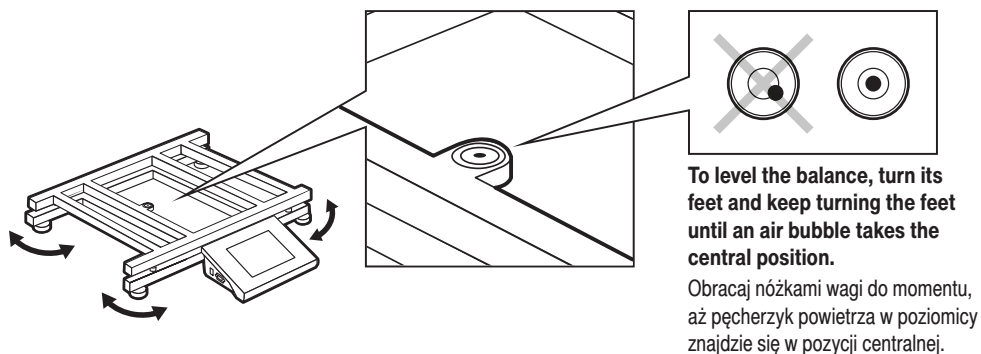
Before levelling the balance, remove the weighing pan. After levelling put the weighing pan again. The operation described in pt. 4

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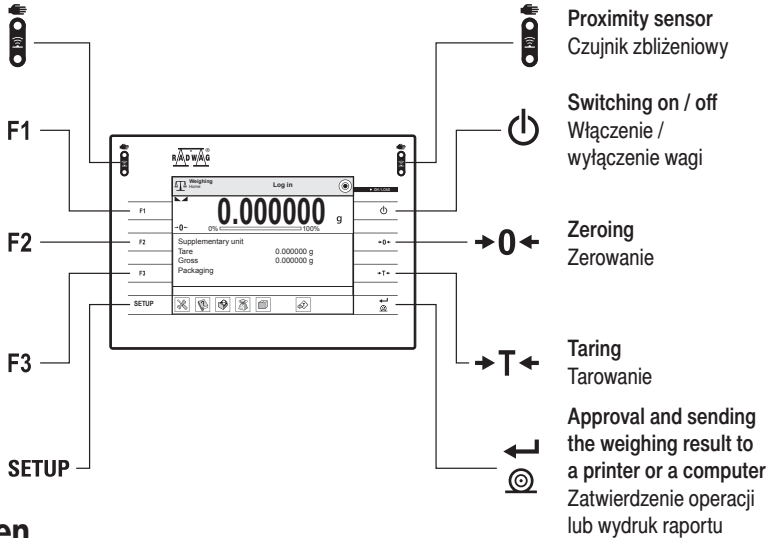
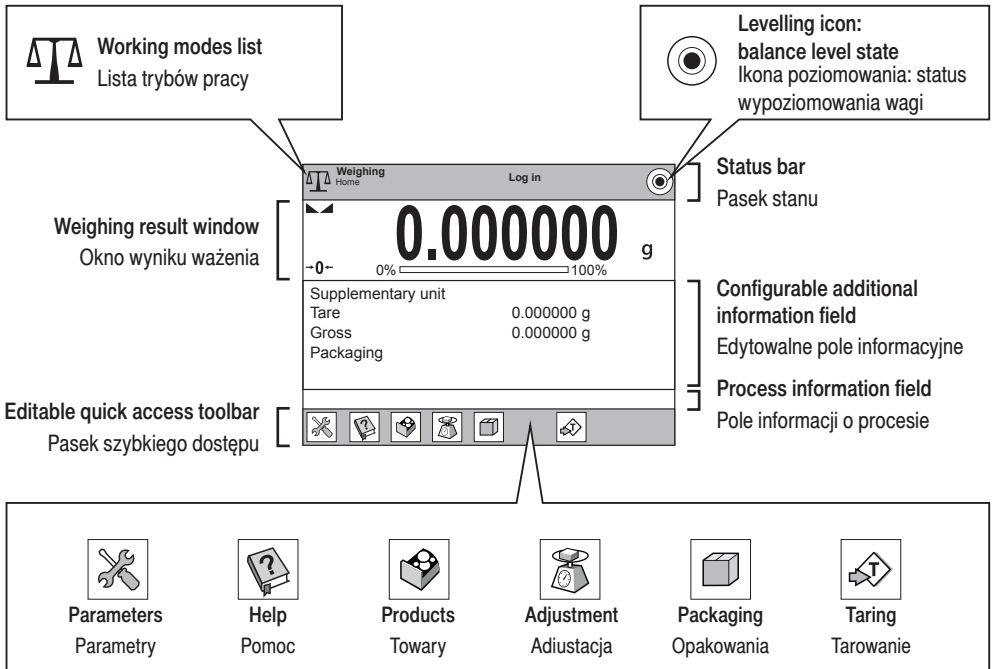
Ekran główny

8. BASIC OPERATIONS

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8.1. Adjustment (for models with external adjustment)

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External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on balance type and capacity.

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The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use a mass standard of class F1 or F2. View the table and select the mass standard you need.

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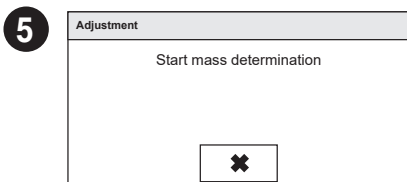
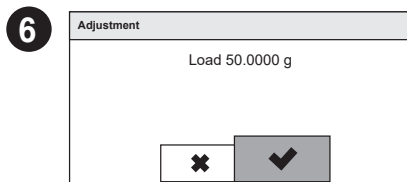
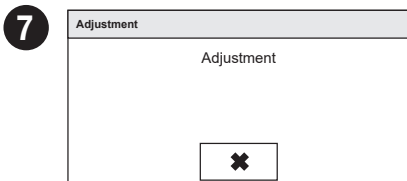
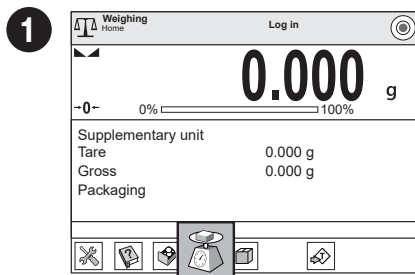
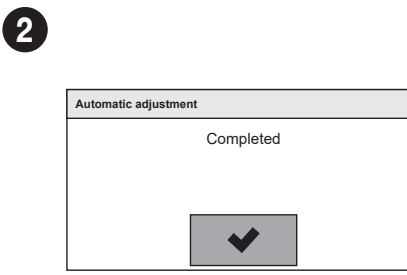
8.2. Adjustment (for models with automatic adjustment)

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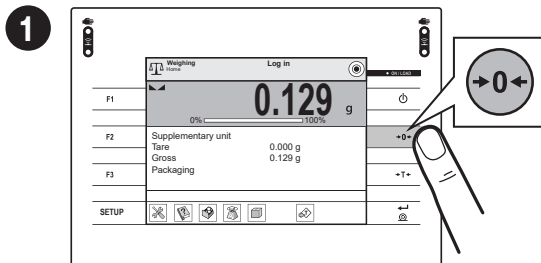
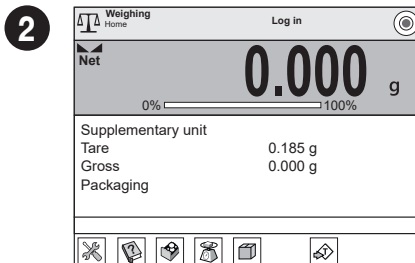
8.3. Zeroing

Zerowanie

Make sure that the weighing pan is empty and press “Zeroing” button.

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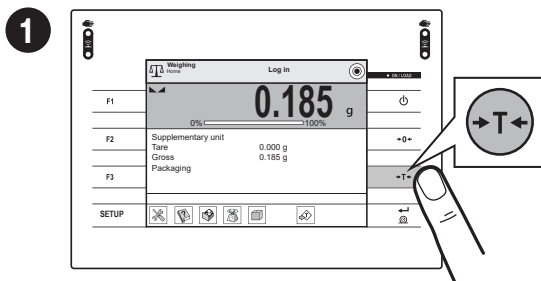
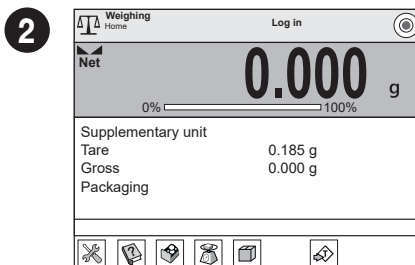
8.4. Taring

Tarowanie

With loaded weighing pan: upon stabilisation of the weighing result, press “Tare” button.

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8.5. Working modes / units selection

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Working modes

1	Weighing
2	Parts counting
3	Checkweighing
4	Dosing

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0.000 g

Supplementary unit
 Tare 0.000 g
 Gross 0.000 g
 Packaging

Press the current weight unit icon to display list of available units.

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8.6. Working modes settings

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1

Press the icon of the current working mode, to display the list of available working modes.

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2

Press information field to access current working mode options.

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3

Press selected button to view the list of available options.

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9. SETTINGS

USTAWIENIA

Some settings are accessible for Administrator exclusively.

Prior to balance parameters setup, log in as the Administrator.

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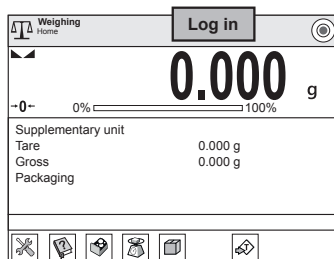
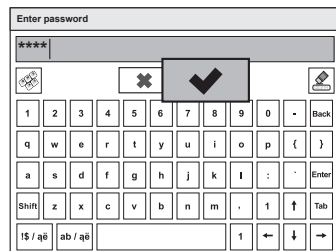
3

Enter "1111" password and press button to confirm.

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9.2. User settings

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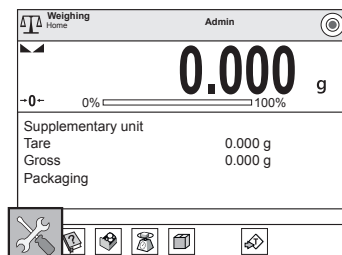
Prior to users settings, log in as Administrator (point 9.1.).

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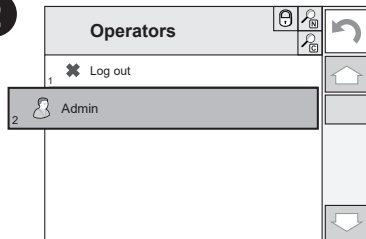
Press "Parameters" button.

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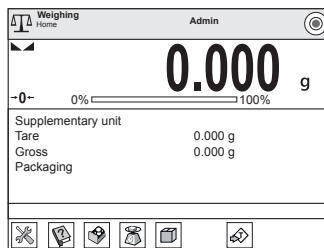
2

Users list is displayed, select "Admin" option.

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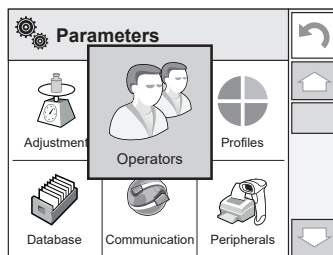
4

You have logged in as the Administrator.

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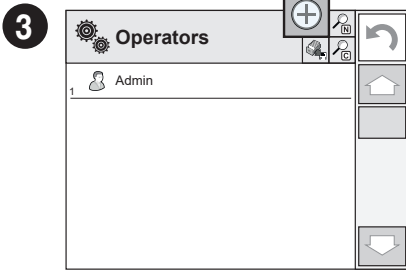
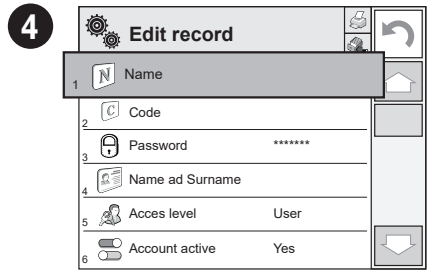
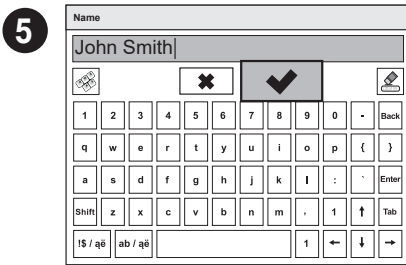
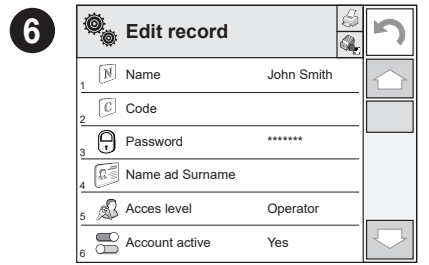
Password: Enter user password

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First and last name: Enter user's first and last name

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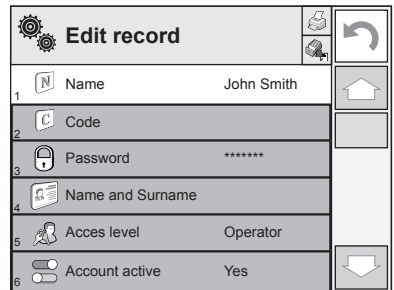
Access levels: Set respective access level
(user, advanced, admin)

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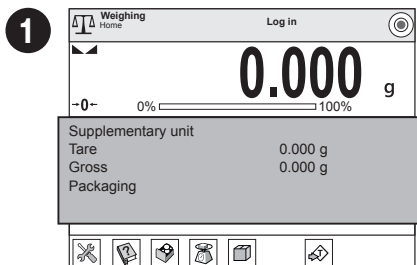
Prior to proximity sensors setup, log in as Administrator (point 9.1).

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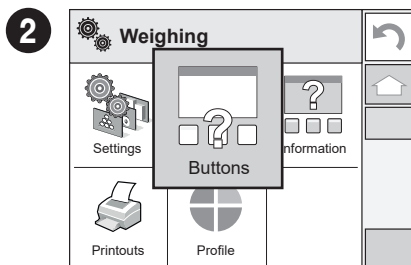
You can assign the proximity sensors with different functions for each working mode (the sensors will trigger different operation for each working mode). See the below procedure of sensors setup for the "Weighing" mode.

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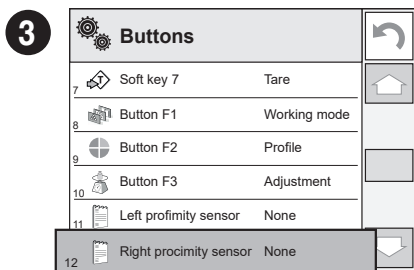
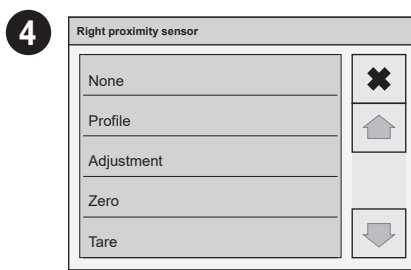
Press "Buttons" button.

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Select the appropriate sensor (right or left).

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Select function that is to be assigned to the chosen sensor.

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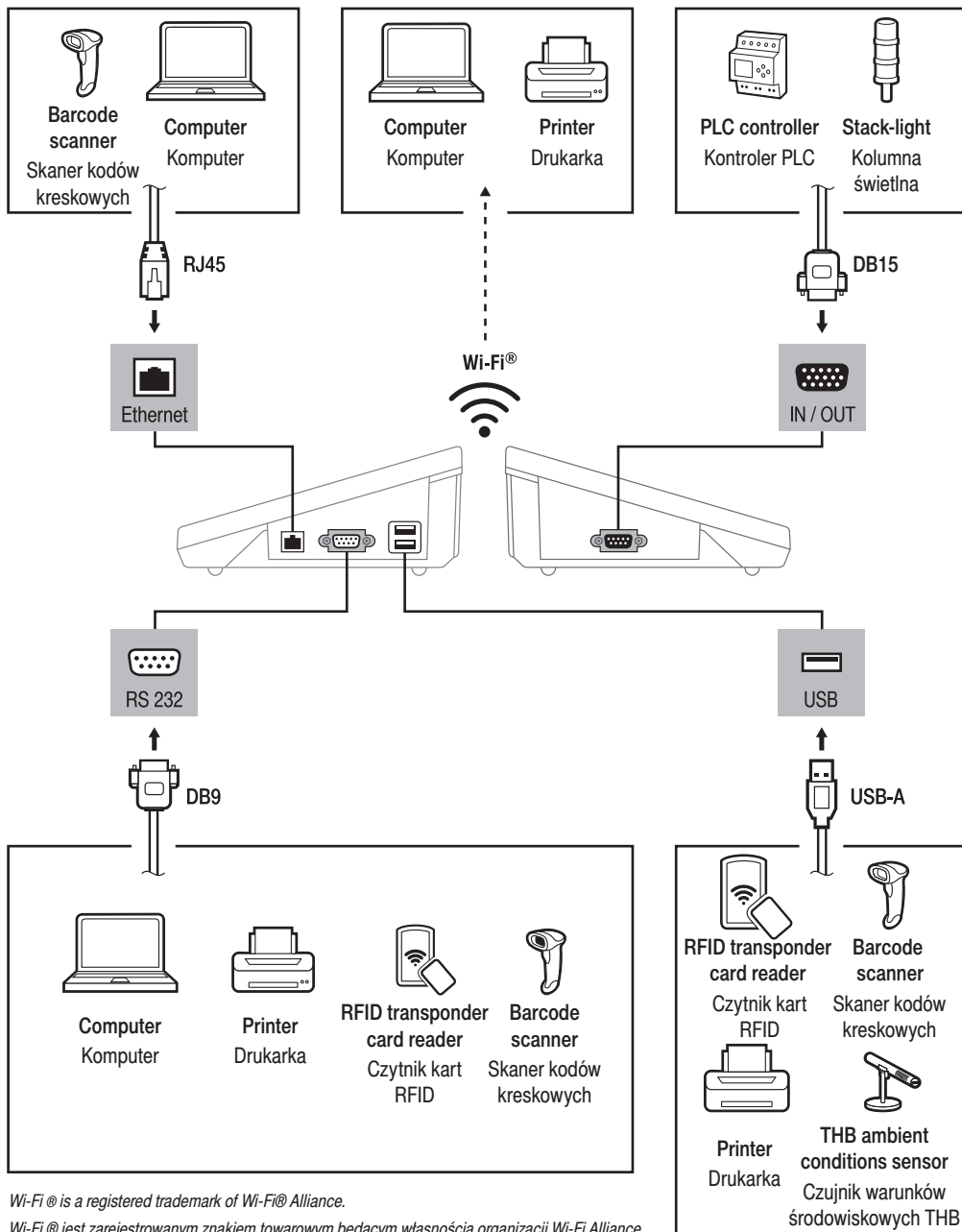
Following the above procedure you can configure sensors for other working modes.

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Wi-Fi® is a registered trademark of Wi-Fi® Alliance.

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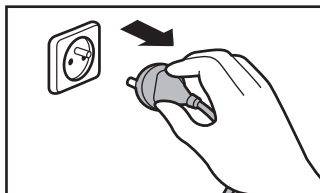
11. DEVICE CLEANING

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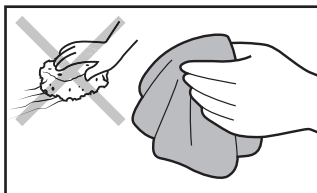
11.1. Guidelines and precautions

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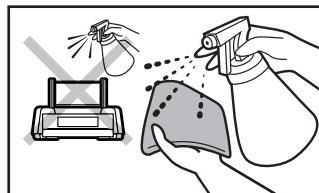
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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Use soft cloth made of either natural or synthetic fiber exclusively. Avoid using cloth made of stiff/hard fabric, this could cause scratches on the surface!

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Apply the cleanser onto the cloth. Mind not to apply the cleanser directly onto the device!

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Shall it be necessary to disassemble the weighing chamber, follow the description provided in section 11.3.

- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).

- Dry sample leftovers can be removed using either brush or handheld vacuum cleaner.
- Clean components can be dried using soft cloth or dust-free paper towel, with this the remaining moist will be absorbed.
- Install all the components providing they are completely dry.

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