IMMB-11-02-05-22-EN-PL

Start up Guide

Szybki start

Moisture Analyzer

Wagosuszarki

- MA X2.A
- MA X2.IC.A

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

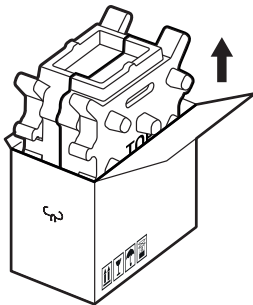
2

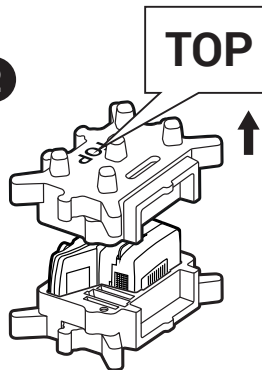3

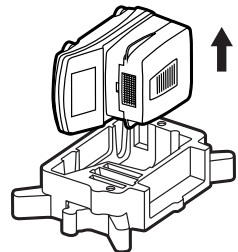
Keep the packaging should a warranty claim or service be required.

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3. RECOMMENDATIONS AND WARNINGS

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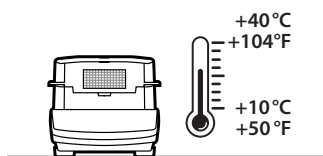
Operate the device in a room where the temperature ranges between 10–40 °C (50–104 °F) and where the relative humidity is below 80%.


Place the moisture analyzer on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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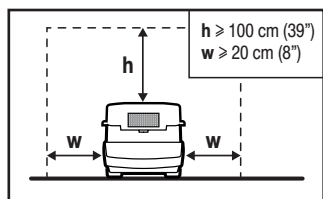
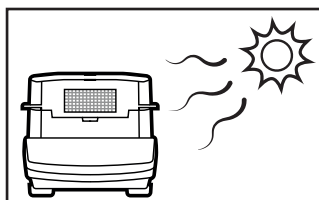
Place the device away from heat sources. Avoid exposing the device to the sunlight.

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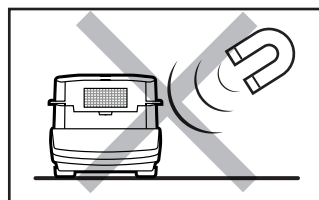
Avoid exposing the device to a magnetic field. Do not weigh magnetic substances.

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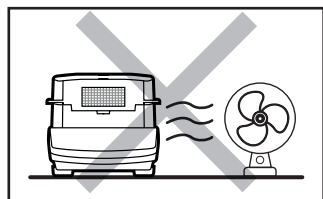
Avoid air drafts and air movements at the workstation.

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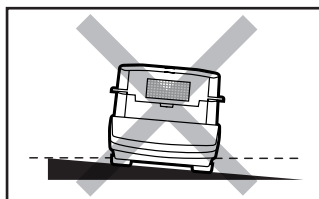
Make sure that the device is placed on an even surface.

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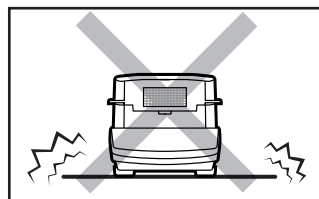
Do not place the device on an unstable surface exposed to shocks and vibrations.

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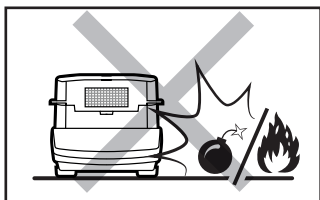
Do not place the device in a hazardous area. Do not weigh explosive or flammable materials.

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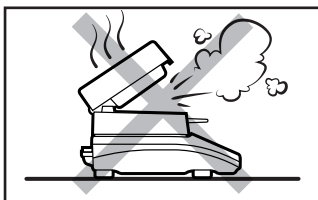Do not open in the course of operation since the heating components reach very high temperature.

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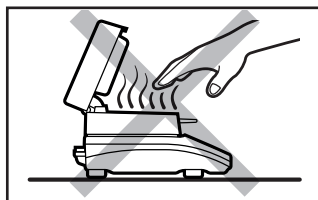Be extremely careful while taking the sample out. The sample and the moisture analyzer elements may be hot.

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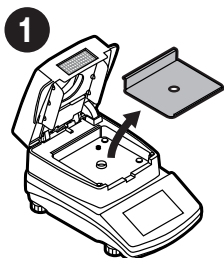
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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
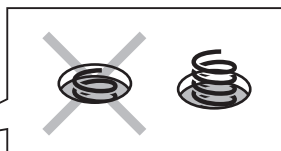
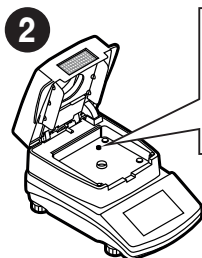
Check whether the grounding spring takes its intended location. Make sure that the grounding spring juts slightly out of the hole.

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


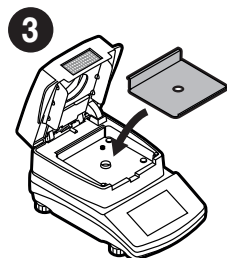
Install the base insert.

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

5. COMPONENTS ASSEMBLY

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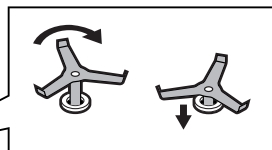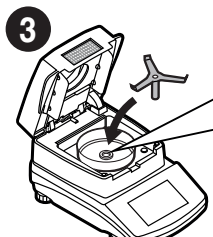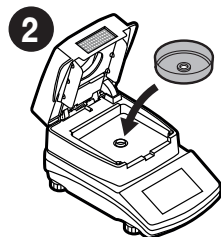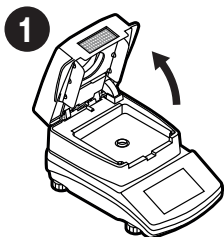Slowly rotate the cross-shaped holder, keep rotating until the holder sinks deeper.

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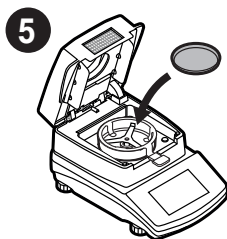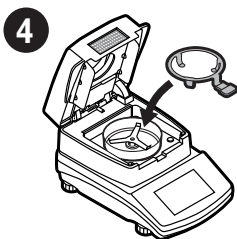6. PREPARING FOR WORK

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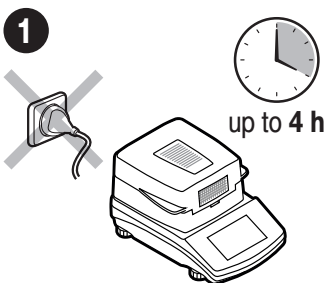
Place the moisture analyzer at the workstation, wait until it reaches room temperature.

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
6.1. Temperature stabilisation

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
Prior to switching the device on let it reach room temperature (estimated stabilisation time: up to 4 hours).

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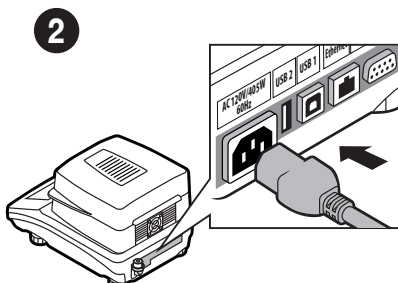Connect the power cord to the AC socket located at the back of the device.

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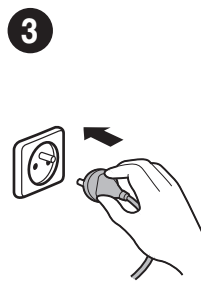
6.2. Levelling

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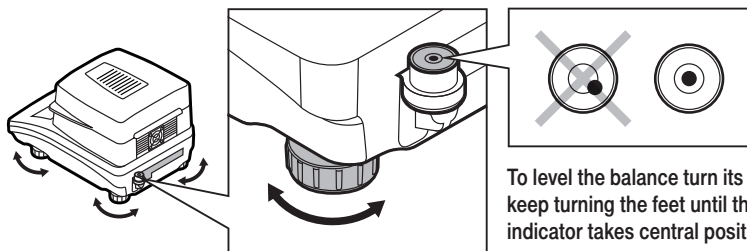
To level the balance turn its feet, keep turning the feet until the level indicator takes central position.

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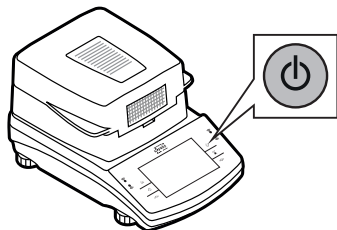Before using WEIGHING mode, carry out instrument adjustment: refer to section 8.1. (adjustment is not required for DRYING mode).

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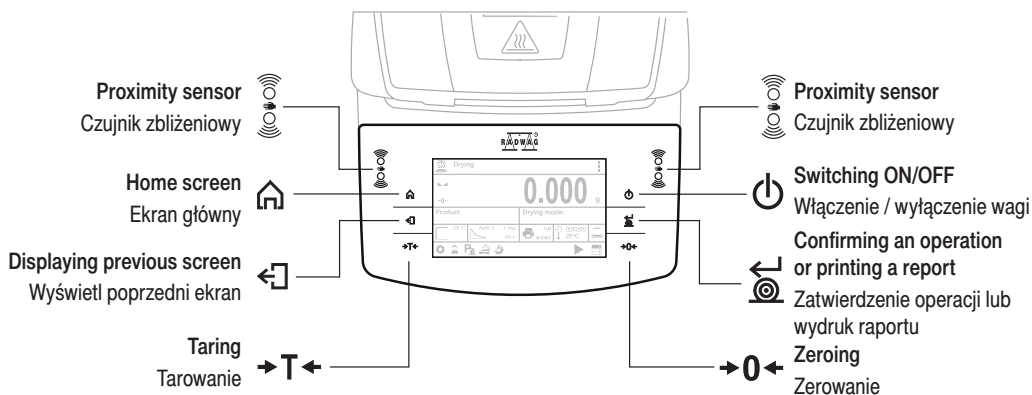
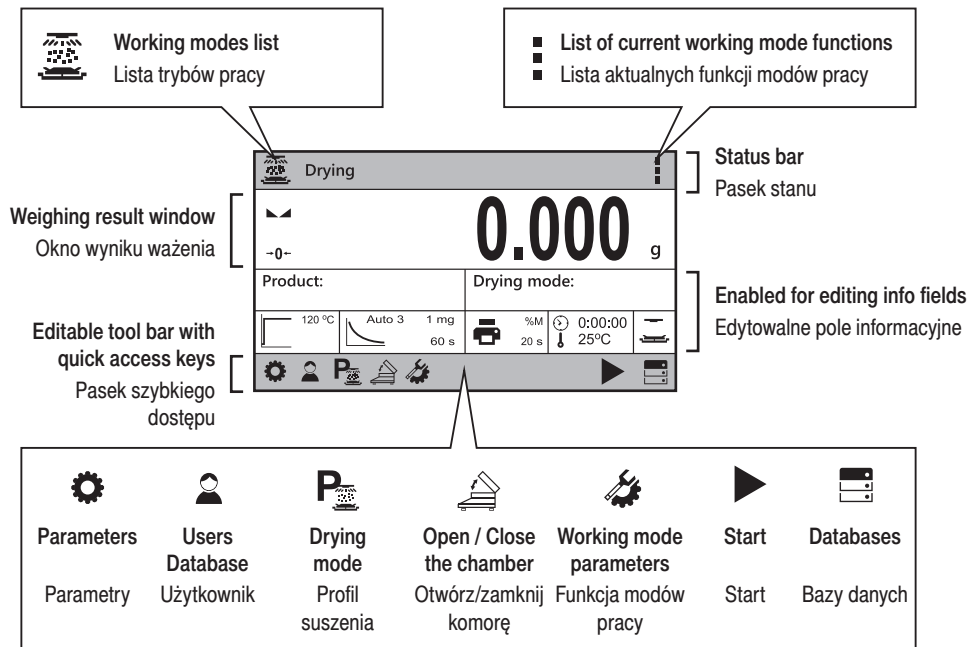
Ekran główny


8. BASIC OPERATIONS

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
8.1.1. Adjustment (for models with an internal adjustment only)

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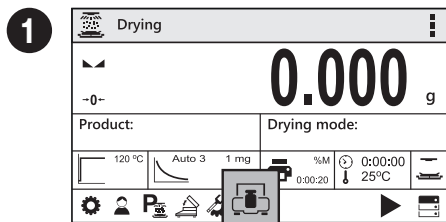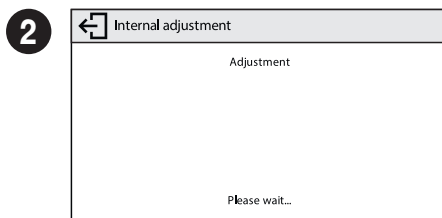8.1.2. Adjustment (for models with an external adjustment only)

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It is recommended to carry out adjustment operation prior to running the WEIGHING mode (adjustment is not required for DRYING). External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on device type and capacity.

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

The adjustment mass standard does not come with the moisture analyzer. Prior to adjustment prepare the right mass standard. You are recommended to use mass standard of class F1 or F2. View the table and select the mass standard you need. **Caution!** Make sure that the disposable weighing pan has been installed (read point 5.5.). Do not close the drying chamber in the course of adjustment.

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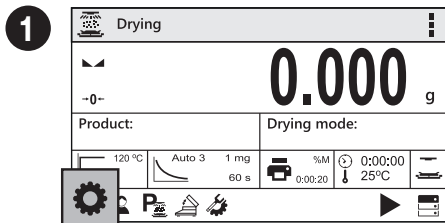
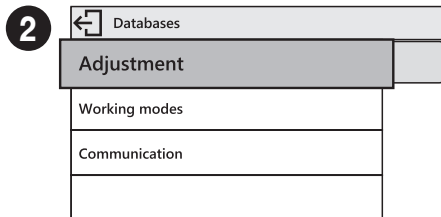
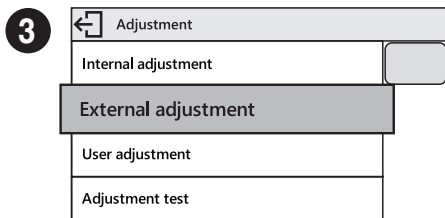
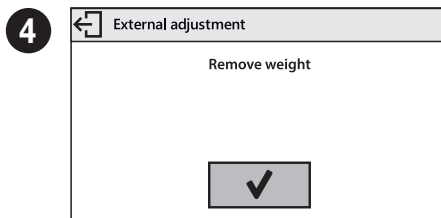
Make sure that the weighing pan is empty and press button for confirmation.

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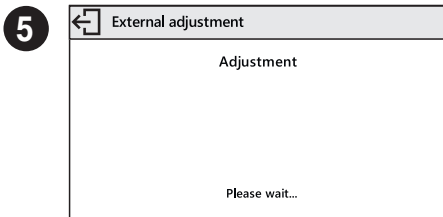
Wait for completion of the start mass determination process.

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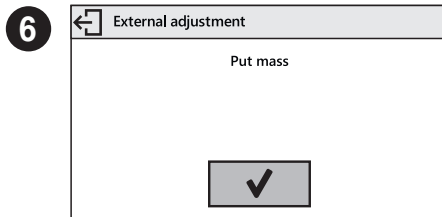
Load the weighing pan with mass standard and press button for confirmation.

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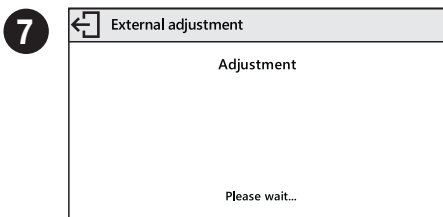
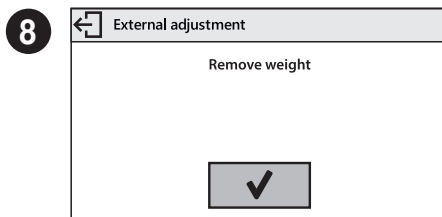
8.2. Zeroing

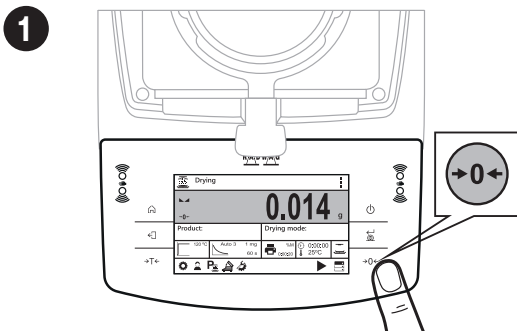Zerowanie

Make sure that the weighing pan is empty and press "Zeroing" button.

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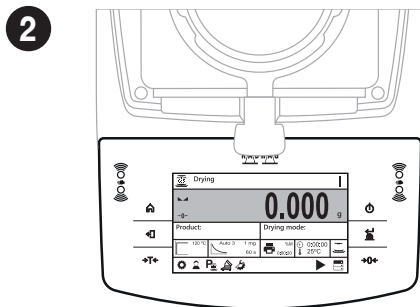
The moisture analyzer has been zeroed.

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
8.3. Taring

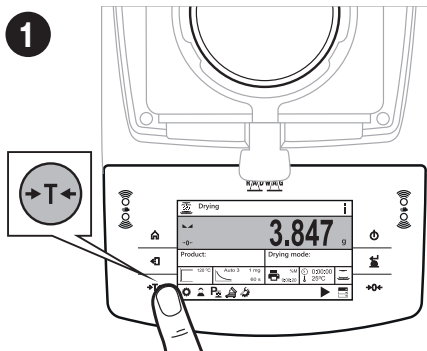Taring

With loaded weighing pan: upon stabilisation of the weighing result, press “Tare” button.

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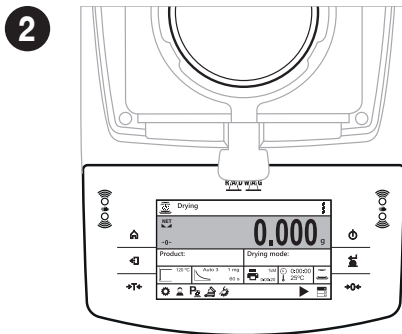
The moisture analyzer has been tared.

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8.4. Drying profiles / drying process finish modes

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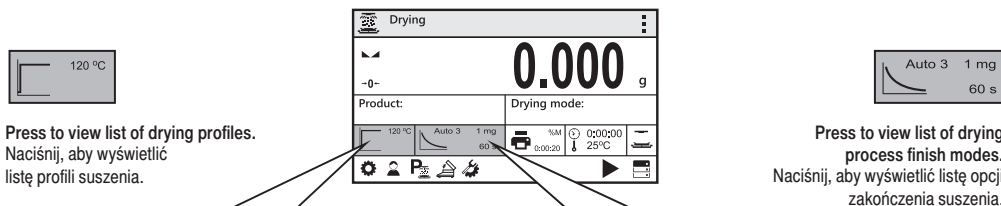Press to view list of drying profiles.

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Press to view list of drying process finish modes.

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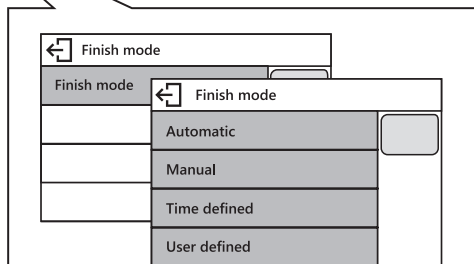
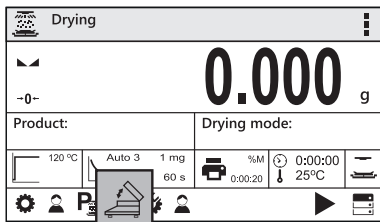
8.5. Automatically opened and closed drying chamber

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In order to open/close the drying chamber, press an on-screen button: "Open / Close the chamber"

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9. SETTINGS

USTAWIENIA

Some balance settings are accessible for Administrator exclusively.

Prior to balance parameters setup, log in as Administrator.

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
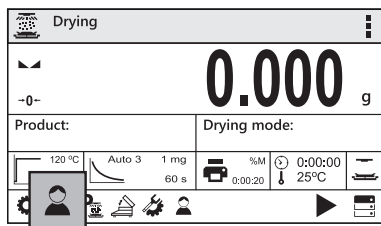
Press "Operators" button.

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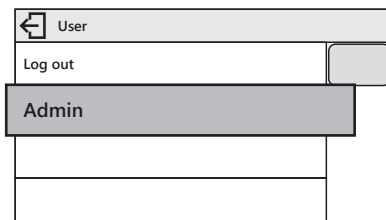
2

Users list is displayed, select "Admin" option.

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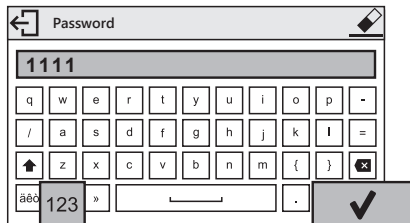3


Press **123** button to activate numeric keyboard. Enter

„1111” password and press **✓** button to confirm.

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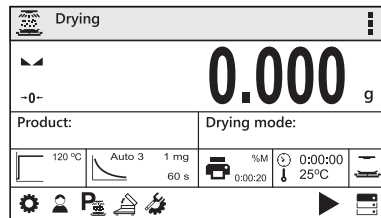4

You have logged as the Administrator.

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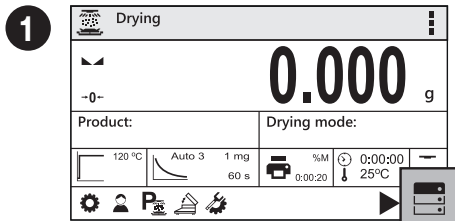
9.2. User settings

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
Prior to balance users setup, log in as Administrator (point 9.1).

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
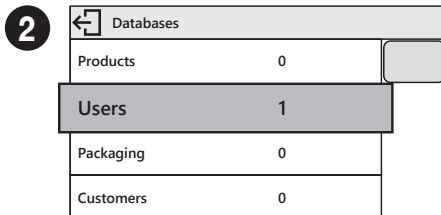
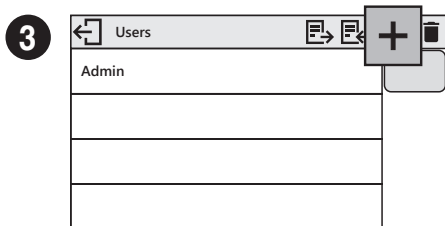
Press **+** button.

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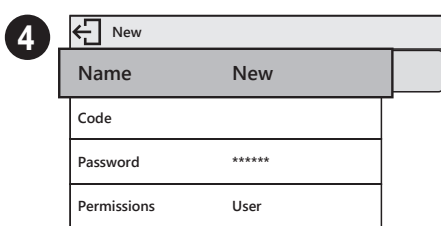
Select "Name" parameter.

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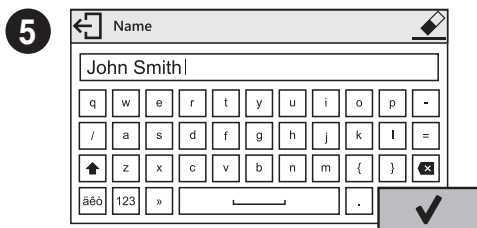Enter user name and press **✓** button to confirm.

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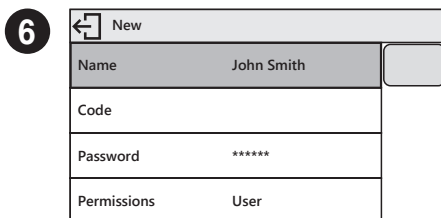7**Following the above procedure, set remaining parameters:**

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New	
Name	John Smith
Code	
Password	*****
Permissions	User

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**Prior to balance users setup, log in as Administrator (point 9.1.).**

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You can assign the proximity sensors with different functions for each working mode (the sensors will trigger different operation for each working mode). See the below procedure of sensors setup for the weighing mode.

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Press "Parameters" button.

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2

Parameters	
Adjustment	>
Working modes	
Communication	>
Peripherals	>

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3

Working modes	
Availability	>
Weighing	>
Drying	

Select "Weighing" parameter.

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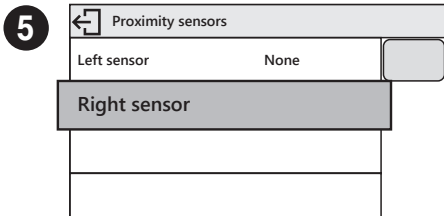
4


Weighing	
Readout	>
Proximity sensors	
Autotare	✓
Print mode	When stable

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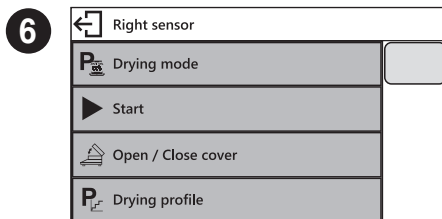10. SAMPLES AND DRYING METHODS

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10.1. Preparing samples

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Optimal sample structure and size are crucial features influencing measurements speed, accuracy and repeatability.

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Solids:

- sample crushing necessary,
- sample of low weight required,
- large sample surface.

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- large sample surface.

Ciecze:

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Some sample types require particular caution!

They may pose risk for people and objects. It is the moisture analyzer operator who is responsible for potential loss being a result of such samples use.

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Substances that in the course of heating may emit vapours causing corrosion (e.g. acids).

Apply small amount of samples, this helps to prevent risk of settling of vapour over cool housing components.

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Explosive and flammable and materials that in the course of heating emit flammable vapours.

Apply small amount of samples and lower drying temperature, this helps to prevent flames from rising and explosion. Use glasses and protective gloves.

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Toxic and corrosive substances and substances that in the course of heating emit toxic vapours.

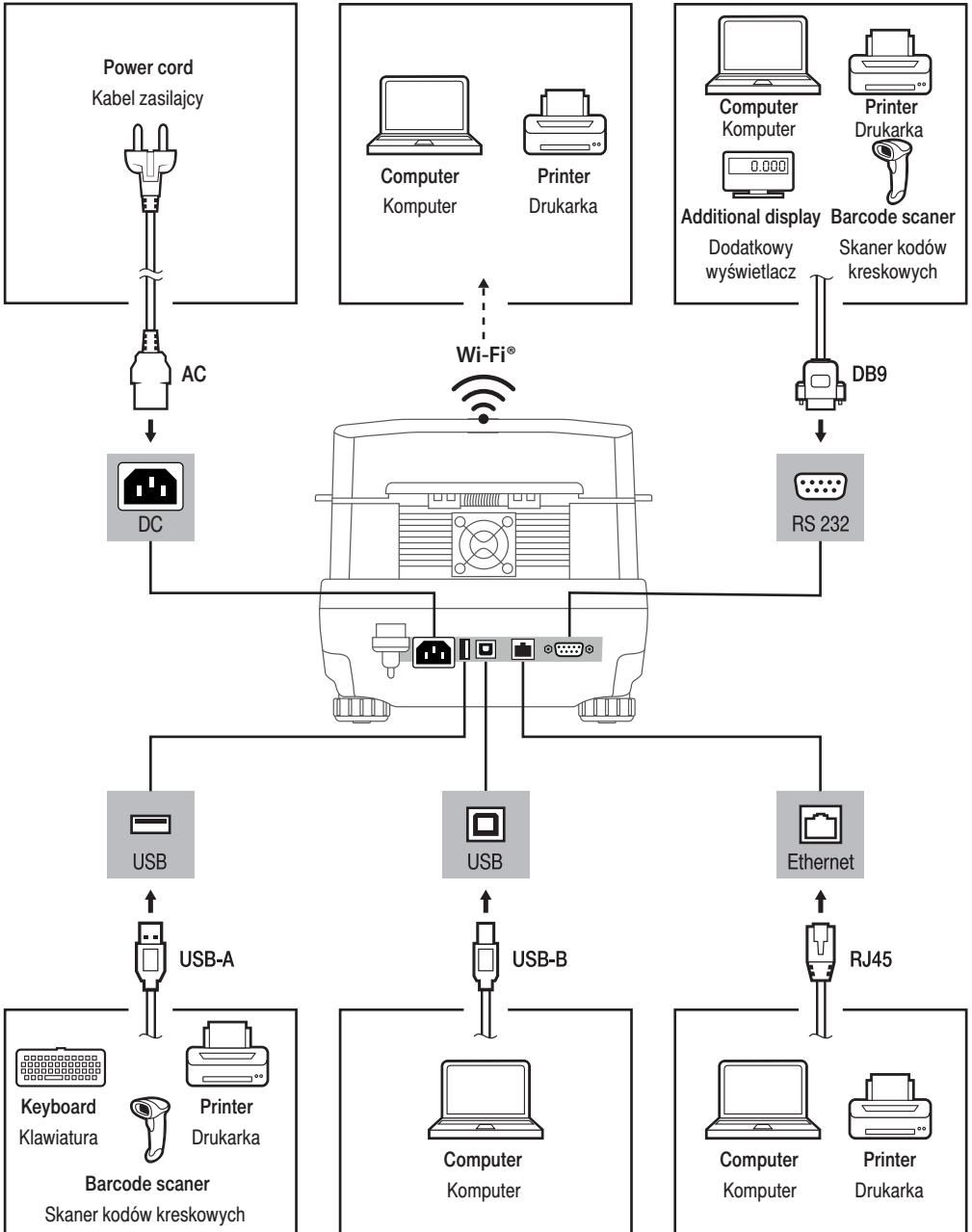
Due to risk of eyes, skin and lungs irritation, and due to risk of death, such samples must be dried in fume cupboard.

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Wi-Fi® is a registered trademark of Wi-Fi Alliance.

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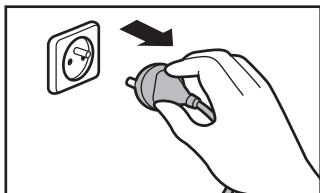
12. DEVICE CLEANING

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
12.1. General recommendations

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
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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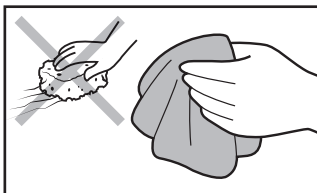Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface!

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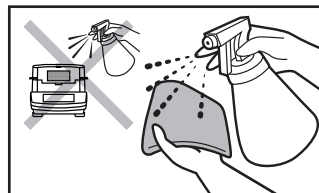Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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