IMMB-03-03-04-22-EN-PL

Start up Guide Szybki start

Analytical and Precision Balances

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Analytical balances:

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• AS R2 PLUS

Precision balances:

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- PS R1
- PS R1.M
- PS R2
- PS R2.M
- PS R2.H
- PS R2.M.H

To view a complete User Manual go to the website or scan QR code:

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3. RECOMMENDATIONS AND WARNINGS

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Operate the device in a room where the temperature ranges between 10–40 $^\circ C$ (50–104 $^\circ F)$ and where the relative humidity is below 80%.

Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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Avoid air drafts and air movements at the workstation.

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Make sure that the balance is placed on an even surface.

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Do not place the balance on an unstable ground exposed to shocks and vibrations.

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Do not place the balance in a hazardous area. Do not weigh explosive materials.

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4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Transport lock removal

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AS R2 Series

Keep the transport lock should a warranty claim or service be required.

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6.2. Leveling

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To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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7. PANEL AND SCREEN

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7.1. Panel keys

8.1.1. Adjustment (for models with an internal adjustment only)

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8.1.2. Adjustment (for models with an external adjustment only)

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It is recommended to carry out adjustment operation prior to running the WEIGHING mode. External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on device type and capacity.

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The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use mass standard of class F1 or F2. View the table and select the mass standard you need.

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Wait for completion of the start mass determination process.

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Wait for the adjustment completion.

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Load the weighing pan with mass standard and press $\left[\begin{array}{c} \textcircled{\bullet} \\ \hline \end{array} \right]$ key.

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Upon process completion unload the weighing pan and press $\left[\begin{array}{c} \textcircled{a} \\ \hline \end{array}
ight]$ key.

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8.2. Zeroing

Zerowanie

Make sure that the weighing pan is empty and press $\frac{1}{2 \text{ def}}$ key.

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RADWAG Net Mode + On/Off Cal Database TARE 3.24 Setup Function +T+ t © +0+ Delete Units Esc Insert

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To configure other working modes, follow the above procedure.

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The menu features the following working modes:

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- Weiahina
- Parts Counting •
- Checkweighing
- Dosing •
- Percent Weighing •
- Density Determination of Solids ٠
- Density Determination of Liquids •
- Statistics
- Totalising
- Peak Hold
- · Pipettes Calibration

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4

Detailed overview of the settings of the working modes is to be found in balance's main user manual.

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"Jednostki" i naciśnij przycisk

Using keys select unit for which accessibility is to be set and press key.

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8.7. Weighing unit selection Wybór jednostek

Wciśnij przycisk wy to select the weighing diff.

To configure other mass units, follow the above procedure.

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The balance enables selecting the following weighing units:

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2

- Troy ounce [ozt]Pennyweight [dwt]
- Newton [N]
- * Units available for verified balances exclusively

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Kod

Hasło

Using keys enter user name and press

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11. DEVICE CLEANING

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11.1. Guidelines and precautions

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Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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Use soft cloth made of either natural or synthetic fiber exclusively. Avoid using cloth made of stiff/hard fabric, this could cause scratches on the surface!

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Apply the cleanser onto the cloth. Mind not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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Shall it be necessary to disassemble the weighing chamber, follow the description provided in section 11.3.

- GLASS PANES AND COMPONENTS can be cleaned using window cleaner.
- WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).

- OPERATION PANEL AND HOUSING can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- Dry sample leftovers can be removed using either brush or handheld vacuum cleaner.
- Clean components can be dried using soft cloth or dust-free paper towel, with this the remaining moist will be absorbed.
- · Install all the components providing they are completely dry.

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11.3. Weighing chamber elements disassembly

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5

Take out the back glass. The housing has recesses to facilitate removing the glass. Wyciągnij tylną szybę do góry.

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