IMMB-07-02-05-22-EN-PL

Start up Guide

Szybki start

Precision Balances

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- WLC F1
- WLC C2

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

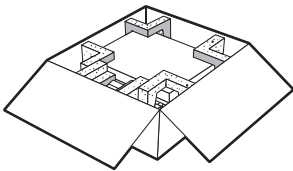
2

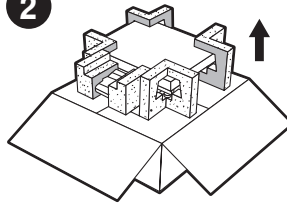3

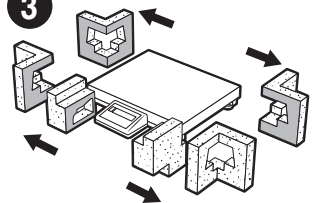
Keep the packaging should a warranty claim or service be required.

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3. RECOMMENDATIONS AND WARNINGS

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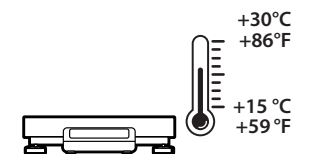
Operate the device in a room where the temperature ranges between 15–30 °C (59–86 °F) and where the relative humidity is below 80%.

Place the balance on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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
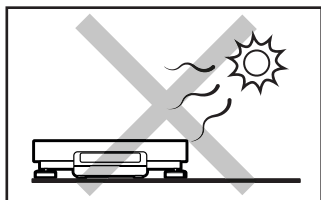
Place the balance away from heat sources. Avoid exposing the balance to the sunlight.

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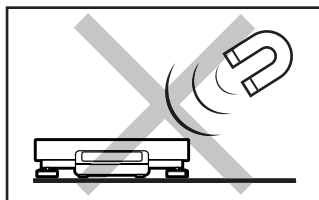
Avoid exposing the balance to a magnetic field. Do not weigh magnetic substances.

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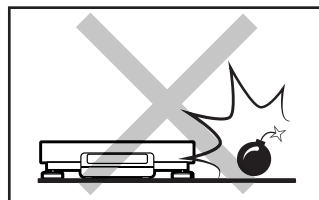
Do not place the balance in a hazardous area. Do not weigh explosive materials.

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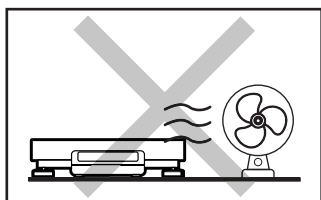
Avoid air drafts and air movements at the workstation.

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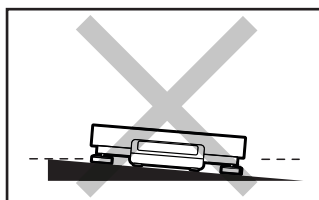Make sure that the balance is placed on an even surface.

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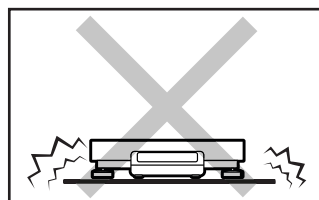Do not place the balance on an unstable ground exposed to shocks and vibrations.

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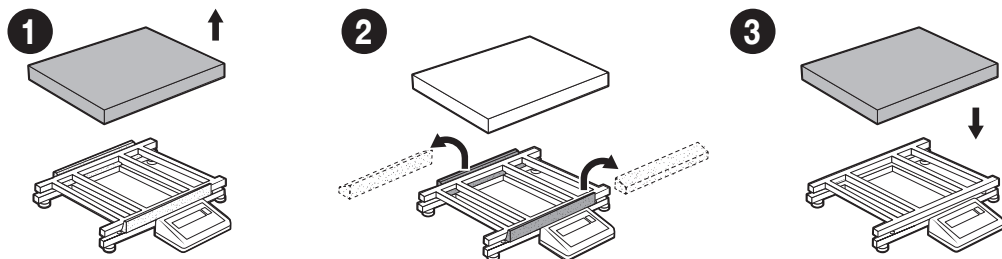
4. ACTIVITIES TO BE DONE PRIOR TO OPERATION

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4.1. Remove transport lock

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
5. PREPARING FOR WORK

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5.1. Temperature stabilization

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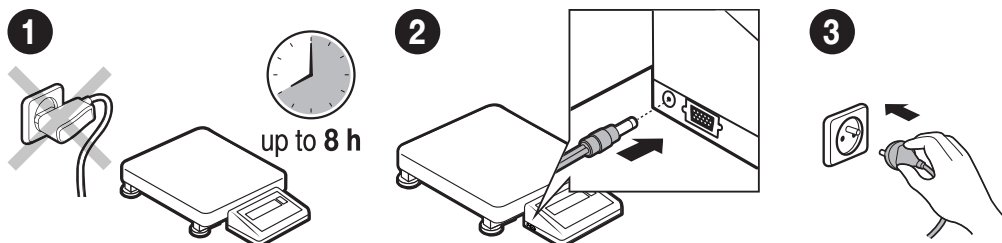Prior to switching the device on, let it reach room temperature (estimated stabilisation time: up to 8 hours).

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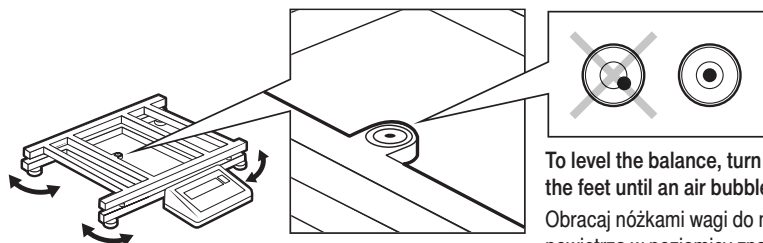
5.2. Levelling

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
To level the balance, turn its feet and keep turning the feet until an air bubble takes the central position.

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5.3. Switching the balance on

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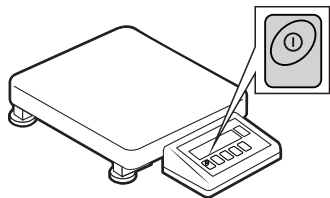Your balance is ready to use!

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To switch the device on, press ON / OFF key.

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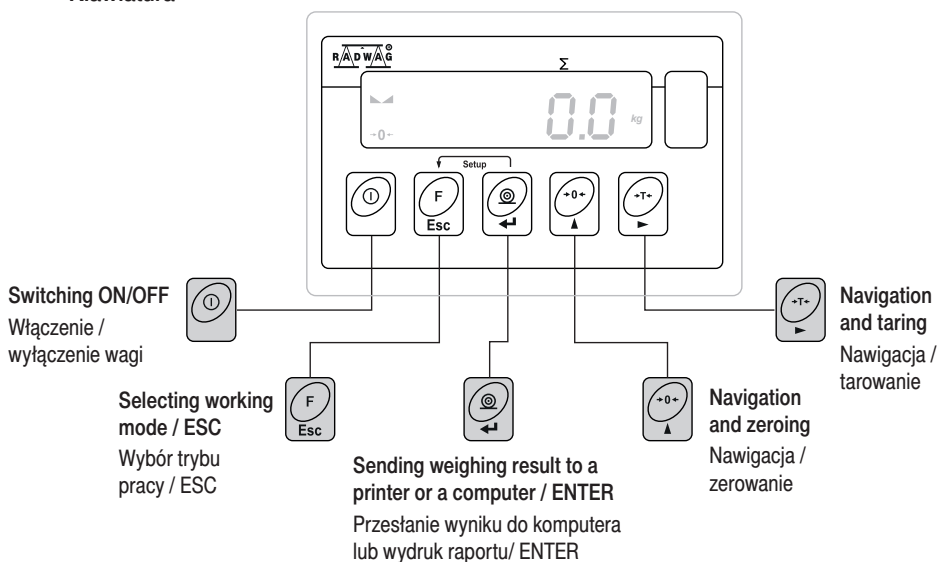
Stable result pictogram

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Balance zeroing pictogram

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Weighing result window and communication and information section

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7.1. Adjustment

Adiustacja

External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on balance type and capacity.

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
The adjustment mass standard does not come with the balance. Prior to adjustment prepare the right mass standard. You are recommended to use a mass standard of class F1 or F2. View the table and select the mass standard you need.

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
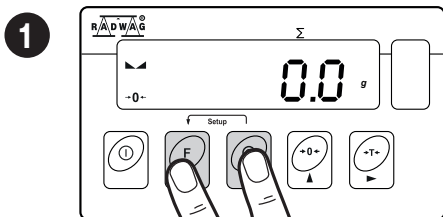
Simultaneously press **Esc F** + **→** keys to enter home screen.

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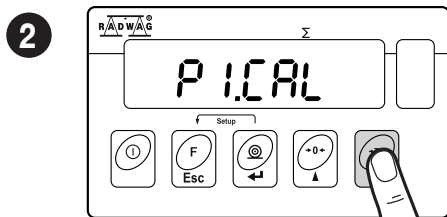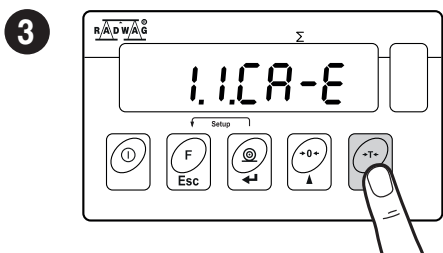
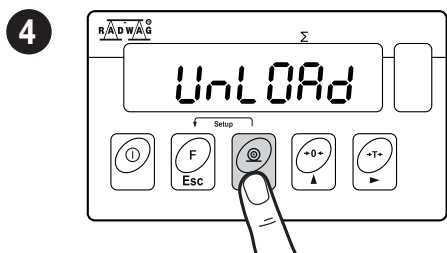
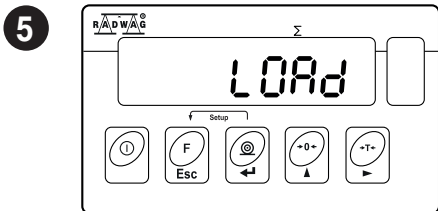
Unload the weighing pan and press **→** key.

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
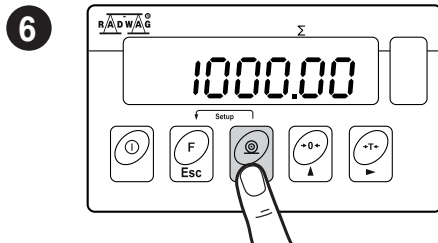
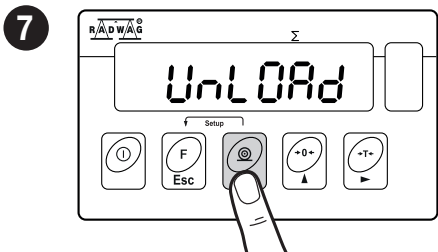
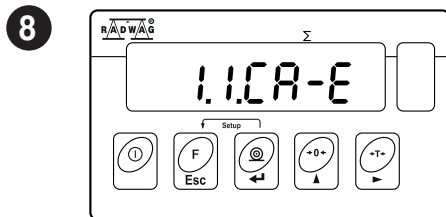
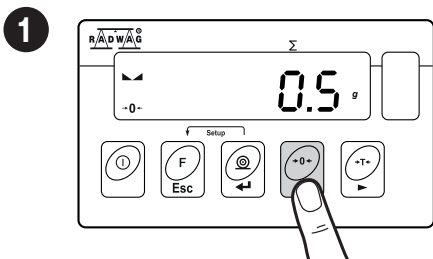
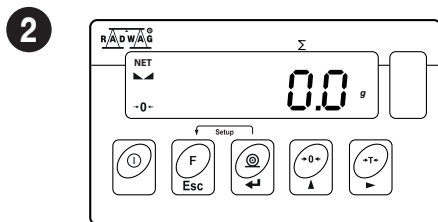
7.3. Taring

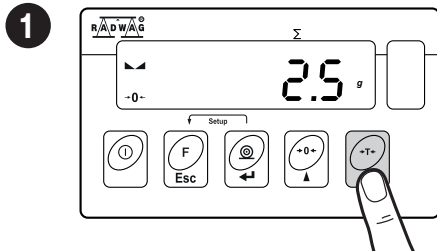Tarowanie

With loaded weighing pan: upon stabilisation of the weighing result, press **+T-** key.

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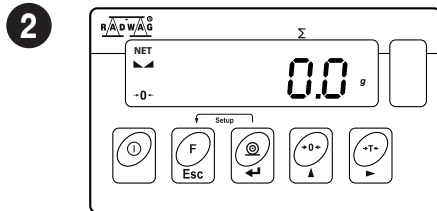
The balance has been tared.

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
7.4. Working modes accessibility

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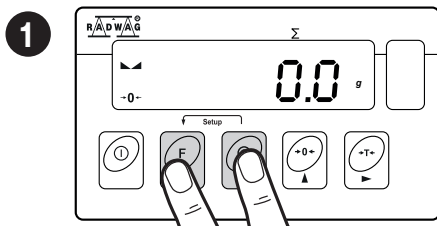
Simultaneously press **Esc F** + **↻** keys to enter home screen.

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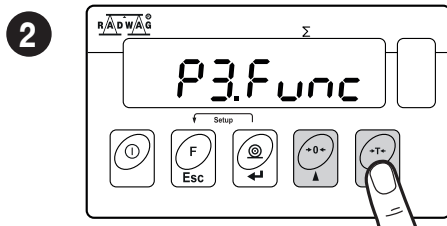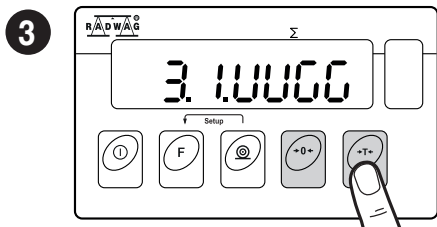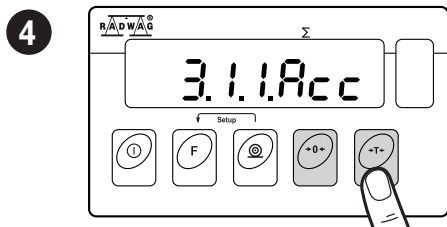
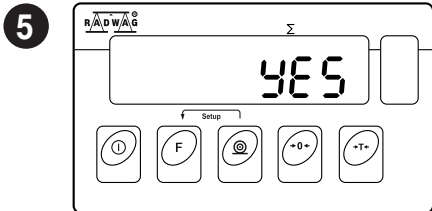
Using key select "YES" value and press key to confirm.

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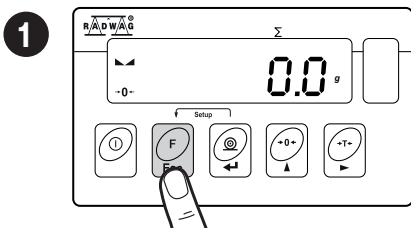
To configure other working modes, follow the above procedure.

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
Tryby pracy

Using key enter working modes menu.

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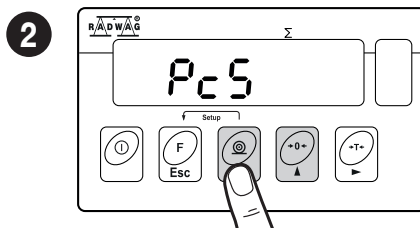Using key select working mode and press .

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The menu features the following working modes:

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Detailed overview of the settings of the working modes is to be found in balance's main user manual.

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

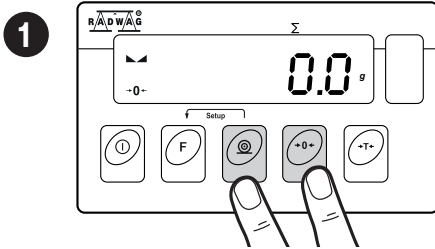
Simultaneously press + keys to enter home screen.

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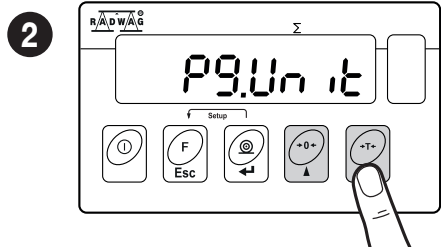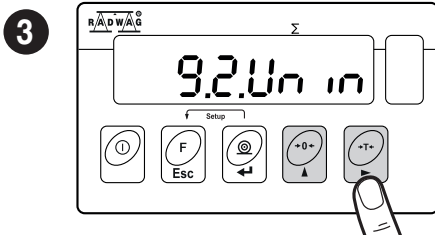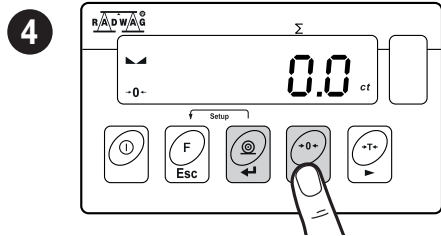
Using key select weighing unit and press key to confirm.

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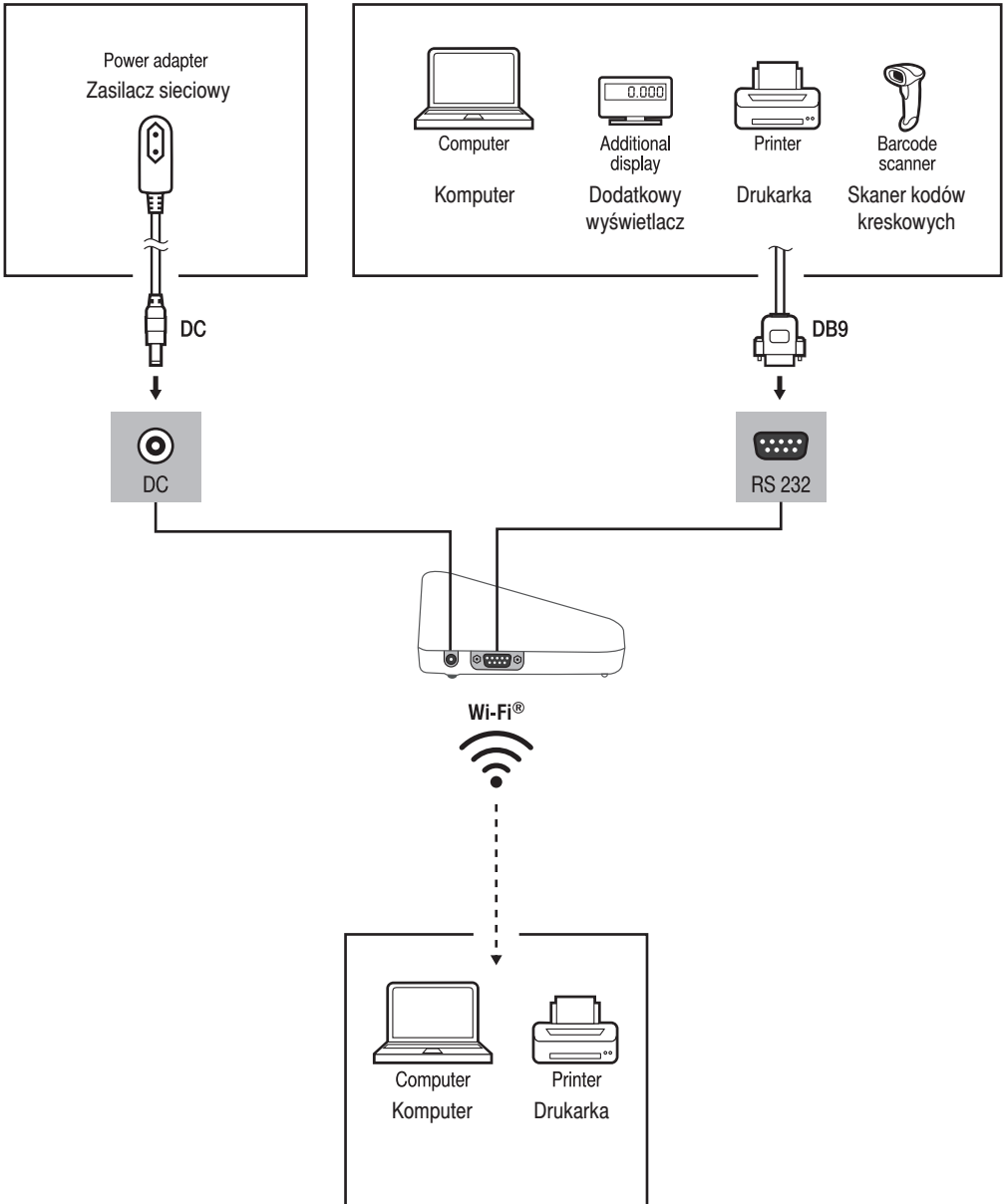
* (lb) pound unit unavailable for verified balances.

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Wi-Fi® is a registered trademark of Wi-Fi® Alliance.

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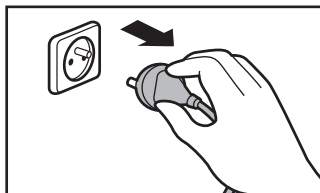
9. DEVICE CLEANING

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
9.1. General recommendations

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
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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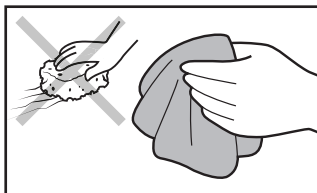Use soft cloth made of either natural or synthetic fiber exclusively. Avoid using cloth made of stiff/hard fabric, this could cause scratches on the surface!

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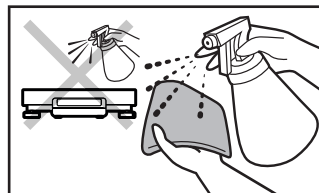Apply the cleanser onto the cloth. Mind not to apply the cleanser directly onto the device!

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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