IMMB-20-01-01-24-EN-PL

Start up Guide

Szybki start

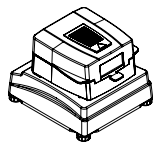
Moisture Analyzer

Wagosuszarka

- MA 5Y

1. CONTENT

ZAWARTOŚĆ

Moisture analyzer

Wagosuszarka

x1

Disposable pan

Szalka jednorazowa

x10

Cross-shaped holder

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x1

Drying pan handle

Uchwyt szalki

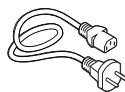
x1

Drying pan shield

Ostona szalki

x1

Power cord*

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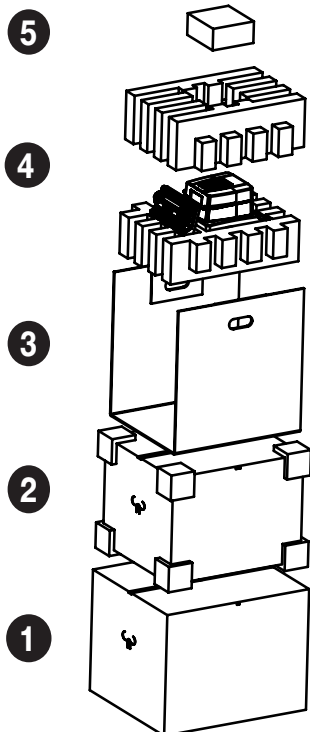
x1

**The plug type may vary by country.*

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2. UNPACKING

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
Keep the packaging should a warranty claim or service be required.

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3. RECOMMENDATIONS AND WARNINGS

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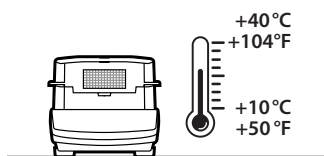
Operate the device in a room where the temperature ranges between 10–40 °C (50–104 °F) and where the relative humidity is below 80%.


Place the moisture analyzer on a solid surface to ensure stability. To obtain stable and repeatable weighing results, an anti-vibration table is recommended.

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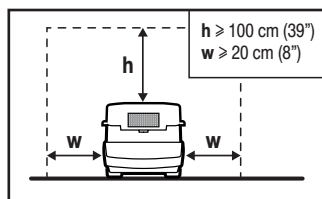
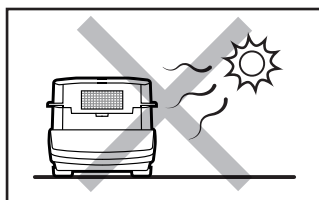
Place the device away from heat sources. Avoid exposing the device to the sunlight.

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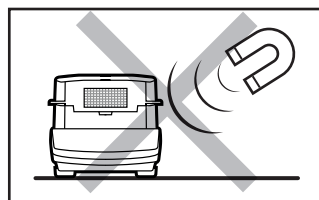
Avoid exposing the device to a magnetic field. Do not weigh magnetic substances.

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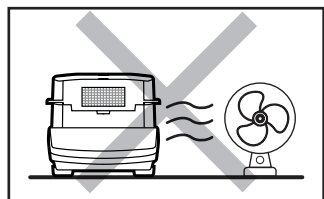
Avoid air drafts and air movements at the workstation.

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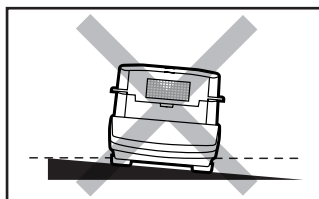
Make sure that the device is placed on an even surface.

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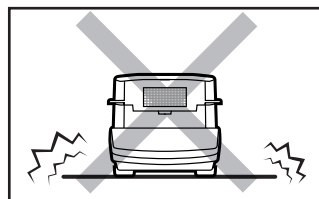
Do not place the device on an unstable surface exposed to shocks and vibrations.

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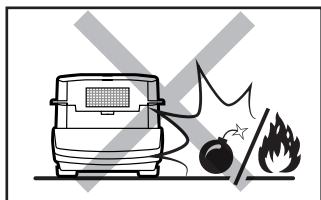
Do not place the device in a hazardous area. Do not weigh explosive or flammable materials.

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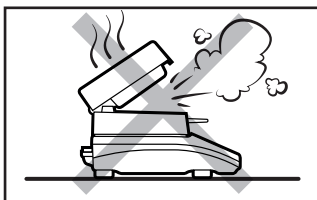Do not open in the course of operation since the heating components reach very high temperature.

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



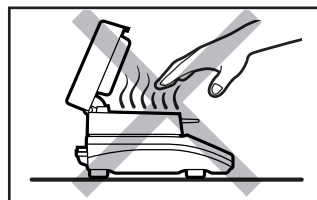
Be extremely careful while taking the sample out. The sample and the moisture analyzer elements may be hot.

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4. COMPONENTS ASSEMBLY

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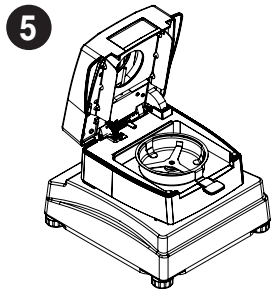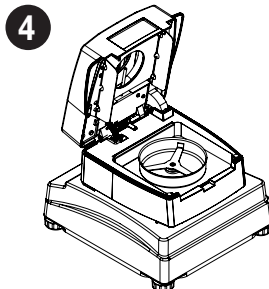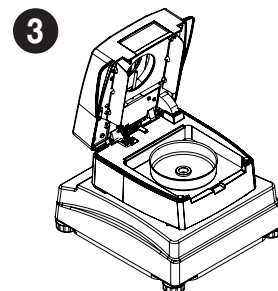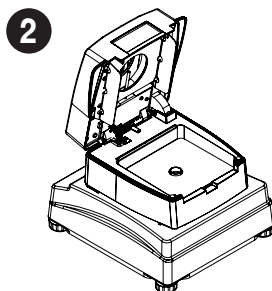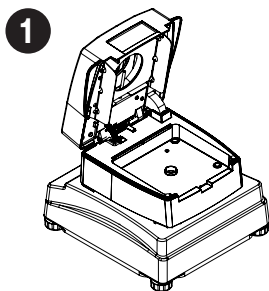5. PREPARING FOR WORK

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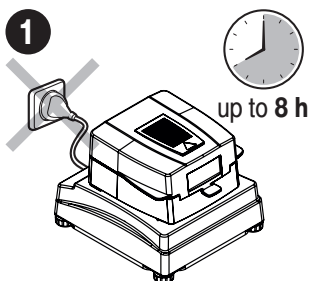
Place the device at the workstation, wait until it reaches room temperature.

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
5.1. Temperature stabilization

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
Prior to switching the device on let it reach room temperature (estimated stabilisation time: up to 8 hours).

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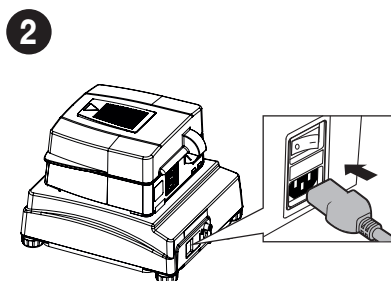Connect the power cord to the AC socket located at the back of the device.

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
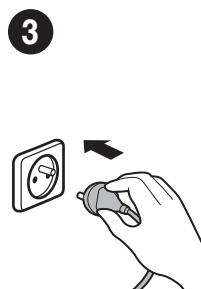
5.2. Switching the device on

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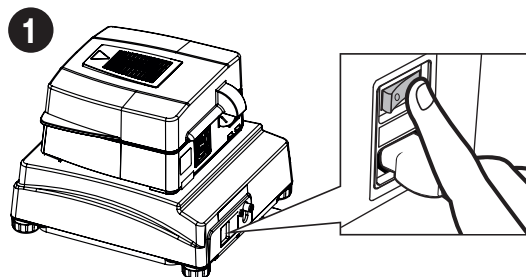
Press the switch located next to the power supply socket at the back of the housing to switch on the device.

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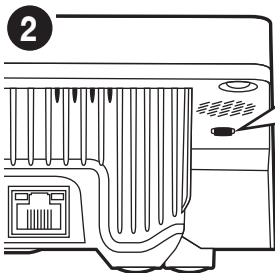
The balance switches on automatically when connected to the mains. If the balance is manually switched off using the top menu, it must be switched on using the button underneath the display housing on the left.

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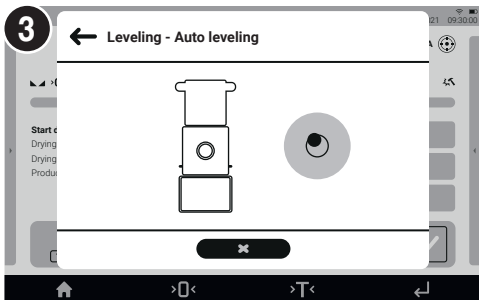The balance with automatic levelling starts the program and the levelling process runs automatically.

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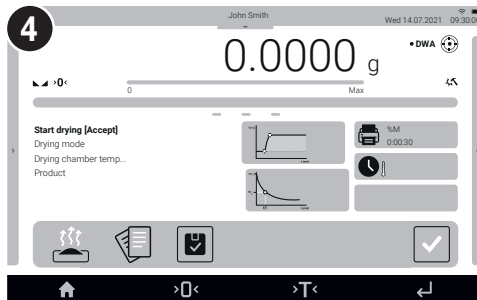When autolevelling is complete, the home screen is displayed.

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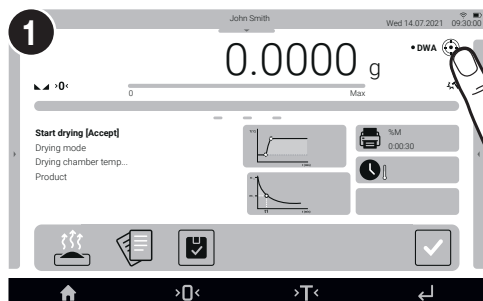
5.3 Manual Levelling

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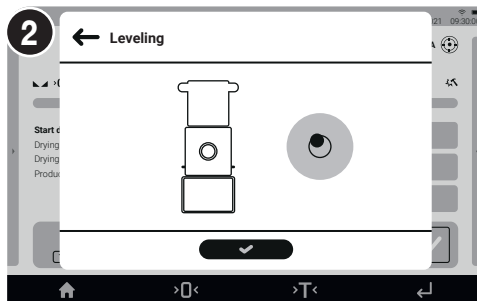
Ultra-microbalances and microbalances are equipped with the Reflex Level System, which monitors and automatically corrects the balance level.

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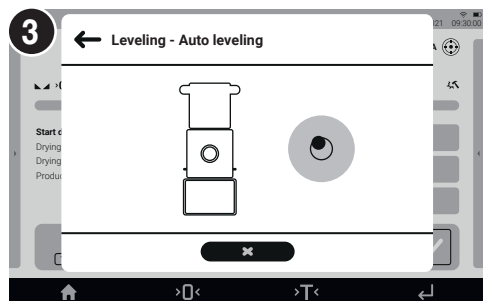
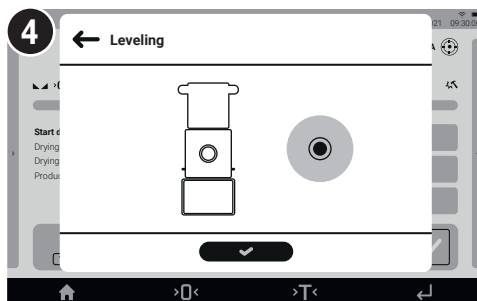
The moisture analyzer is ready to use!

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Before using WEIGHING mode, carry out instrument adjustment: refer to section 7.1.
(adjustment is not required for DRYING mode).

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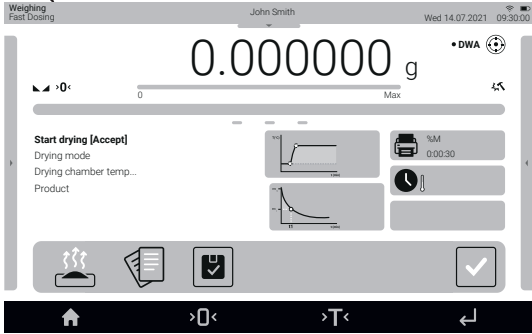
6. HOME SCREEN

EKRAN GŁÓWNY

information about the currently used working mode, logged-in operator, date, time, active connection with a computer.

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7. BASIC OPERATIONS

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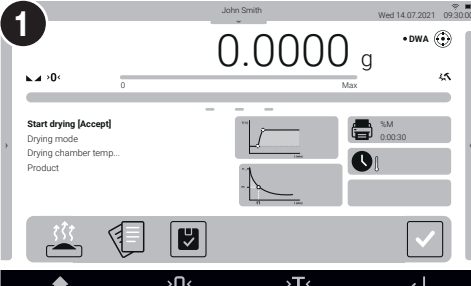
It is recommended to carry out adjustment operation prior to running the WEIGHING mode (adjustment is not required for DRYING). External adjustment is carried out using an external mass standard of the right accuracy and weight value, which value depends on device type and capacity.

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

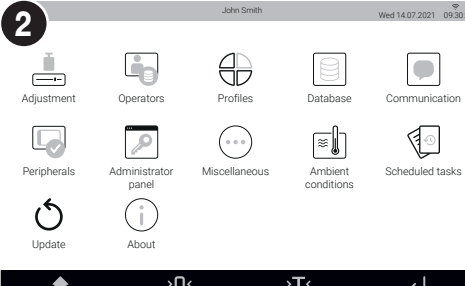
2

Press „Adjustment”.

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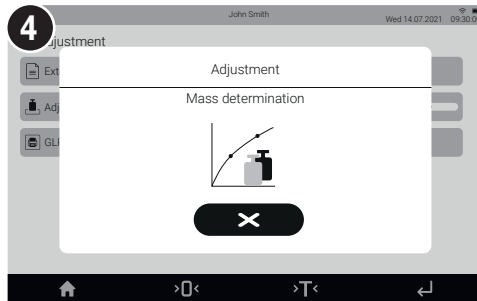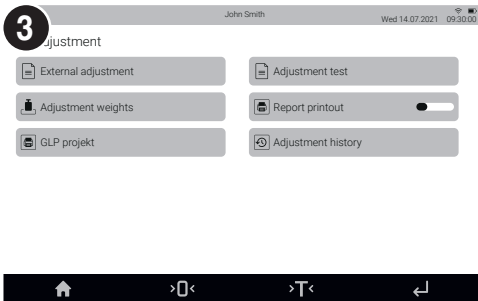Press „External adjustment” and follow the instructions on display.

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Wait for the process to be completed.

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The adjustment mass standard does not come with the moisture analyzer. Prior to adjustment prepare the right mass standard. You are recommended to use mass standard of class F1 or F2. View the table and select the mass standard you need. **Caution!** Make sure that the disposable weighing pan has been installed (read point 4.). Do not close the drying chamber in the course of adjustment.

Balance model Model wagi	Mass standard [g] Wzorzec masy [g]
PMV 50	50

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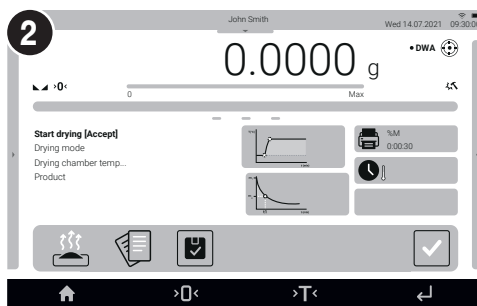
7.2. Zeroing

Zerowanie

Make sure that the weighing pan is empty and press “Zeroing” button.

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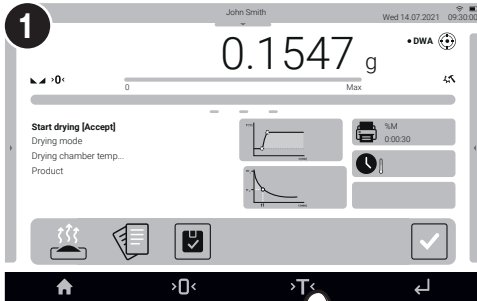
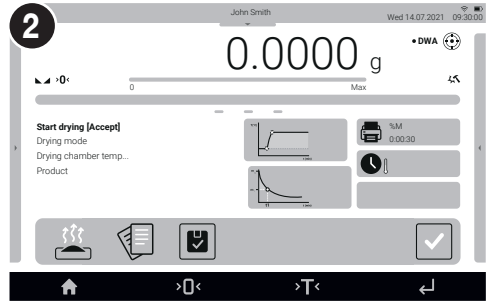
7.4. Drying profiles / drying process finish modes

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Prior to drying profiles / drying process finish modes setup, log in as Administrator (point 8.1.).

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
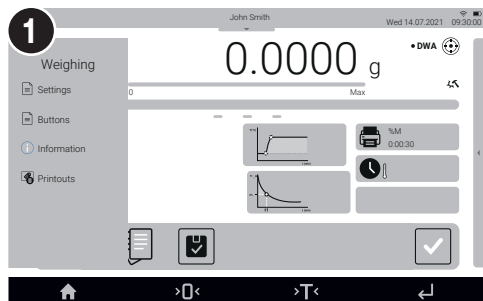
You can assign a function of your choice to each proximity sensor. In addition, you can assign various functions to the same sensor in different working modes. See the exemplary procedure for sensors configuration for the “Weighing” mode.

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Slide out the menu.

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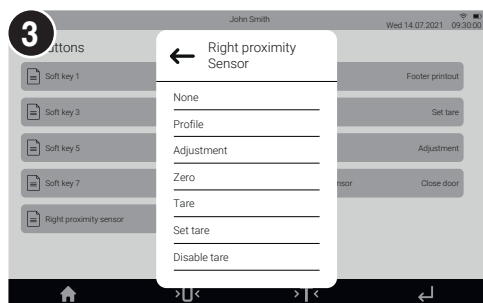
Select a respective sensor (right or left) from the list.

Select a function to be assigned to the sensor.

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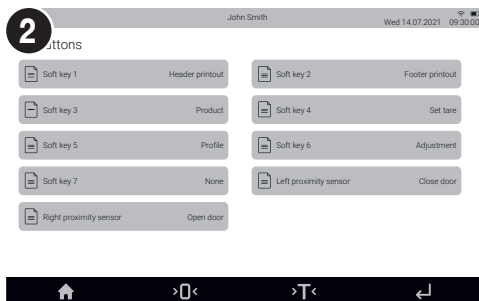9. SAMPLES AND DRYING METHODS

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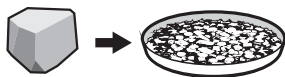
9.1. Preparing samples

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Optimal sample structure and size are crucial features influencing measurements speed, accuracy and repeatability.

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Solids:

- sample crushing necessary,
- sample of low weight required,
- large sample surface.

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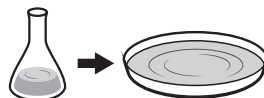
Some sample types require particular caution!

They may pose risk for people and objects. It is the moisture analyzer operator who is responsible for potential loss being a result of such samples use.

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Substances that in the course of heating may emit vapours causing corrosion (e.g. acids).

Apply small amount of samples, this helps to prevent risk of settling of vapour over cool housing components.

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Explosive and flammable and materials that in the course of heating emit flammable vapours.

Apply small amount of samples and lower drying temperature, this helps to prevent flames from rising and explosion. Use glasses and protective gloves.

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Toxic and corrosive substances and substances that in the course of heating emit toxic vapours.

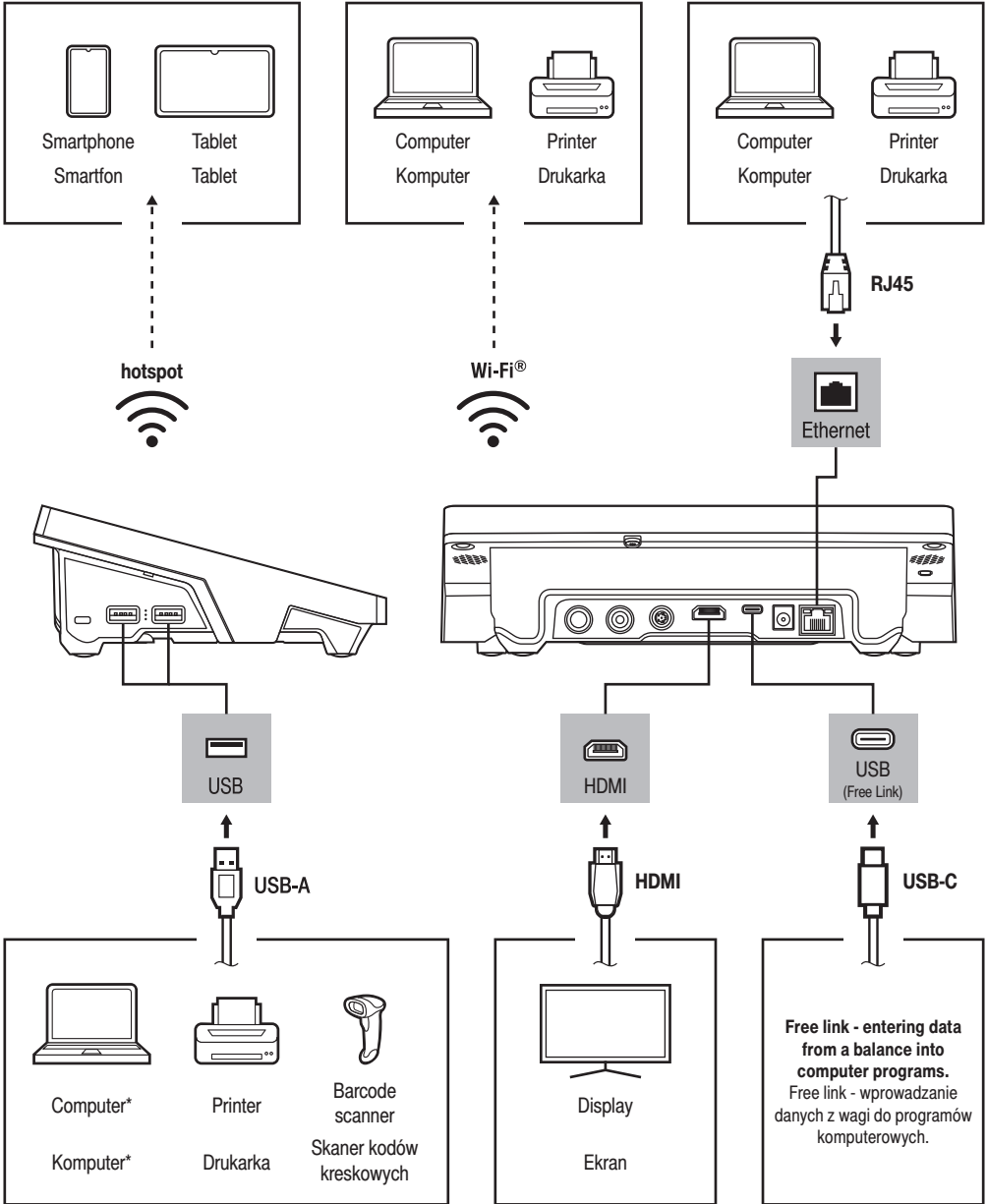
Due to risk of eyes, skin and lungs irritation, and due to risk of death, such samples must be dried in fume cupboard.

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Wi-Fi ® is a registered trademark of Wi-Fi Alliance.

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*A USB to RS232 converter must be used for proper connection.

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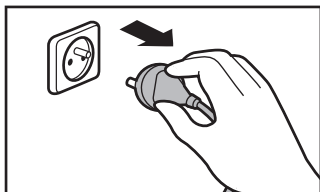
11. DEVICE CLEANING

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
11.1. General recommendations

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
Prior to cleaning it is necessary to disconnect the device from the mains! Remove the plug from the outlet, disconnect all cables.

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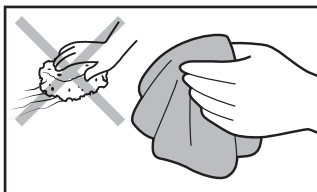Use soft cloths made of microfiber, natural fiber or man-made fiber. Avoid using abrasive cloths or cloths that might scratch the surface.

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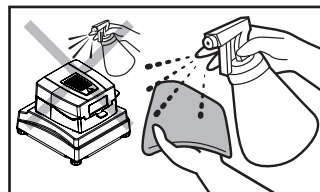Apply the cleanser onto the cloth first. Avoid applying the cleanser directly onto the device.

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Do not use aggressive cleaning agents (e.g. solvents, chlorine preparations, corrosive chemicals, bleach).

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Neither use substances of sharp structure nor intended for scrubbing.

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Make sure that neither dust nor liquid gets inside the weighing system (device inside).

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- **GLASS PANES AND COMPONENTS** can be cleaned using window cleaner.
- **WEIGHING PANS AND STAINLESS STEEL OR ALUMINIUM COMPONENTS** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).
- **OPERATION PANEL AND HOUSING** can be cleaned using a cloth slightly soaked in a solution of water and gentle cleanser (e.g. soap or dishwashing liquid).

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